



ETHANOL GAS STANDARD CYLINDER REPORT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59282 (08/2018)

Chemical Test Operator Name (Print) <i>Elmore, William</i>		
Location <i>BE PD</i>	Intoxilyzer® Serial Number <i>80-007096</i>	
Gas Lot Number <i>26021080A1</i>	Gas Cylinder Number <i>36</i>	Gas Expiration Date <i>10-5-2023</i>

Check When Done:

- 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 2. Perform an ACA Test (Level 1, Function C).
 Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 A. *0080AC*
 B. *0081AC*
 C. *0080AC*
- 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 B. ACA Test Record.

Chemical Test Operator Signature <i>William Elmore</i>	Date <i>11-8-21</i>
Reviewed By (Crime Laboratory Use Only) <i>Charles E. Ehl</i>	Date <i>11-8-2021</i>


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-007096
Location = BEPD 8164.16.00 09/18
11/08/2021 09:08

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	09:09
02 Std. Gas	0.080	09:09
03 Room Air	0.000	09:09
04 Std. Gas	0.081	09:10
05 Room Air	0.000	09:10
06 Std. Gas	0.080	09:11
07 Room Air	0.000	09:11

Lot No = 26021080A1
Cyl No = 36
Exp Date = 10/05/2023
County = 29 Oper No. = 130634



Operator Signature
WILLIAM ELMORE

Remarks:

Form 106-I8000