

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-006513

Location: MOTT
Hettinger CO SD

- Check When Done
- A. Is the breath tube heated? X
 - B. Is the diagnostic check complete? X
 - C. Is the time, date, and year correct (re-set if necessary)? X
 - D. Print test. (Attach test record.) X
 - E. Low AC. Use ≤ 0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5113 Lot No. 201507F Rep AC 0.020AC X
 - F. High AC. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR3378 Lot No. 16060 Rep AC 0.300AC X
 - G. Interferent check. Use 0.05% acetone in ethanol solution in ABA mode.
(Attach test record.)
Sim SN DR3803 Lot No. 1055 Solution 0.100AC + 0.05% Acetone X
 - H. RFI check. Run CMS mode. Key radio during CMS test.
(Attach test record.) X
 - I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.) 0.080AC

Lot No. 19615080AS Cyl No. 25 Exp. Date 5 Sep 17

Test 1 0.080 Test 4 0.080 Test 7 0.082

Test 2 0.080 Test 5 0.081 Test 8 0.081

Test 3 0.080 Test 6 0.081 Test 9 0.080 + Rew

Average 0.081

J. Tank pressure, Level 3 Func D Sub G P: 535 psi Reg 525 psi

K. Remarks/Maintenance record: Annual inspection.

30 May 17
Date

Roberta Mudge-Remmo
Field Inspector's Signature

31 May 17
Date

Deb Shanaver
Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:04

***** Printer Test *****

abcdefghijklmnopqrstuvwxy1234567890-=
ABCDEFGHIJKLMNopqrstuVwxyz!@#%\$^&*()_+?

abcdefghijklmnopqrstuvwxy1234567890-=
ABCDEFGHIJKLMNopqrstuVwxyz!@#%\$^&*()_+?

Current Instrument Setup

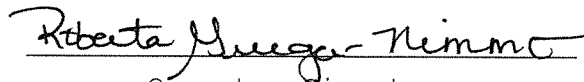
Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 19615080A5
Standard Cyl #? 25
Standard Expiration? 09/05/2017
Oper No? 888888

Flow Cal. Date: 08/20/2015
Slope 680
Intercept -696037

IR Calibration Date: 08/21/2015

	3um	9um
0th Coef(*100):	-14436	-13610
1st Coef(*100):	265490	136147
2nd Coef(*100):	4422	1710
H2O adj(mg/l*10k):	672	448

***** Printer Test End *****


Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: *Print Test*

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

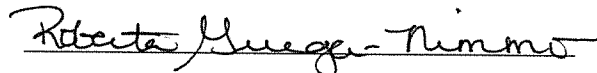
CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:29

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:30
02 Std. Sol.	0.021	08:31
03 Room Air	0.000	08:31
04 Std. Sol.	0.021	08:32
05 Room Air	0.000	08:32
06 Std. Sol.	0.021	08:33
07 Room Air	0.000	08:34

08 Sim Temp = 34.0°C

Simul Ser No = DR5113
Std Sol No = 201507F
County = 21 Oper No. = 888888


Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: low AC
0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:42

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:42
02 Std. Sol.	0.306	08:43
03 Room Air	0.000	08:44
04 Std. Sol.	0.305	08:45
05 Room Air	0.000	08:45
06 Std. Sol.	0.306	08:46
07 Room Air	0.000	08:46

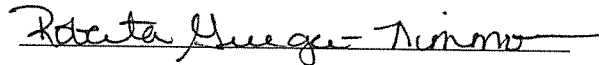
08 Sim Temp = 34.0°C

Simul Ser No = DR3378

Std Sol No = 16060

County = 21

Oper No. = 888888



Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: High AC
0.300 AC

Form 106-I8000

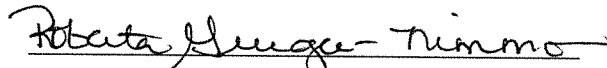
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:50

Test	AC	Time
01 Room Air	0.000	08:51
02 *Subject Test	INT*	08:51
03 Room Air	0.000	08:52

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS821456
Lot No = 19615080A5
Cyl No = 25
Expiration Date = 09/05/2017
County = 21 Oper No. = 888888



Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Interferent Check
0.100AC + 0.05% Acetone

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

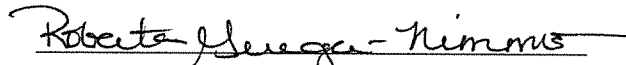
CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:24

Test	AC	Time
01 Diagnostic	OK	08:25
02 Room Air	0.000	08:25
03 Subject Test 1	0.000	08:26
04 Room Air	RFI*	08:27
05 Room Air	0.000	08:27

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = RFI CHECK
Dr. Lic. = ND/DIS821456
Lot No = 19615080A5
Cyl No = 25
Expiration Date = 09/05/2017
County = 21 Oper No. = 888888

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.


Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: RFI Check

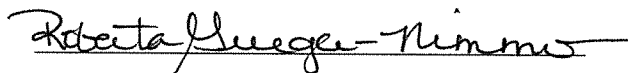
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:04

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:05
02 Std. Gas	0.080	08:05
03 Room Air	0.000	08:06
04 Std. Gas	0.080	08:06
05 Room Air	0.000	08:07
06 Std. Gas	0.080	08:07
07 Room Air	0.000	08:08

Lot No = 19615080A5
Cyl No = 25
Exp Date = 09/05/2017
County = 21 Oper No. = 888888



Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check
0.080 AC

Form 106-I8000

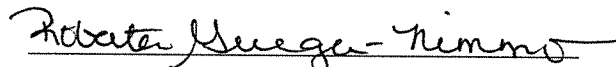
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:12

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:13
02 Std. Gas	0.080	08:13
03 Room Air	0.000	08:13
04 Std. Gas	0.081	08:14
05 Room Air	0.000	08:14
06 Std. Gas	0.081	08:15
07 Room Air	0.000	08:15

Lot No = 19615080A5
Cyl No = 25
Exp Date = 09/05/2017
County = 21 Oper No. = 888888


Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check

0.080 AC

Form 106-I8000

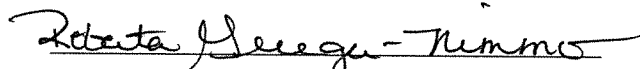
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006513
Location = MOTT 8164.13.00 06/09
05/30/2017 08:17

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:18
02 Std. Gas	0.082	08:18
03 Room Air	0.000	08:19
04 Std. Gas	0.081	08:19
05 Room Air	0.000	08:20
06 Std. Gas	0.080	08:20
07 Room Air	0.000	08:20

Lot No = 19615080A5
Cyl No = 25
Exp Date = 09/05/2017
County = 21 Oper No. = 888888


Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks:

Calibration Check
0.080 AC

Form 106-I8000