



**INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT**  
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL  
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM  
 SFN 59281 (06/2018)

Serial Number <i>80-006682</i>	Instrument Location <i>BCCC</i>
Reason for Install/Repair <input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
  - A. Print Test (Level 1, Function P).
  - B. ACA Test (Level 1, Function C).
  - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): \_\_\_\_\_
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>[Signature]</i>	Date <i>5/13/2020</i>
---	--------------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By <i>[Signature]</i>	Certified Date <i>5/14/2020</i>
---	------------------------------------

Intoxilyzer Test Record and Checklist  
 NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer          Alcohol Analyzer  
 North Dakota Model 8000      SN 80-006682  
 Location = BCCC                  8164.14.00 09/16  
 05/13/2020                          20:43

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=  
 ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=  
 ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#%^&\*()\_+?

Current Instrument Setup


Data Entry Mode:                  Enabled  
 Start Test Sequence:              DABACABA  
 Display Prelim Rslt?              Yes  
 Display Third Digit?              Yes  
 Inhib Printer(Y/N)?              No  
 Display Volume?                  No  
 Disable On Memfull?              Yes  
 # of Print Copies?                1  
 Select Std (D/W/I)?              Dry  
 Standard Value?                  0.080  
 Standard Lot #?                  05620080A1  
 Standard Cyl #?                  47  
 Standard Expiration?              04/05/2022  
 Oper No?                          100012  
  
 Flow Cal. Date:                  09/06/2016  
 Slope                              701  
 Intercept                         -706193

IR Calibration Date:              09/06/2016  
    3um                              9um  
 -----  
 0th Coef(\*100):                  -17150                              -23247  
 1st Coef(\*100):                  271726                              134617  
 2nd Coef(\*100):                  831                                  1215  
 H2O adj(mg/l\*10k):              393                                  518

\*\*\*\*\* Printer Test End \*\*\*\*\*

*CYL #47 USED AT  
 CRIME LAB.  
 FIELD INSPECTOR  
 SCANNED IN GAS  
 CYLINDER #45  
 ASSIGNED TO BCCC  
 BEFORE RUNNING  
 ACA & RFI TESTS.*

*GOOD INSTALL.  
 Charles E. Ehl  
 5/14/2020*

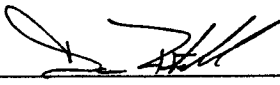  
 \_\_\_\_\_  
 Operator Signature  
 DANA RUSTEBAKKE

Remarks:

Form 106-I8000


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-006682  
Location = BCCC      8164.14.00 09/16  
05/13/2020      20:48

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	20:49
02 Std. Gas	0.082	20:49
03 Room Air	0.000	20:50
04 Std. Gas	0.081	20:50
05 Room Air	0.000	20:51
06 Std. Gas	0.082	20:51
07 Room Air	0.000	20:52

Lot No = 05620080A1  
Cyl No = 45  
Exp Date = 04/05/2022  
County = 02      Oper No. = 100012

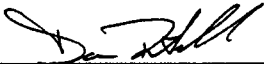

Operator Signature  
DANA RUSTEBAKKE

Remarks:

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-006682  
Location = BCCC      8164.14.00 09/16  
05/13/2020      20:52

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	20:53
02 Std. Gas	0.082	20:53
03 Room Air	RFI*	20:53
04 Room Air	0.000	20:54

\*Invalid Test  
Inhibited - RFI

Lot No = 05620080A1  
Cyl No = 45  
Exp Date = 04/05/2022  
County = 02      Oper No. = 100012



Operator Signature  
DANA RUSTEBAKKE

Remarks:

Form 106-I8000

*Install after Annual Ins.*