Toxicology Section/Breath Alcohol Program Intoxilyzer® 8000 Inspection

## INTOXILYZER® 8000 INSPECTION

Intoxilyzer<sup>®</sup> 8000 Serial Number: 80-006687

Location: TOXL

- Α. General Setup and Checks:
  - X Diagnostic Tests Pass and Instrument in Ready Mode 1.
  - Breath tube heated 2.
  - Date, time and location code (Level 2,E). Re-set if necessary. 3.
  - Print test (Level 1,P). Sign and attach test record. 4.
  - X Tank monitor (Level 3,D,G). Display and Regulator within 50 psi. 5.
    - Display: 100 A psi Regulator: 1000 psi a.
- Β. Tests (Sign and attach test records):

 $\mathbf{K}$  Configure simulator for the following tests (Level 1,S).

- Wet Calibration Check 1.
  - а.
  - A Low AC (Level 1,C): Known Value  $\leq 0.03$  AC: 0.030 AC Sim. Ser #: MP5289 Lot #:201911E Exp. Date: 1.19.21W High AC (Level 1,C): Known Value  $\geq 0.25$  AC: 0.4900 AC Sim. Ser #: MP3062 Lot #:202103E Exp. Date: 3.29.29.23b.

Configure dry gas standard for the following tests (Level 1,S).

- X Interferent Check (Level 1,B):\_Known Value: 0.10 AC + 0.05% Acetone 2. Sim. Ser #: DR3877Lot #: ICS 6 Exp. Date: NA Display reads "Interferent Detect"
- Display reads "RFI Detect" RFI Check (CMS Mode) 3.
- Dry Calibration Check (Level 1,C): Known Value 0.080 AC 4. Gas Cyl Lot #:05620080 A1 Cyl #: 48 Exp. Date: 4.5.22 Test 10.082Test 40.082Test 70.082Test 20.081Test 50.082Test 80.082Test 30.082Test 60.083Test 90.082Average 0.082
- Remarks/Maintenance: ANNUAL JUSPECTION C.

Instrument is acceptable to be used in the field.

Analyst Signature

Reviewed by

02/18 Issuing Authority - CEE

<u>5.5-202</u> Date 5-11-21

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501 CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 SN 80-006687 Location = TOXL 8164.14.00 09/16 05/05/2021 10:35 abcdefghijklmnopqrstuvwxyz1234567890-= ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#\$%^&\*() +? abcdefghijklmnopqrstuvwxyz1234567890-=| ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#\$%^&\*() +? Current Instrument Setup Data Entry Mode: Enabled Start Test Sequence: DABACABA Display Prelim Rslt? Yes Display Third Digit? Yes Inhib Printer(Y/N)? No Display Volume? No Disable On Memfull? Yes # of Print Copies? 1 Select Std (D/W/I)? Dry Standard Value? 0.080 Standard Lot #? 02621080A1 Standard Cyl #? 28 Standard Expiration? 03/05/2023 Oper No? 666666 Flow Cal. Date: 10/05/2018 Slope 696 -609977 Intercept IR Calibration Date: 10/05/2018 3um 9um 0th Coef(\*100): -11478 -18771 255256 4432 1st Coef(\*100): 132135 2nd Coef(\*100): 1676 H2O adj(mg/l\*10k): 604575 

Operator Signature CHARLES EDER

Remarks: PRINT TEST

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. IntoxilyzerAlcohol AnalyzerNorth Dakota Model 8000SN 80-006687Location = TOXL8164.14.00 09/1605/05/202111:00

	WET CAL CHECK	
Test	AC	Time
01 Room Air 02 Std. Sol. 03 Room Air 04 Std. Sol. 05 Room Air 06 Std. Sol. 07 Room Air	0.000 0.029 0.000 0.029 0.000 0.030 0.030	11:00 11:01 11:02 11:02 11:03 11:04 11:04

 $08 \text{ Sim Temp} = 34.0^{\circ}\text{C}$ 

Simul Ser No = MP5289 Std Sol No = 201911E County = 08

Oper No. = 666666

Uð 6

Operator Signature CHARLES EDER

Low AC 0.030AC

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. IntoxilyzerAlcohol AnalyzerNorth Dakota Model 8000SN 80-006687Location = TOXL8164.14.00 09/1605/05/202111:05

	WET CAL CHECK	]
Test	AC	Time
01 Room Air	0.000	11:06
02 Std. Sol.	0.395	11:07
03 Room Air	0.000	11:07
04 Std. Sol.	0.395	11:08
05 Room Air	0.000	11:09
06 Std. Sol.	0.397	11:09
07 Room Air	0.000	11:10

08 Sim Temp = 34.0°C

Simul Ser No = MP3062 Std Sol No = 2022103E County = 08

Oper No. = 666666

LARL Cat

Operator Signature CHARLES EDER

HIGH AC 0.400 AC

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501 CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 SN 80-006687 Location = TOXL 8164.14.00 09/16 05/05/2021 11:15 Test AC Time 01 Room Air 0.000 11:16 02 \*Subject Test INT\* 11:17 03 Room Air 0.000 11:17 \*Invalid Test Interferent Detected Sub Name = DISCOVER, THE SPIRIT Sub DOB = 01/08/1987Sub Sex = Male Weight = 150Test = DUI Cit = INTERFERENT Dr. Lic. = ND/DIS871456Lot No = 05620080A1Cyl No = 48Expiration Date = 04/05/2022County = 08Oper No. = 666666

Operator Signature CHARLES EDER

c + 0.05% ACETONE ler 0.10AC +

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501 CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 SN 80-006687 Location = TOXL 8164.14.00 09/16 05/05/2021 11:18 Test AC Time 01 Diagnostic OK 11:19 02 Room Air 0.000 11:20 03 \*Subject Test RFI\* 11:20 04 Room Air 0.000 11:20 \*Invalid Test Inhibited - RFI Sub Name = DISCOVER, THE SPIRIT Sub DOB = 01/08/1987Sub Sex = Male Weight = 150Test = DUI Cit = RFI CHECK Dr. Lic. = ND/DIS871456Lot No = 05620080A1Cyl No = 48Expiration Date = 04/05/2022County = 08Oper No. = 666666

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

Operator Signature CHARLES EDER

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 SN 80-006687 Location = TOXL 8164.14.00 09/16 05/05/2021 11:22

	DRY CAL CHECK	
Test	AC	Time
01 Room Air 02 Std. Gas 03 Room Air 04 Std. Gas 05 Room Air 06 Std. Gas 07 Room Air	0.000 0.082 0.000 0.081 0.000 0.082 0.000	11:22 11:23 11:23 11:23 11:24 11:24 11:24 11:25

Lot No = 05620080A1Cyl No = 48Exp Date = 04/05/2022County = 08

Oper No. = 666666

Operator Signature

CHARLES EDER ALIBRATION CHECK 0.080 AC

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 SN 80-006687 Location = TOXL 8164.14.00 09/16 05/05/2021 11:25

	DRY CAL CHECK	
Test	AC	Time
01 Room Air 02 Std. Gas 03 Room Air 04 Std. Gas 05 Room Air 06 Std. Gas 07 Room Air	0.000 0.082 0.000 0.082 0.000 0.083 0.000	11:26 11:26 11:27 11:27 11:27 11:28 11:28

Lot No = 05620080A1Cyl No = 48Exp Date = 04/05/2022County = 08

Oper No. = 666666

Ó

Operator Signature CHARLES EDER

AECK ALIBRATION 0.080 AC

Form 106-18000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. IntoxilyzerAlcohol AnalyzerNorth Dakota Model 8000SN 80-006687Location = TOXL8164.14.00 09/1605/05/202111:29

	DRY CAL CHECK	
Test	AC	Time
01 Room Air 02 Std. Gas 03 Room Air 04 Std. Gas 05 Room Air 06 Std. Gas 07 Room Air	0.000 0.082 0.000 0.082 0.000 0.082 0.000	11:29 11:29 11:30 11:30 11:31 11:31 11:32

Lot No = 05620080A1 Cyl No = 48 Exp Date = 04/05/2022 County = 08

Oper No. = 666666

Operator Signature CHARLES EDER CALIBRATTON CHECK

0,080 AC

Form 106-18000