



ETHANOL GAS STANDARD CYLINDER REPORT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59282 (08/2018)

Chemical Test Operator Name (Print) <i>Shawn Holm</i>		
Location <i>Divide Co So Intake / Processing</i>	Intoxilyzer® Serial Number <i>80-00161682</i>	
Gas Lot Number <i>26021 080A1</i>	Gas Cylinder Number <i>024</i>	Gas Expiration Date <i>10/5/2023</i>

Check When Done:

- 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 2. Perform an ACA Test (Level 1, Function C).
 Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 A. 0.079 AC
 B. 0.080 AC
 C. 0.079 AC
- 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

Chemical Test Operator Signature <i>Shawn Holm</i>	Date <i>08/19/22</i>
Reviewed By (Crime Laboratory Use Only) <i>Janelle Putscheller</i>	Date <i>23 Aug 22</i>

