



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION
 Toxicology Section/Breath Alcohol Program
 SFN 59281 (1-2010)

Serial Number: <u>80-004936</u>	Instrument Location: <u>Mint AFB</u>
Reason for Install/Repair:	
<input type="checkbox"/> Install After Lab Repair/Inspection <input checked="" type="checkbox"/> Other (Specify) <u>New Install</u>	

Check When Done:

- A. Surge Protector Installed/Properly Grounded.
- B. Telephone Line Connected to Intoxilyzer®.
- C. Breath Tube Heated.
- D. Review/Enter Preliminary Data Entry (Level 2, Function E).
- E. Review/Enter Gas Setup (Level 1, Function S).
- F. Print Test (Level 1, Function P). (Attach Test Record).
- G. ACA Test (Level 1, Function C). (Attach Test Record).
- H. RFI Test (Use CMS Mode-First Room Air). (Attach Test Record).
- I. Repair and/or Maintenance Performed: None Required

W. [Signature]
 Field Inspector's Signature

17 FEB 10
 Date

Deb Kasher
 Reviewed By

10 May 2010
 Date

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = MAFB 8164.13.00 06/09
02/17/2010 08:08

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 659358
Standard Cyl #? 49
Standard Expiration? 09/03/2011
Oper No? 020809

Flow Cal. Date: 01/18/2010
Slope 732
Intercept -767769

IR Calibration Date: 02/09/2010
 3um 9um

0th Coef(*100): -33294 -38146
1st Coef(*100): 263151 133532
2nd Coef(*100): -1136 209
H2O adj(mg/l*10k): 633 600

***** Printer Test End *****

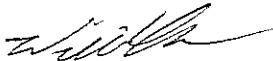

Operator Signature
WILLIAM CLARK

Remarks:


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = MAFB 8164.13.00 06/09
02/17/2010 08:35

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:36
02 Std. Gas	0.079	08:36
03 Room Air	0.000	08:37
04 Std. Gas	0.078	08:37
05 Room Air	0.000	08:38
06 Std. Gas	0.079	08:38
07 Room Air	0.000	08:39

Lot No = 659358
Cyl No = 49
Exp Date = 09/03/2011
County = 51 Oper No. = 020809

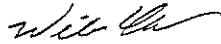

Operator Signature
WILLIAM CLARK

Remarks:

Form 106-18000

80-004936
P. 37/40 KM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = MAFB 8164.13.00 06/09
02/17/2010 08:30

Test	AC	Time
01 Diagnostic	OK	08:32
02 Room Air	RFI*	08:32
03 Room Air	RFI*	08:32

*Invalid Test
Inhibited - RFI

Sub Name = CLARK, WILLIAM NA
Sub DOB = 08/13/1983
Sub Sex = Male Weight = 210
Test = OTH Cit = NA
Dr. Lic. = ND/CLA831711
Lot No = 659358
Cyl No = 49
Expiration Date = 09/03/2011
County = 51 Oper No. = 020809

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

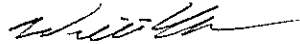

Operator Signature
WILLIAM CLARK

Remarks:

Form 106-I8000

80-004936
P. 40/40 JH