



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT

OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION
Toxicology Section/Breath Alcohol Program
SFN 59281 (1-2010)

Serial Number: ^{DK} 80-004943	Instrument Location: STAN ^{DK}
Reason for Install/Repair: <input checked="" type="checkbox"/> Install After Lab Repair/Inspection <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- A. Surge Protector Installed/Properly Grounded.
- B. Telephone Line Connected to Intoxilyzer® 8000.
- C. Breath Tube Heated.
- D. Review/Enter Preliminary Data Entry (Level 2, Function E).
- E. Review/Enter Gas Setup (Level 1, Function S).
- F. Print Test (Level 1, Function P). (Attach Test Record.)
- G. ACA Test (Level 1, Function C). (Attach Test Record.)
- H. RFI Test (Use CMS Mode-First Room Air). (Attach Test Record.)
- I. Repair and/or Maintenance Performed: _____

<u><i>[Signature]</i></u> Field Inspector Signature	<u>03/19/11</u> Date
<u>Deb Kashner</u> Reviewed By	<u>22 Mar 11</u> Date

Does not scan!

80-004943
P. 65 of 68
FM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004943
Location = STAN 8164.13.00 06/09
03/19/2011 19:07

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPNOPQRSTUVWXYZ!@#%&^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPNOPQRSTUVWXYZ!@#%&^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? No
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 709352
Standard Cyl #? 34
Standard Expiration? 12/23/2011
Oper No? 100019

Flow Cal. Date: 01/18/2010
Slope 693
Intercept -665387

IR Calibration Date: 01/18/2010
 3um 9um

0th Coef(*100):	-20351	-24446
1st Coef(*100):	248542	131200
2nd Coef(*100):	2180	1190
H2O adj(mg/l*10k):	594	450

***** Printer Test End *****

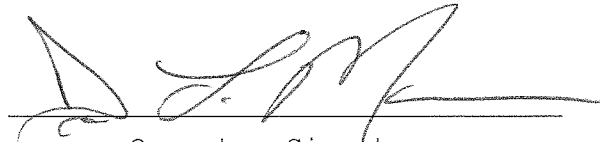

Operator Signature
DUANE MARMON

Remarks:


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004943
Location = STAN 8164.13.00 06/09
03/19/2011 19:00

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	19:01
02 Std. Gas	0.078	19:01
03 Room Air	0.000	19:02
04 Std. Gas	0.079	19:02
05 Room Air	0.000	19:03
06 Std. Gas	0.079	19:03
07 Room Air	0.000	19:03

Lot No = 709352
Cyl No = 34
Exp Date = 12/23/2011
County = 31 Oper No. = 100019

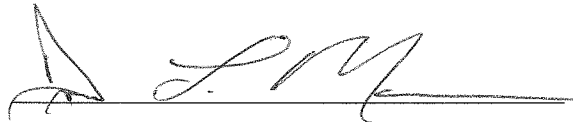

Operator Signature
DUANE MARMON

Remarks:

Form 106-I8000

80-004943
P. 67 of 68
JM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004943
Location = STAN 8164.13.00 06/09
03/19/2011 19:11

Test	AC	Time
01 Diagnostic	OK	19:14
02 Room Air	RFI*	19:14
03 Room Air	0.000	19:14

*Invalid Test
Inhibited - RFI

Sub Name = SUBJECT, TEST P
Sub DOB = 01/02/1989
Sub Sex = Female Weight = 350
Test = OTH Cit = 1234
Dr. Lic. = ND/SUB891234
Lot No = 709352
Cyl No = 34
Expiration Date = 12/23/2011
County = 31 Oper No. = 100019

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

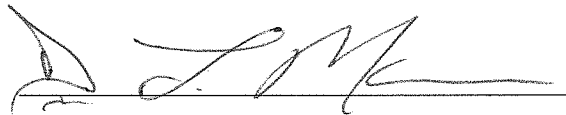

Operator Signature
DUANE MARMON

Remarks:

Form 106-I8000

80-004943
P. 68 of 68
SM