

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-003059

Location: TOXL

- Check When Done
- A. Is the breath tube heated? ✓
 - B. Is the diagnostic check complete? ✓
 - C. Is the time, date, and year correct (re-set if necessary)? ✓
 - D. Print test. (Attach test record.) ✓
 - E. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR3377 Lot No. 13270 Rep AC 0.020 ✓
 - F. High AC. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5189 Lot No. 13300 Rep AC 0.300 ✓
 - G. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.)
Sim SN DR5135 Lot No. 10542 Rep AC 0.100AC+0.05% ACETONE ✓
 - H. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
 - I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 16512080A2 Cyl No. 034 Exp. Date 08/01/2014
Test 1 0.080 Test 4 0.081 Test 7 0.081
Test 2 0.081 Test 5 0.082 Test 8 0.081
Test 3 0.081 Test 6 0.081 Test 9 0.080
Average 0.081

J. Tank pressure, Level 3 Func D Sub G P: 68 psi Reg 75 psi

K. Remarks/Maintenance record: REPLACED SIMULATOR RETURN O-RING,

ANNUAL INSPECTION

15 APR 2014
Date

15 Apr 2014
Date

Charles E. Eden
Field Inspector's Signature

Deb Kashner
Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 14:49

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&*()_+?

Current Instrument Setup

Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 16512080A2
Standard Cyl #? 34
Standard Expiration? 08/01/2014
Oper No? 666666

Flow Cal. Date: 04/07/2008
Slope 713
Intercept -692349

IR Calibration Date: 07/24/2013
 3um 9um

0th Coef(*100): -16999 -10479
1st Coef(*100): 275659 137940
2nd Coef(*100): 2781 1186
H2O adj(mg/l*10k): 674 466

***** Printer Test End *****



Operator Signature
CHARLES EDER

Remarks:

PRINT TEST

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 14:51

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:52
02 Std. Sol.	0.021	14:53
03 Room Air	0.000	14:53
04 Std. Sol.	0.020	14:54
05 Room Air	0.000	14:54
06 Std. Sol.	0.020	14:55
07 Room Air	0.000	14:56

08 Sim Temp = 34.0°C

Simul Ser No = DR3377
Std Sol No = 13270
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

Low AC
0.020 AC

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 14:57

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:58
02 Std. Sol.	0.299	14:58
03 Room Air	0.000	14:59
04 Std. Sol.	0.301	15:00
05 Room Air	0.000	15:00
06 Std. Sol.	0.302	15:01
07 Room Air	0.000	15:02


08 Sim Temp = 34.0°C

Simul Ser No = DR5189

Std Sol No = 13300

County = 08

Oper No. = 666666

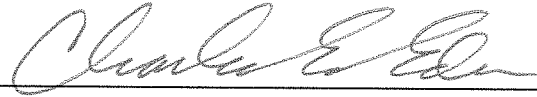

Operator Signature
CHARLES EDER

Remarks:

HIGH AC
0.300 AC

Form 106-I8000


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 15:03

Test	AC	Time
01 Room Air	0.000	15:04
02 *Subject Test	INT*	15:04
03 Room Air	0.000	15:05

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS821456
Lot No = 16512080A2
Cyl No = 34
Expiration Date = 08/01/2014
County = 08 Oper No. = 666666

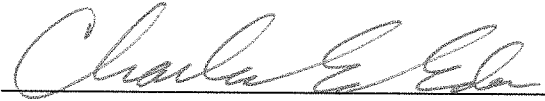

Operator Signature
CHARLES EDER

Remarks: *INTERFERENT CHECK*
0.100 AC + 0.05% ACETONE

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 15:07

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:08
02 Std. Gas	0.080	15:08
03 Room Air	0.000	15:08
04 Std. Gas	0.081	15:09
05 Room Air	0.000	15:09
06 Std. Gas	0.081	15:10
07 Room Air	0.000	15:10

Lot No = 16512080A2
Cyl No = 34
Exp Date = 08/01/2014
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 15:10

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:11
02 Std. Gas	0.081	15:11
03 Room Air	0.000	15:12
04 Std. Gas	0.082	15:12
05 Room Air	0.000	15:13
06 Std. Gas	0.081	15:13
07 Room Air	0.000	15:13

Lot No = 16512080A2

Cyl No = 34

Exp Date = 08/01/2014

County = 08

Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
04/15/2014 15:14

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:14
02 Std. Gas	0.081	15:15
03 Room Air	0.000	15:15
04 Std. Gas	0.081	15:15
05 Room Air	0.000	15:16
06 Std. Gas	0.080	15:16
07 Room Air	0.000	15:17

Lot No = 16512080A2
Cyl No = 34
Exp Date = 08/01/2014
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Form 106-I8000