



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION
 Toxicology Section/Breath Alcohol Program
 SFN 59281 (03-2014)

Serial Number <i>80-006504</i>	Instrument Location <i>Bott</i>
Reason for Install/Repair <input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Properly Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. RFI Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if Any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer®8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>Tyler Wolfe</i>	Date <i>05-12-16</i> ✓
Reviewed By (Crime Laboratory Use Only) <i>Deb Shanaver</i>	Date <i>20 May 16</i>

This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Certified By: *Charles E. Edin* Certified Date: *10 Nov. 16*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006504
Location = BOTT 8164.13.00 06/09
05/12/2016 10:09

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:09
02 Std. Gas	0.079	10:10
03 Room Air	0.000	10:10
04 Std. Gas	0.079	10:11
05 Room Air	0.000	10:11
06 Std. Gas	0.079	10:11
07 Room Air	0.000	10:12

Lot No = 01316080A2
Cyl No = 29
Exp Date = 02/05/2018
County = 05 Oper No. = 060901

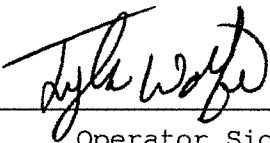

Operator Signature
TYLER WOLFE

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006504
Location = BOTT 8164.13.00 06/09
05/12/2016 10:37

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:38
02 Std. Gas	0.079	10:38
03 Room Air	RFI*	10:38
04 Room Air	0.000	10:39

*Invalid Test
Inhibited - RFI

Lot No = 01316080A2
Cyl No = 29
Exp Date = 02/05/2018
County = 05 Oper No. = 060901

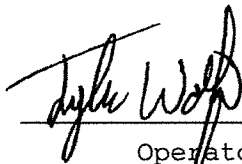

Operator Signature
TYLER WOLFE

Remarks: