



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION
 Toxicology Section/Breath Alcohol Program
 SFN 59281 (5-2011)

Serial Number 80-004193	Instrument Location FYPD - Breath Alcohol Testing Vehicle
Reason for Install/Repair	
<input type="checkbox"/> Install After Receiving From Crime Laboratory <input checked="" type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Properly Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P). ✓
 - B. ACA Test (Level 1, Function C). ✓
 - C. RFI Test (Use CMS Mode; Key Radio During First Room Air). ✓
- 7. Repair and/or Maintenance Performed (if Any): N/A
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer®8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature Sterling Reel	Date 05-16-13
Reviewed By (Crime Laboratory Use Only) Deb Kashur	Date 20 May 13

This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Certified By: **Charles E. Eden** Certified Date: **08 Nov. 16**

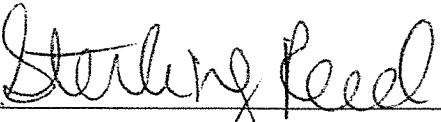
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = FYPD 8164.13.00 06/09
05/16/2013 11:41

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:41
02 Std. Gas	0.078	11:42
03 Room Air	0.000	11:42
04 Std. Gas	0.078	11:42
05 Room Air	0.000	11:43
06 Std. Gas	0.078	11:43
07 Room Air	0.000	11:44

Lot No = 16512080A2
Cyl No = 25
Exp Date = 08/01/2014
County = 61 Oper No. = 020717


Operator Signature
STERLING R. REED

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = FYPD 8164.13.00 06/09
05/16/2013 11:44

Test	AC	Time
01 Room Air	RFI*	11:46
02 Room Air	0.000	11:47

*Invalid Test
Inhibited - RFI

Sub Name = TEST, TEST NA
Sub DOB = 12/12/1975
Sub Sex = Male Weight = 200
Test = MIC Cit = 0001
Dr. Lic. = ND/TES750001
Lot No = 16512080A2
Cyl No = 25
Expiration Date = 08/01/2014
County = 61 Oper No. = 020717

Sterling Reed

Operator Signature
STERLING R. REED

Remarks:

RFI TEST

Form 106-I8000

RFI test done in ABA mode. DK 20 May 13