

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-005937

Location: STEE (KIDDER COUNTY S.O.)

- Check When Done
- A. Is the breath tube heated? ✓
 - B. Is the diagnostic check complete? ✓
 - C. Is the time, date, and year correct (re-set if necessary)? ✓
 - D. Print test. (Attach test record.) ✓
 - E. Low AC. Use ≤ 0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5191 Lot No. 15160 Rep AC 0.020 ✓
 - F. High AC. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR3847 Lot No. 16060 Rep AC 0.300 ✓
 - G. Interferent check. Use 0.05% acetone in ethanol solution in ABA mode.
(Attach test record.)
Sim SN DR5190 Lot No. IC54 Solution 0.10AC + 0.05% ACETONE ✓
 - H. RFI check. Run CMS mode. Key radio during CMS test.
(Attach test record.) ✓
 - I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.) 0.080 AC

Lot No. 19615080A5 Cyl No. 27 Exp. Date 09/05/2017

Test 1 <u>0.079</u>	Test 4 <u>0.078</u>	Test 7 <u>0.079</u>
Test 2 <u>0.080</u>	Test 5 <u>0.079</u>	Test 8 <u>0.079</u>
Test 3 <u>0.080</u>	Test 6 <u>0.080</u>	Test 9 <u>0.080</u>

Average 0.079

J. Tank pressure, Level 3 Func D Sub G P: 808 psi Reg 800 psi

K. Remarks/Maintenance record: ANNUAL INSPECTION, REPLACED BREATH HOSE O-RING.

5/13/2016
Date

Charles E. [Signature]
Field Inspector's Signature

He May 16 ✓
Date

Deb Shanaver
Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:17

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:18
02 Std. Sol.	0.018	15:19
03 Room Air	0.000	15:19
04 Std. Sol.	0.018	15:20
05 Room Air	0.000	15:21
06 Std. Sol.	0.019	15:21
07 Room Air	0.000	15:22


08 Sim Temp = 34.0°C

Simul Ser No = DR5191

Std Sol No = 15160

County = 22

Oper No. = 666666

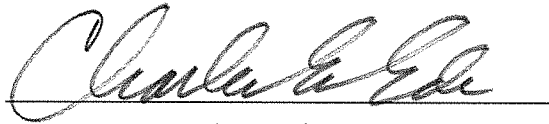

Operator Signature
CHARLES EDER

Remarks:

LOW AC
0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:22

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:23
02 Std. Sol.	0.306	15:24
03 Room Air	0.000	15:24
04 Std. Sol.	0.305	15:25
05 Room Air	0.000	15:26
06 Std. Sol.	0.305	15:26
07 Room Air	0.000	15:27

08 Sim Temp = 34.0°C

Simul Ser No = DR3847

Std Sol No = 16060

County = 22

Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

HIGH AC
0.300 AC

Form 106-I8000


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 14:59

Test	AC	Time
01 Room Air	0.000	15:00
02 *Subject Test	INT*	15:02
03 Room Air	0.000	15:02

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS821456
Lot No = 19615080A5
Cyl No = 27
Expiration Date = 09/05/2017
County = 22 Oper No. = 666666

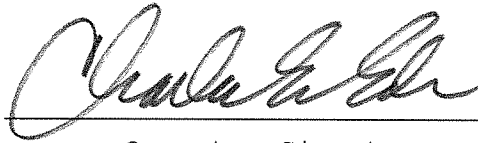

Operator Signature
CHARLES EDER

Remarks:

Interferent Check
0.10AC + 0.05% ACETONE

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:03

Test	AC	Time
01 Diagnostic	OK	15:04
02 Room Air	0.000	15:04
03 *Subject Test	RFI*	15:05
04 Room Air	0.000	15:05

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = RFI CHECK
Dr. Lic. = ND/DIS821456
Lot No = 19615080A5
Cyl No = 27
Expiration Date = 09/05/2017
County = 22 Oper No. = 666666

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

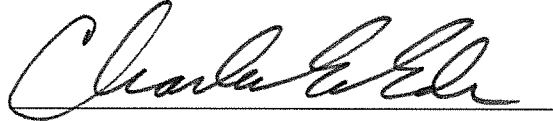

Operator Signature
CHARLES EDER

Remarks:

RFI CHECK


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:07

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:07
02 Std. Gas	0.079	15:08
03 Room Air	0.000	15:08
04 Std. Gas	0.080	15:09
05 Room Air	0.000	15:09
06 Std. Gas	0.080	15:10
07 Room Air	0.000	15:10

Lot No = 19615080A5
Cyl No = 27
Exp Date = 09/05/2017
County = 22 Oper No. = 666666


Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:28

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:29
02 Std. Gas	0.078	15:29
03 Room Air	0.000	15:29
04 Std. Gas	0.079	15:30
05 Room Air	0.000	15:30
06 Std. Gas	0.080	15:31
07 Room Air	0.000	15:31

Lot No = 19615080A5
Cyl No = 27
Exp Date = 09/05/2017
County = 22 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = STEE 8164.13.00 06/09
05/13/2016 15:31

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:32
02 Std. Gas	0.079	15:32
03 Room Air	0.000	15:33
04 Std. Gas	0.079	15:33
05 Room Air	0.000	15:34
06 Std. Gas	0.080	15:34
07 Room Air	0.000	15:34

Lot No = 19615080A5
Cyl No = 27
Exp Date = 09/05/2017
County = 22 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: *CALIBRATION CHECK*
0.080 AC