

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-004196

Location: Minot Air Force Base

- Check When Done
- A. Is the breath tube heated? ✓
 - B. Is the diagnostic check complete? ✓
 - C. Is the time, date, and year correct (re-set if necessary)? ✓
 - D. Print test. (Attach test record.) ✓
 - E. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5115 Lot No. 12111 Rep AC 0.020 ✓
 - F. High AC. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5143 Lot No. 13080 Rep AC 0.300 ✓
 - G. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.)
Sim SN DR3375 Lot No. 105#2 Rep AC 0.10 AC + 0.05% acetone ✓
 - H. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
 - I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 16512080A2 Cyl No. 38 Exp. Date 08/01/2014

Test 1 <u>0.078</u>	Test 4 <u>0.078</u>	Test 7 <u>0.079</u>
Test 2 <u>0.079</u>	Test 5 <u>0.079</u>	Test 8 <u>0.079</u>
Test 3 <u>0.079</u>	Test 6 <u>0.079</u>	Test 9 <u>0.079</u>

Average 0.079

J. Tank pressure, Level 3 Func D Sub G P: 920 psi Reg 930 psi

K. Remarks/Maintenance record: Annual inspection

21 May 2014
Date

22 MAY 2014 ✓
Date CEE

Deb Kasler
Field Inspector's Signature

Charles E. Ed
Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:23

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-|=|
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-|=|
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 16512080A2
Standard Cyl #? 38
Standard Expiration? 08/01/2014
Oper No? 777777

Flow Cal. Date: 06/04/2009
Slope 727
Intercept -600779

IR Calibration Date: 04/25/2013
 3um 9um

0th Coef(*100): -31440 -19588
1st Coef(*100): 280397 135844
2nd Coef(*100): -1584 629
H2O adj(mg/l*10k): 540 419

***** Printer Test End *****

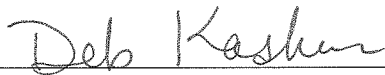

Operator Signature
DEB KASHUR

Remarks:

Print test

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:42

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:43
02 Std. Sol.	0.020	10:44
03 Room Air	0.000	10:44
04 Std. Sol.	0.020	10:45
05 Room Air	0.000	10:45
06 Std. Sol.	0.020	10:46
07 Room Air	0.000	10:47

08 Sim Temp = 34.0°C

Simul Ser No = DR5115

Std Sol No = 12111

County = 64

Oper No. = 777777

Deb Kasher

Operator Signature
DEB KASHUR

Remarks:

Low AC 0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:49

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:49
02 Std. Sol.	0.302	10:50
03 Room Air	0.000	10:50
04 Std. Sol.	0.302	10:51
05 Room Air	0.000	10:52
06 Std. Sol.	0.302	10:53
07 Room Air	0.000	10:53

08 Sim Temp = 34.0°C

Simul Ser No = DR5143

Std Sol No = 13080

County = 64

Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

High AC 0.300 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:28

Test	AC	Time
01 Room Air	0.000	10:28
02 *Subject Test	INT*	10:29
03 Room Air	0.000	10:30

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS921456
Lot No = 16512080A2
Cyl No = 38
Expiration Date = 08/01/2014
County = 64 Oper No. = 777777

Deb Kasher

Operator Signature
DEB KASHUR

Remarks:

Interferent check 0.10 Ae + 0.05% acetone

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:40

Test	AC	Time
01 Diagnostic	OK	10:41
02 Room Air	RFI*	10:41
03 Room Air	0.000	10:41

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS921456
Lot No = 16512080A2
Cyl No = 38
Expiration Date = 08/01/2014
County = 64 Oper No. = 777777

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

Debs Kashur

Operator Signature
DEB KASHUR

Remarks:

RFI check

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:24

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:24
02 Std. Gas	0.078	10:25
03 Room Air	0.000	10:25
04 Std. Gas	0.079	10:26
05 Room Air	0.000	10:26
06 Std. Gas	0.079	10:27
07 Room Air	0.000	10:27

Lot No = 16512080A2
Cyl No = 38
Exp Date = 08/01/2014
County = 64 Oper No. = 777777

Deb Kasher
Operator Signature
DEB KASHUR

Remarks:

Calibration check 0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:31

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:32
02 Std. Gas	0.078	10:32
03 Room Air	0.000	10:33
04 Std. Gas	0.079	10:33
05 Room Air	0.000	10:34
06 Std. Gas	0.079	10:34
07 Room Air	0.000	10:35

Lot No = 16512080A2

Cyl No = 38

Exp Date = 08/01/2014

County = 64

Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Calibration check 0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004196
Location = MNAF 8164.13.00 06/09
05/21/2014 10:36

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:36
02 Std. Gas	0.079	10:37
03 Room Air	0.000	10:37
04 Std. Gas	0.079	10:37
05 Room Air	0.000	10:38
06 Std. Gas	0.079	10:38
07 Room Air	0.000	10:39

Lot No = 16512080A2
Cyl No = 38
Exp Date = 08/01/2014
County = 64 Oper No. = 777777

Deb Kashur
Operator Signature
DEB KASHUR

Remarks:

Calibration check 0.080 AC

Form 106-I8000