

INTOXILYZER 8000 INSPECTION

Serial No.: 80-004205

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)  
Sim SN DL3455 Lot No.      Rep AC 0.021 ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)  
Sim SN DL5174 Lot No.      Rep AC 0.294 ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.  
(Attach test record.)  
Sim SN DL3452 Lot No.      Rep AC 0.10AC + Acetone ✓
- J. RFI check. Run CMS mode. Key radio on first room air.  
(Attach test record.) ✓

K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.  
(Attach test record.)

Lot No. <u>15410080A1</u>	Cyl No. <u>21</u>	Exp. Date <u>7-1-12</u>
Test 1 <u>0.080</u>	Test 4 <u>0.080</u>	Test 7 <u>0.080</u>
Test 2 <u>0.080</u>	Test 5 <u>0.079</u>	Test 8 <u>0.080</u>
Test 3 <u>0.080</u>	Test 6 <u>0.080</u>	Test 9 <u>0.080</u>
Average <u>0.080</u>		

L. Tank pressure, Level 3 Func D Sub G P: 58 psi Reg 60 psi

M. Remarks/Maintenance record: Annual Inspection

6-1-11  
Date

[Signature]  
Field Inspector's Signature

02 June 2011  
Date

Debi Kashan  
Reviewed by

80-004205  
P. 114 of 122  
FM

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-004205  
Location = TOXL                    8164.13.00 06/09  
06/01/2011                            13:25

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopQRSTUVWXYZ!@#\$\$%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopQRSTUVWXYZ!@#\$\$%^&\*()\_+?

Current Instrument Setup

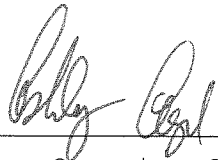
Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?        No  
Display Volume?            No  
Disable On Memfull?        Yes  
# of Print Copies?         1  
Select Std (D/W/I)?        Dry  
Standard Value?            0.080  
Standard Lot #?            15410080A1  
Standard Cyl #?            23<sup>rd</sup> 21  
Standard Expiration?       07/01/2012  
Oper No?                    555555

Flow Cal. Date:            06/04/2009  
Slope                      704  
Intercept                  -610062

IR Calibration Date:        06/03/2009

	3um	9um
0th Coef(*100):	-16944	-29673
1st Coef(*100):	265939	132317
2nd Coef(*100):	1696	900
H2O adj(mg/l*10k):	490	583

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
ASHLEY AYD

Remarks: *Print test*

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-004205  
Location = TOXL                    8164.13.00 06/09  
06/01/2011                            14:00

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:00
02 Std. Sol.	0.019	14:01
03 Room Air	0.000	14:02
04 Std. Sol.	0.019	14:02
05 Room Air	0.000	14:03
06 Std. Sol.	0.019	14:04
07 Room Air	0.000	14:04

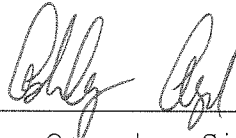
08 Sim Temp = 34.0°C

Simul Ser No = 0

Std Sol No = 0

County = 08

Oper No. = 555555



Operator Signature  
ASHLEY AYD

Remarks:

Low AC 0.021 AC

Form 106-I8000

80-004205  
P 116 of 122  
JFM

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004205  
Location = TOXL      8164.13.00 06/09  
06/01/2011      14:07

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:07
02 Std. Sol.	0.283	14:08
03 Room Air	0.000	14:09
04 Std. Sol.	0.285	14:09
05 Room Air	0.000	14:10
06 Std. Sol.	0.286	14:11
07 Room Air	0.000	14:11

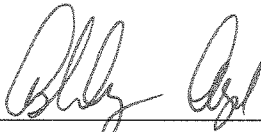
08 Sim Temp = 34.0°C

Simul Ser No = 0

Std Sol No = 0

County = 08

Oper No. = 555555

  
\_\_\_\_\_  
Operator Signature  
ASHLEY AYD

Remarks: *Linearity test 0.294AC*

Form 106-I8000

80-004205  
P. 117 of 122  
KEM

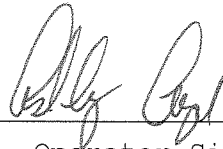
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004205  
Location = TOXL      8164.13.00 06/09  
06/01/2011      13:56

Test	AC	Time
01 Room Air	0.000	13:56
02 *Subject Test	INT*	13:57
03 Room Air	0.000	13:57

\*Invalid Test  
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 01/01/1982  
Sub Sex = Female      Weight = 150  
Test = DUI      Cit = NA  
Dr. Lic. = ND/DIS821456  
Lot No = 15410080A1  
Cyl No = 21  
Expiration Date = 07/01/2012  
County = 08      Oper No. = 555555



Operator Signature  
ASHLEY AYD

Remarks: *Interferent check 0.10AC + Acetone*

Form 106-I8000

80-004205  
P 11 P of 122  
JEM

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-004205  
Location = TOXL                 8164.13.00 06/09  
06/01/2011                        13:58

Test	AC	Time
01 Diagnostic	OK	13:58
02 Room Air	RFI*	13:58
03 Room Air	0.000	13:59

\*Invalid Test  
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 01/01/1982  
Sub Sex = Female                    Weight = 150  
Test = DUI                            Cit = NA  
Dr. Lic. = ND/DIS821456  
Lot No = 15410080A1  
Cyl No = 21  
Expiration Date = 07/01/2012  
County = 08                          Oper No. = 555555

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature  
ASHLEY AYD

Remarks: RFI

Form 106-I8000

80-004205  
P. 119 of 122  
JEM

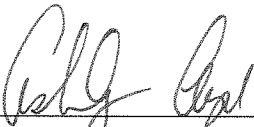
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-004205  
Location = TOXL                    8164.13.00 06/09  
06/01/2011                            13:27

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:27
02 Std. Gas	0.080	13:28
03 Room Air	0.000	13:28
04 Std. Gas	0.080	13:29
05 Room Air	0.000	13:29
06 Std. Gas	0.080	13:29
07 Room Air	0.000	13:30

Lot No = 15410080A1  
Cyl No = 21  
Exp Date = 07/01/2012  
County = 08                            Oper No. = 555555

  
\_\_\_\_\_  
Operator Signature  
ASHLEY AYD

Remarks: Calibration check 0.080AC

Form 106-I8000

80-004205  
01200102  
JPM

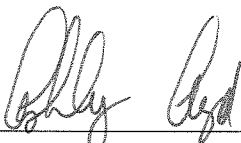
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004205  
Location = TOXL      8164.13.00 06/09  
06/01/2011      13:30

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:31
02 Std. Gas	0.080	13:31
03 Room Air	0.000	13:32
04 Std. Gas	0.079	13:32
05 Room Air	0.000	13:33
06 Std. Gas	0.080	13:33
07 Room Air	0.000	13:33

Lot No = 15410080A1  
Cyl No = 21  
Exp Date = 07/01/2012  
County = 08      Oper No. = 555555

  
\_\_\_\_\_  
Operator Signature  
ASHLEY AYD

Remarks: Calibration check 0.080 AC

Form 106-I8000

80-004205  
P. 12/07/22  
JAM



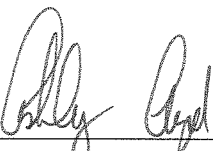
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-004205  
Location = TOXL                 8164.13.00 06/09  
06/01/2011                        13:34

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:35
02 Std. Gas	0.080	13:35
03 Room Air	0.000	13:35
04 Std. Gas	0.080	13:36
05 Room Air	0.000	13:36
06 Std. Gas	0.080	13:37
07 Room Air	0.000	13:37

Lot No = 15410080A1  
Cyl No = 21  
Exp Date = 07/01/2012  
County = 08                        Oper No. = 555555

  
\_\_\_\_\_  
Operator Signature  
ASHLEY AYD

Remarks: Calibration check 0.080 AC

80-004205  
R.1224122  
YPM