

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-003065

Location: GFSO CORR.

- Check When Done
- A. Is the breath tube heated? ✓
  - B. Is the diagnostic check complete? ✓
  - C. Is the time, date, and year correct (re-set if necessary)? ✓
  - D. Print test. (Attach test record.) ✓
  - E. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)  
Sim SN DR5115 Lot No. 12111 Rep AC 0.020 ✓
  - F. High AC. Use ≥ 0.25 AC in ACA mode. (Attach test record.)  
Sim SN DR5143 Lot No. 13080 Rep AC 0.300 ✓
  - G. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.  
(Attach test record.)  
Sim SN DR3375 Lot No. ICS#2 Rep AC 0.100 AC + 0.05% ACETONE ✓
  - H. RFI check. Run CMS mode. Key radio on first room air.  
(Attach test record.) ✓
  - I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.  
(Attach test record.)

Lot No. 33913080A4 Cyl No. 047 Exp. Date 01/01/2016

Test 1 <u>0.079</u>	Test 4 <u>0.078</u>	Test 7 <u>0.079</u>
Test 2 <u>0.079</u>	Test 5 <u>0.079</u>	Test 8 <u>0.078</u>
Test 3 <u>0.079</u>	Test 6 <u>0.078</u>	Test 9 <u>0.079</u>

Average 0.079

J. Tank pressure, Level 3 Func D Sub G P: 716 psi Reg 725 psi

K. Remarks/Maintenance record: REPLACED SIMULATOR RETURN O-RING,  
ANNUAL INSPECTION

03 JUN 14  
Date  
04 June 14 ✓  
Date

Charles E. Eden  
Field Inspector's Signature  
Deb Kashan  
Reviewed by

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      10:31

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxy1234567890-=  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%^&\*()\_+?

abcdefghijklmnopqrstuvwxy1234567890-=  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%^&\*()\_+?

Current Instrument Setup

Data Entry Mode:      Enabled  
Start Test Sequence:      DABACABA  
Display Prelim Rslt?      Yes  
Display Third Digit?      Yes  
Inhib Printer(Y/N)?      No  
Display Volume?      No  
Disable On Memfull?      Yes  
# of Print Copies?      1  
Select Std (D/W/I)?      Dry  
Standard Value?      0.080  
Standard Lot #?      33913080A4  
Standard Cyl #?      47  
Standard Expiration?      01/01/2016  
Oper No?      666666

Flow Cal. Date:      04/08/2008  
Slope      694  
Intercept      -539075

IR Calibration Date:      06/05/2008

	3um	9um
0th Coef(*100):	-9383	-9235
1st Coef(*100):	288128	138520
2nd Coef(*100):	957	781
H2O adj(mg/l*10k):	391	239

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
CHARLES EDER

Remarks:

PRINT  
TEST

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      10:38

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:38
02 Std. Sol.	0.022	10:39
03 Room Air	0.000	10:40
04 Std. Sol.	0.021	10:40
05 Room Air	0.000	10:41
06 Std. Sol.	0.022	10:42
07 Room Air	0.000	10:42

08 Sim Temp = 34.0°C

Simul Ser No = DR5115

Std Sol No = 12111

County = 18

Oper No. = 666666



Operator Signature  
CHARLES EDER

Remarks:

*Low AC*

*0.020 AC*

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      10:43

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:44
02 Std. Sol.	0.304	10:45
03 Room Air	0.000	10:45
04 Std. Sol.	0.305	10:46
05 Room Air	0.000	10:47
06 Std. Sol.	0.303	10:47
07 Room Air	0.000	10:48


08 Sim Temp = 34.0°C

Simul Ser No = DR5143

Std Sol No = 13080

County = 18

Oper No. = 666666



Operator Signature  
CHARLES EDER

Remarks:

HIGH AC  
0.300 AC

Form 106-I8000



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-003065  
Location = GFSO                    8164.13.00 06/09  
06/03/2014                            10:52

Test	AC	Time
01 Diagnostic	OK	10:53
02 Room Air	RFI*	10:53
03 Room Air	0.000	10:54

\*Invalid Test  
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 01/01/1982  
Sub Sex = Female                    Weight = 150  
Test = DUI                            Cit = RFI CHECK  
Dr. Lic. = ND/DIS821456  
Lot No = 33913080A4  
Cyl No = 47  
Expiration Date = 01/01/2016  
County = 18                            Oper No. = 666666

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

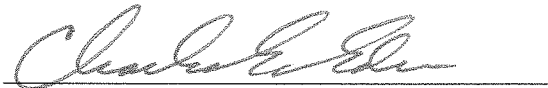

Operator Signature  
CHARLES EDER

Remarks:

RFI CHECK


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      10:54

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:55
02 Std. Gas	0.079	10:55
03 Room Air	0.000	10:56
04 Std. Gas	0.079	10:56
05 Room Air	0.000	10:57
06 Std. Gas	0.079	10:57
07 Room Air	0.000	10:58

Lot No = 33913080A4  
Cyl No = 47  
Exp Date = 01/01/2016  
County = 18      Oper No. = 666666

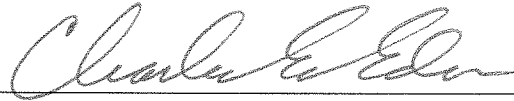

Operator Signature  
CHARLES EDER

Remarks: CALIBRATION  
CHECK  
0.080AC

Form 106-I8000


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      10:58

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:58
02 Std. Gas	0.078	10:59
03 Room Air	0.000	10:59
04 Std. Gas	0.079	11:00
05 Room Air	0.000	11:00
06 Std. Gas	0.078	11:00
07 Room Air	0.000	11:01

Lot No = 33913080A4  
Cyl No = 47  
Exp Date = 01/01/2016  
County = 18      Oper No. = 666666

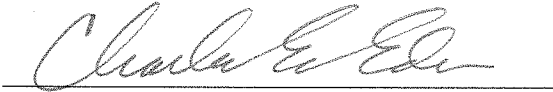

Operator Signature  
CHARLES EDER

Remarks: CALIBRATION  
CHECK  
0.080 AC

Form 106-I8000



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003065  
Location = GFSO      8164.13.00 06/09  
06/03/2014      11:01

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:02
02 Std. Gas	0.079	11:02
03 Room Air	0.000	11:03
04 Std. Gas	0.078	11:03
05 Room Air	0.000	11:03
06 Std. Gas	0.079	11:04
07 Room Air	0.000	11:04

Lot No = 33913080A4  
Cyl No = 47  
Exp Date = 01/01/2016  
County = 18      Oper No. = 666666



Operator Signature  
CHARLES EDER

Remarks:

CALIBRATION  
CHECK  
0.080 AC

Form 106-I8000