



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION
Toxicology Section/Breath Alcohol Program
SFN 59281 (5-2011)

Serial Number <i>80-004936</i>	Instrument Location <i>Burleigh County Sheriff's Dept</i>
Reason for Install/Repair <input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Properly Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. RFI Test (Use CMS Mode; Key Radio During First Room Air).
- 7. Repair and/or Maintenance Performed (if Any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer®8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>M. Gall</i>	Date <i>6-19-13</i>
Reviewed By (Crime Laboratory Use Only) <i>Deb Kashner</i>	Date <i>21 June 13</i>

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = BCSO 8164.13.00 06/09
06/19/2013 13:25

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 16512080A2
Standard Cyl #? 1
Standard Expiration? 08/01/2014
Oper No? 029802

Flow Cal. Date: 01/18/2010
Slope 732
Intercept -767769

IR Calibration Date: 04/26/2013
 3um 9um

	3um	9um
0th Coef(*100):	-20526	-23537
1st Coef(*100):	262392	130145
2nd Coef(*100):	2006	1425
H2O adj(mg/l*10k):	482	626

***** Printer Test End *****

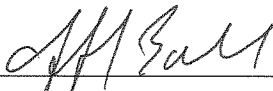

Operator Signature
JEFF BALL

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = BCSO 8164.13.00 06/09
06/19/2013 15:44

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:44
02 Std. Gas	0.078	15:45
03 Room Air	0.000	15:45
04 Std. Gas	0.078	15:45
05 Room Air	0.000	15:46
06 Std. Gas	0.078	15:46
07 Room Air	0.000	15:47

Lot No = 16512080A2

Cyl No = 1

Exp Date = 08/01/2014

County = 08

Oper No. = 029802



Operator Signature
JEFF BALL

Remarks:

ACA Test

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004936
Location = BCSO 8164.13.00 06/09
06/19/2013 16:11

Test	AC	Time
01 Diagnostic	OK	16:14
02 Room Air	RFI*	16:14
03 Room Air	0.000	16:14

*Invalid Test
Inhibited - RFI

Sub Name = SPIRIT, THE DISCOVER
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = 0000
Dr. Lic. = ND/DIS821456
Lot No = 16512080A2
Cyl No = 1
Expiration Date = 08/01/2014
County = 08 Oper No. = 029802

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature
JEFF BALL

Remarks:

Form 106-I8000

RFI Test