



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT

OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION
Toxicology Section/Breath Alcohol Program
SFN 59281 (03-2014)

Serial Number <i>80-005950</i>	Instrument Location <i>FTOT</i>
Reason for Install/Repair <input type="checkbox"/> Install After Receiving From Crime Laboratory <input checked="" type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Properly Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. RFI Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if Any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer®8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>Jacquie Demore</i>	Date <i>7-26-14</i>
Reviewed By (Crime Laboratory Use Only) <i>Deb Kashur</i>	Date <i>15 Sept 14</i>

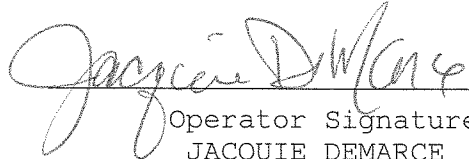
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005950
Location = FTOT 8164.13.00 06/09
07/26/2014 23:32

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	23:33
02 Std. Gas	0.080	23:33
03 Room Air	0.000	23:34
04 Std. Gas	0.081	23:34
05 Room Air	0.000	23:35
06 Std. Gas	0.081	23:35
07 Room Air	0.000	23:36

Lot No = 33913080A4
Cyl No = 46
Exp Date = 01/01/2016
County = 03 Oper No. = 100103


Operator Signature
JACQUIE DEMARCE

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

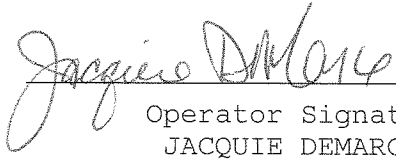
CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005950
Location = FTOT 8164.13.00 06/09
07/26/2014 23:29

DRY CAL CHECK

Test	AC	Time
01 Room Air	RFI*	23:30
02 Room Air	0.000	23:30

*Invalid Test
Inhibited - RFI

Lot No = 33913080A4
Cyl No = 46
Exp Date = 01/01/2016
County = 03 Oper No. = 100103


Operator Signature
JACQUIE DEMARCE

Remarks: