

*✓ emm*

**INTOXILYZER 5000 G and 8000 INSPECTION**

Serial No.: 80-004197

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓

---

- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.) ✓  
Sim SN DR5144 Lot No. --- Rep AC 0.021
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.) ✓  
Sim SN DR5187 Lot No. --- Rep AC 0.290
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode. (Attach test record.) ✓  
Sim SN DR5114 Lot No. --- Rep AC 0.100
- J. RFI check. Run CMS mode. Key radio on first room air. (Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode. (Attach test record.)

Lot No. <u>617774</u>	Cyl No. <u>38</u>	Exp. Date <u>28 May 2011</u>
Test 1 <u>0.681</u>	Test 4 <u>0.080</u>	Test 7 <u>0.080</u>
Test 2 <u>0.679</u>	Test 5 <u>0.080</u>	Test 8 <u>0.080</u>
Test 3 <u>0.080</u>	Test 6 <u>0.080</u>	Test 9 <u>0.080</u>
Average <u>0.080</u>		

L. Remarks/Maintenance record: Annual Inspection

<u>26 Aug 09</u> Date	<u>Deb Keshu</u> Field Inspector's Signature
<u>4 Sept 2009</u> Date	<u>Mary Pearson</u> Reviewed by

08/09

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      16:37

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

Current Instrument Setup  
Data Entry Mode:      Enabled  
Start Test Sequence:      DABACABA  
Display Prelim Rslt?      Yes  
Display Third Digit?      Yes  
Inhib Printer(Y/N)?      No  
Display Volume?      No  
Disable On Memfull?      Yes

---

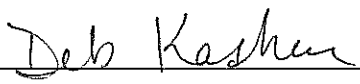
# of Print Copies?      1  
Select Std (D/W/I)?      Wet  
Standard Value?      0.000  
Standard Solution #?      000  
Sim Serial #?      DR0000  
Oper No?      777777

Flow Cal. Date:      06/04/2009  
Slope      697  
Intercept      -597078

IR Calibration Date:      06/03/2009

	3um	9um
0th Coef(*100):	-18382	-32622
1st Coef(*100):	278299	134602
2nd Coef(*100):	929	921
H2O adj(mg/l*10k):	352	627

\*\*\*\*\* Printer Test End \*\*\*\*\*

  
Operator Signature  
DEB KASHUR

Remarks:

*Print test*

Form 106-I8000

80-004197  
P 2 of 9 FR

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      17:07

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:07
02 Std. Sol.	0.016	17:08
03 Room Air	0.000	17:09
04 Std. Sol.	0.017	17:10
05 Room Air	0.000	17:10
06 Std. Sol.	0.017	17:11
07 Room Air	0.000	17:11

08 Sim Temp = 34.0°C

Simul Ser No = DR000

Std Sol No = 000

County = 08

Oper No. = 777777

Deb Kasher

Operator Signature  
DEB KASHUR

Remarks:

Low AC

Form 106-18000

80-004197  
P. 3 of 9 PK

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      16:58

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:59
02 Std. Sol.	0.289	17:01
03 Room Air	0.000	17:01
04 Std. Sol.	0.295	17:02
05 Room Air	0.000	17:02
06 Std. Sol.	0.294	17:03
07 Room Air	0.000	17:04

08 Sim Temp = 34.0°C

Simul Ser No = DR000

Std Sol No = 000

County = 08

Oper No. = 777777

*Deb Kasher*

Operator Signature  
DEB KASHUR

Remarks:

*Linearity test*

Form 106-18000

80-004197  
P. 4 of 9 JCH

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      17:13

Test	AC	Time
01 Room Air	0.000	17:14
02 *Subject Test	INT*	17:14
03 Room Air	0.000	17:15

\*Invalid Test  
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 02/01/1992  
Sub Sex = Female      Weight = 150  
Test = DUI      Cit = NA  
Dr. Lic. = ND/DIS921456  
Simul Ser No = DR000  
Std Sol No = 000  
County = 08      Oper No. = 777777

*Deb Kasher*

Operator Signature  
DEB KASHUR

Remarks:

*Interferent check*

Form 106-18000

80-004197  
P. 5 of 7 AK

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-004197  
Location = TOXL                    8164.12.00 06/09  
08/26/2009                            16:51

Test	AC	Time
01 Diagnostic	OK	16:52
02 Room Air	RFI*	16:52
03 Room Air	0.000	16:52

\*Invalid Test  
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 02/01/1992  
Sub Sex = Female                    Weight = 150  
Test = DUI                            Cit = NA  
Dr. Lic. = ND/DIS921456  
Simul Ser No = DR000  
Std Sol No = 000  
County = 08                            Oper No. = 777777

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

Deb Kasher  
Operator Signature  
DEB KASHUR

Remarks:

*RFI test*

Form 106-18000

*80-004197  
P. Conf? AH*

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      16:38

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:39
02 Std. Gas	0.081	16:39
03 Room Air	0.000	16:40
04 Std. Gas	0.079	16:40
05 Room Air	0.000	16:41
06 Std. Gas	0.080	16:41
07 Room Air	0.000	16:41

Lot No = 617774

Cyl No = 38

Exp Date = 05/28/2011

County = 08

Oper No. = 777777

*Deb Kashur*

Operator Signature  
DEB KASHUR

Remarks:

*Calibration check*

Form 106-I8000

*80-004197  
P. 2 of 8 PK*

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      16:42

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:43
02 Std. Gas	0.080	16:43
03 Room Air	0.000	16:44
04 Std. Gas	0.080	16:44
05 Room Air	0.000	16:44
06 Std. Gas	0.080	16:45
07 Room Air	0.000	16:45

---

Lot No = 617774  
Cyl No = 38  
Exp Date = 05/28/2011  
County = 08      Oper No. = 777777

Deb Kashur  
Operator Signature  
DEB KASHUR

Remarks:  
*Calibration check*

Form 106-I8000

80-004197  
P. 8/29/09



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004197  
Location = TOXL      8164.12.00 06/09  
08/26/2009      16:46

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:46
02 Std. Gas	0.080	16:47
03 Room Air	0.000	16:47
04 Std. Gas	0.080	16:48
05 Room Air	0.000	16:48
06 Std. Gas	0.080	16:48
07 Room Air	0.000	16:49

Lot No = 617774

Cyl No = 38

Exp Date = 05/28/2011

County = 08

Oper No. = 777777

*Deb Kashur*

Operator Signature  
DEB KASHUR

Remarks:

*Calibration check*

Form 106-I8000

80-004197  
P. 9 of 9 PM