

✓ e m l

INTOXILYZER 5000 G and 8000 INSPECTION

Serial No.: 80-004189

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓

- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.) ✓
Sim SN DR5190 Lot No. - Rep AC 0.021
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.) ✓
Sim SN DR5113 Lot No. - Rep AC 0.297
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode. (Attach test record.) ✓
Sim SN DR5135 Lot No. - Rep AC 0.10 AC + Acetone
- J. RFI check. Run CMS mode. Key radio on first room air. (Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode. (Attach test record.)

Lot No. 617774 Cyl No. 41 Exp. Date 5/28/2011

Test 1 <u>0.078</u>	Test 4 <u>0.079</u>	Test 7 <u>0.078</u>
Test 2 <u>0.079</u>	Test 5 <u>0.079</u>	Test 8 <u>0.078</u>
Test 3 <u>0.078</u>	Test 6 <u>0.078</u>	Test 9 <u>0.079</u>

Average 0.078

L. Remarks/Maintenance record: Annual Inspection

27 August 2009 Date Maryn Peann Field Inspector's Signature

04 Sept 2009 Date Deb Kashner Reviewed by

08/09

80-004189
P149 KM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 09:08

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Wet
Standard Value? 0.000
Standard Solution #? 0000
Sim Serial #? DR00000
Oper No? 107501

Flow Cal. Date: 06/03/2009
Slope 666
Intercept -550210

IR Calibration Date: 06/03/2009
 3um 9um

0th Coef(*100): -13136 -24026
1st Coef(*100): 256035 131167
2nd Coef(*100): 2113 1064
H2O adj(mg/l*10k): 465 385

***** Printer Test End *****

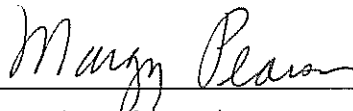

Operator Signature
MARGY PEARSON

Remarks:

Print Test

Form 106-I8000

80-004189
P. 2-19 PM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 08:49

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:49
02 Std. Sol.	0.017	08:50
03 Room Air	0.000	08:51
04 Std. Sol.	0.016	08:51
05 Room Air	0.000	08:52
06 Std. Sol.	0.017	08:52
07 Room Air	0.000	08:53

08 Sim Temp = 34.0°C

Simul Ser No = DR00000
Std Sol No = 0000
County = 08 Oper No. = 107501



Operator Signature
MARGY PEARSON

Remarks:

LOW AC

Form 106-I8000

80-004189
P. 3 of 9 FM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 09:00

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	09:01
02 Std. Sol.	0.285	09:03
03 Room Air	0.000	09:04
04 Std. Sol.	0.291	09:05
05 Room Air	0.000	09:05
06 Std. Sol.	0.290	09:06
07 Room Air	0.000	09:06

08 Sim Temp = 34.0°C

Simul Ser No = DR00000
Std Sol No = 0000
County = 08 Oper No. = 107501

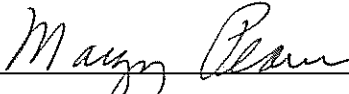

Operator Signature
MARGY PEARSON

Remarks:

Linearity

Form 106-I8000

80-004189
P. 4/9 PM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 08:46

Test	AC	Time
01 Room Air	0.000	08:47
02 *Subject Test	INT*	08:47
03 Room Air	0.000	08:48

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 11/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = DR00000
Std Sol No = 0000
County = 08 Oper No. = 107501



Operator Signature
MARGY PEARSON

Remarks:

Interfer

Form 106-18000

80-004189
P. 549 AM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 09:08

Test	AC	Time
01 Diagnostic	OK	09:09
02 Room Air	0.000	09:10
03 *Subject Test	RFI*	09:11
04 Room Air	0.000	09:11

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 11/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = DR00000
Std Sol No = 0000
County = 08 Oper No. = 107501

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

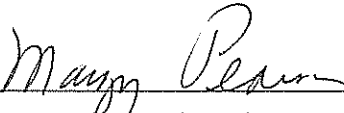

Operator Signature
MARGY PEARSON

Remarks:

RFI

Form 106-I8000

80-004189
8/27/09 JRM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/27/2009 07:57

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	07:58
02 Std. Gas	0.078	07:58
03 Room Air	0.000	07:59
04 Std. Gas	0.078	07:59
05 Room Air	0.000	07:59
06 Std. Gas	0.079	08:00
07 Room Air	0.000	08:00

Lot No = 617774
Cyl No = 41
Exp Date = 05/28/2011
County = 08 Oper No. = 107501

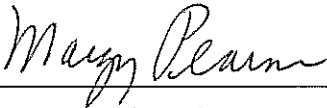
Operator Signature
MARGY PEARSON

Remarks:

AC A

Form 106-18000

80-004189
P. 2/9 KH

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/26/2009 17:14

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:14
02 Std. Gas	0.079	17:15
03 Room Air	0.000	17:15
04 Std. Gas	0.079	17:16
05 Room Air	0.000	17:16
06 Std. Gas	0.078	17:16
07 Room Air	0.000	17:17

Lot No = 617774
Cyl No = 41
Exp Date = 05/28/2011
County = 08 Oper No. = 107501



Operator Signature
MARGY PEARSON

Remarks:

AC A
Form 106-I8000

80-004189
P. 8-19 FM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004189
Location = TOXL 8164.12.00 06/09
08/26/2009 16:59

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:00
02 Std. Gas	0.078	17:00
03 Room Air	0.000	17:00
04 Std. Gas	0.079	17:01
05 Room Air	0.000	17:01
06 Std. Gas	0.078	17:02
07 Room Air	0.000	17:02

Lot No = 617774
Cyl No = 41
Exp Date = 05/28/2011
County = 08 Oper No. = 107501

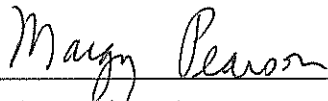

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

80-004189
P. 9/9 KM