

✓ cmn

INTOXILYZER 5000 G and 8000 INSPECTION

Serial No.: 8D-004198

Location: TOXC

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR3845 Lot No. — Rep AC 0.0217 ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5113 Lot No. — Rep AC 0.297 ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.) ✓
Sim SN DR 5135 Lot No. — Rep AC 0.10 AC + ACETONE ✓
- J. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 617774 Cyl No. 36 Exp. Date 5/28/2011
Test 1 0.080 Test 4 0.080 Test 7 0.079
Test 2 0.080 Test 5 0.080 Test 8 0.079
Test 3 0.081 Test 6 0.081 Test 9 0.079
Average 0.080 AC

L. Remarks/Maintenance record: Annual Inspection
31 August 2009 Mary Plann
Date Field Inspector's Signature
04 Sept 2009 Deb Kasher
Date Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 18:04

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	18:05
02 Std. Sol.	0.017	18:06
03 Room Air	0.000	18:06
04 Std. Sol.	0.017	18:07
05 Room Air	0.000	18:08
06 Std. Sol.	0.017	18:08
07 Room Air	0.000	18:09


08 Sim Temp = 34.0°C

Simul Ser No = DR0000

Std Sol No = 0000

County = 08

Oper No. = 107501

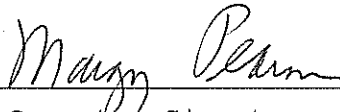

Operator Signature
MARGY PEARSON

Remarks:

LOW AC

Form 106-I8000

P. 3 of 9
FM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 17:06

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:07
02 Std. Sol.	0.292	17:08
03 Room Air	0.000	17:08
04 Std. Sol.	0.297	17:09
05 Room Air	0.000	17:09
06 Std. Sol.	0.296	17:10
07 Room Air	0.000	17:11


08 Sim Temp = 34.0°C

Simul Ser No = DR0000

Std Sol No = 0000

County = 08

Oper No. = 107501

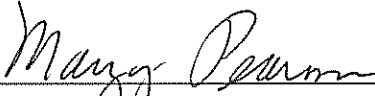

Operator Signature
MARGY PEARSON

Remarks:

Linearity

Form 106-18000

*P. 4 of 9
KM*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 18:14

Test	AC	Time
01 Room Air	0.000	18:15
02 *Subject Test	INT*	18:15
03 Room Air	0.000	18:15

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 11/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Lot No = 617774
Cyl No = 36
Expiration Date = 05/28/2011
County = 08 Oper No. = 107501

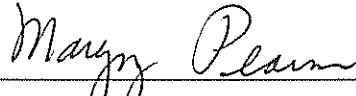

Operator Signature
MARGY PEARSON

Remarks:

AIC + ACETONE

Form 106-I8000

P. 5/9
JAN

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 09:34

Test	AC	Time
01 Diagnostic	OK	09:36
02 Room Air	RFI*	09:36
03 Room Air	0.000	09:37

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 11/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = DR0000
Std Sol No = 0000
County = 08 Oper No. = 107501

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature
MARGY PEARSON

Remarks:

RFI

Form 106-I8000

P. Cat 9
SM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 09:17

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	09:18
02 Std. Gas	0.080	09:18
03 Room Air	0.000	09:19
04 Std. Gas	0.080	09:19
05 Room Air	0.000	09:20
06 Std. Gas	0.081	09:20
07 Room Air	0.000	09:21

Lot No = 617774
Cyl No = 36
Exp Date = 05/28/2011
County = 08 Oper No. = 107501

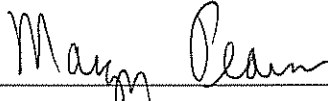

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

P. 7/9
m


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 09:13

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	09:14
02 Std. Gas	0.080	09:14
03 Room Air	0.000	09:15
04 Std. Gas	0.080	09:15
05 Room Air	0.000	09:16
06 Std. Gas	0.081	09:16
07 Room Air	0.000	09:17

Lot No = 617774
Cyl No = 36
Exp Date = 05/28/2011
County = 08 Oper No. = 107501

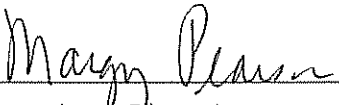

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

P. 8 of 9
FM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004198
Location = TOXL 8164.12.00 06/09
08/31/2009 18:16

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	18:16
02 Std. Gas	0.079	18:17
03 Room Air	0.000	18:17
04 Std. Gas	0.079	18:17
05 Room Air	0.000	18:18
06 Std. Gas	0.079	18:18
07 Room Air	0.000	18:19

Lot No = 617774
Cyl No = 36
Exp Date = 05/28/2011
County = 08 Oper No. = 107501



Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

P. 9 of 9
JPM