

**INTOXILYZER 5000 G and 8000 INSPECTION**

Serial No.: 80-003069

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)  
Sim SN DR5144 Lot No. — Rep AC 0.019 ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)  
Sim SN DR5187 Lot No. — Rep AC 0.285 ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.  
(Attach test record.)  
Sim SN DR5114 Lot No. — Rep AC 0.100 ✓
- J. RFI check. Run CMS mode. Key radio on first room air.  
(Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.  
(Attach test record.)

Lot No. 659358 Cyl No. 50 Exp. Date 03 Sept 2011

Test 1 0.078 Test 4 0.080 Test 7 0.080

Test 2 0.079 Test 5 0.080 Test 8 0.081

Test 3 0.080 Test 6 0.080 Test 9 0.081

Average 0.080

L. Remarks/Maintenance record: New software 8164.13 Annual inspection

12 Nov 09  
Date

Deb Kashur  
Field Inspector's Signature

1 Dec 09  
Date

Margy Pearson  
Reviewed by

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-003069  
Location = TOXL                    8164.13.00 06/09  
11/10/2009                            16:27

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxy1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

abcdefghijklmnopqrstuvwxy1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

Current Instrument Setup

Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?        No  
Display Volume?            No  
Disable On Memfull?        Yes  
# of Print Copies?          1  
Select Std (D/W/I)?        Wet  
Standard Value?            0.000  
Standard Solution #?        000  
Sim Serial #?              00000  
Oper No?                    777777

Flow Cal. Date:            04/08/2008  
Slope                        691  
Intercept                   -599264

IR Calibration Date:        06/04/2008  
                                  3um                            9um

	3um	9um
0th Coef(*100):	-14879	-13696
1st Coef(*100):	280678	136270
2nd Coef(*100):	177	615
H2O adj (mg/l*10k):	391	185

\*\*\*\*\* Printer Test End \*\*\*\*\*

Deb Keshur  
Operator Signature  
DEB KASHUR

Remarks:

*Print test*

Form 106-I8000

80-DC3069  
P. 2 of 9  
JKM

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/10/2009      16:19

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:19
02 Std. Sol.	0.018	16:21
03 Room Air	0.000	16:21
04 Std. Sol.	0.018	16:22
05 Room Air	0.000	16:23
06 Std. Sol.	0.018	16:23
07 Room Air	0.000	16:24

08 Sim Temp = 34.0°C

Simul Ser No = <sup>DR5144</sup>~~00000~~ DK

Std Sol No = ~~000~~ 0.019

County = 08

Oper No. = 777777

Deb Kashur

Operator Signature  
DEB KASHUR

Remarks:

low AC

Form 106-I8000

80-003069  
P. 3 of 9

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/12/2009      10:32

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:33
02 Std. Sol.	0.287	10:34
03 Room Air	0.000	10:34
04 Std. Sol.	0.291	10:35
05 Room Air	0.000	10:36
06 Std. Sol.	0.290	10:36
07 Room Air	0.000	10:37

08 Sim Temp = 34.0°C

Simul Ser No = <sup>DR5187</sup> 00000-01  
Std Sol No = <sup>0.285</sup> 000 0.285  
County = 08      Oper No. = 777777

Deb Kashur

Operator Signature  
DEB KASHUR

Remarks:

Linearity test

Form 106-I8000

80-003069  
P. 4 of 9  
JK

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/12/2009      10:46

Test	AC	Time
01 Room Air	0.000	10:46
02 *Subject Test	INT*	10:47
03 Room Air	0.000	10:47

\*Invalid Test  
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 02/01/1992  
Sub Sex = Female      Weight = 150  
Test = DUI      Cit = NA  
Dr. Lic. = ND/DIS921456  
Lot No = 659358  
Cyl No = 50  
Expiration Date = 09/03/2011  
County = 08      Oper No. = 777777

*Deb Kasher*

Operator Signature  
DEB KASHUR

Remarks:

*Interferent check*

Form 106-I8000

80-003069  
P. Sol 9/12/11

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-003069  
Location = TOXL                8164.13.00 06/09  
11/12/2009                        10:55

Test	AC	Time
01 Diagnostic	OK	10:56
02 Room Air	RFI*	10:56
03 Room Air	0.000	10:56

\*Invalid Test  
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 02/01/1992  
Sub Sex = Female                Weight = 150  
Test = DUI                        Cit = NA  
Dr. Lic. = ND/DIS921456  
Lot No = 659358  
Cyl No = 50  
Expiration Date = 09/03/2011  
County = 08                        Oper No. = 777777

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

Deb Kasher

Operator Signature  
DEB KASHUR

Remarks:

RFI test

Form 106-18000

80-603069  
P. 6 of 9  
JFM

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/12/2009      11:05

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:05
02 Std. Gas	0.078	11:06
03 Room Air	0.000	11:06
04 Std. Gas	0.079	11:06
05 Room Air	0.000	11:07
06 Std. Gas	0.080	11:07
07 Room Air	0.000	11:08

Lot No = 659358  
Cyl No = 35  
Exp Date = 09/03/2011  
County = 08      Oper No. = 777777

Deb Kasher  
Operator Signature  
DEB KASHUR

Remarks:

*Calibration check*

Form 106-I8000

80-003069  
p. 7 of 9

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/12/2009      11:08

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:09
02 Std. Gas	0.080	11:09
03 Room Air	0.000	11:10
04 Std. Gas	0.080	11:10
05 Room Air	0.000	11:11
06 Std. Gas	0.080	11:11
07 Room Air	0.000	11:11

Lot No = 659358  
Cyl No = 35  
Exp Date = 09/03/2011  
County = 08      Oper No. = 777777

*Deb Kashur*

Operator Signature  
DEB KASHUR

Remarks:

*Calibration check*

Form 106-I8000

80-003069  
P. 8 of 9  
JK

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003069  
Location = TOXL      8164.13.00 06/09  
11/12/2009      11:12

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:13
02 Std. Gas	0.080	11:13
03 Room Air	0.000	11:14
04 Std. Gas	0.081	11:14
05 Room Air	0.000	11:15
06 Std. Gas	0.081	11:15
07 Room Air	0.000	11:15

Lot No = 659358  
Cyl No = 35  
Exp Date = 09/03/2011  
County = 08      Oper No. = 777777

Deb Kasher  
Operator Signature  
DEB KASHUR

Remarks:

*Calibration check*

Form 106-I8000

80-003069  
P. 9 of 9 gm