



INTOXILYZER 5000 G and 8000 INSPECTION

Serial No.: 80-003059

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5190 Lot No. - Rep AC 0.021 ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5113 Lot No. - Rep AC 0.285 ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.)
Sim SN DR5135 Lot No. - Rep AC 0.10 AC ✓
- J. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 659358 Cyl No. 34 Exp. Date 9/03/2011
Test 1 0.081 Test 4 0.082 Test 7 0.082
Test 2 0.082 Test 5 0.082 Test 8 0.082
Test 3 0.082 Test 6 0.082 Test 9 0.082
Average 0.082

L. Remarks/Maintenance record: Software 8164.13 (Annual Inspec)
3^{mm} December 2009
Date 08 Dec 2009
Date 08 Dec 2009
Field Inspector's Signature Mary Plann
Reviewed by Deb Kasher

08/09

80-003059
P. 1 of 9
1/22

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:24

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:25
02 Std. Sol.	0.019	11:26
03 Room Air	0.000	11:26
04 Std. Sol.	0.020	11:27
05 Room Air	0.000	11:28
06 Std. Sol.	0.020	11:28
07 Room Air	0.000	11:29


08 Sim Temp = 34.0°C

Simul Ser No = 0000000000

Std Sol No = 0000000

County = 08

Oper No. = 107501

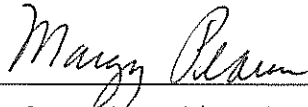

Operator Signature
MARGY PEARSON

Remarks:

Low 0.021 AC.

Form 106-18000

80-003059
P. 3 of 7 JAN

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:42

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:43
02 Std. Sol.	0.281	11:44
03 Room Air	0.000	11:44
04 Std. Sol.	0.299	11:45
05 Room Air	0.000	11:46
06 Std. Sol.	0.296	11:46
07 Room Air	0.000	11:47

08 Sim Temp = 34.0°C

Simul Ser No = 0000000000
Std Sol No = 0000000
County = 08 Oper No. = 107501

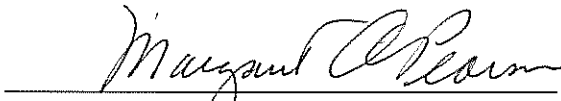

Operator Signature
MARGY PEARSON

Remarks:

Linearity 0.285% C.

Form 106-I8000

80-003059
P. 4 of 9 PM

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:49

Test	AC	Time
01 Diagnostic	OK	11:50
02 Room Air	0.000	11:51
03 *Subject Test	INT*	11:51
04 Room Air	0.000	11:52

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = 0000000000
Std Sol No = 0000000
County = 08 Oper No. = 107501

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

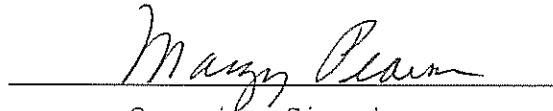

Operator Signature
MARGY PEARSON

Remarks:

0.10 AC Alcohol + 0.05% Acetone

Form 106-18000

80-003059
P. 5 of 9 JFH

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:39

Test	AC	Time
01 Diagnostic	OK	11:40
02 Room Air	RFI*	11:40
03 Room Air	0.000	11:41

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = 0000000000
Std Sol No = 0000000
County = 08 Oper No. = 107501

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

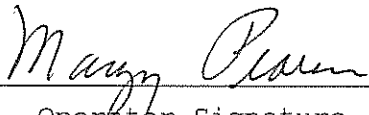

Operator Signature
MARGY PEARSON

Remarks:

RFI

Form 106-I8000

80-003059
P. 6 of 9 JEM


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:16

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:17
02 Std. Gas	0.081	11:17
03 Room Air	0.000	11:18
04 Std. Gas	0.082	11:18
05 Room Air	0.000	11:19
06 Std. Gas	0.082	11:19
07 Room Air	0.000	11:20

Lot No = 659358
Cyl No = 34
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

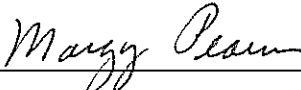

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-18000

80-003059
P. 2 of 9 Jan


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:20

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:20
02 Std. Gas	0.082	11:21
03 Room Air	0.000	11:21
04 Std. Gas	0.082	11:22
05 Room Air	0.000	11:22
06 Std. Gas	0.082	11:23
07 Room Air	0.000	11:23

Lot No = 659358
Cyl No = 34
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

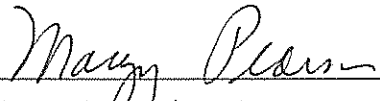

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

80-003059
P. 8-19-11


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003059
Location = TOXL 8164.13.00 06/09
12/03/2009 11:57

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:57
02 Std. Gas	0.082	11:58
03 Room Air	0.000	11:58
04 Std. Gas	0.082	11:58
05 Room Air	0.000	11:59
06 Std. Gas	0.082	11:59
07 Room Air	0.000	12:00

Lot No = 659358
Cyl No = 34
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

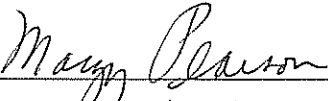

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

80-003059
P. 9 of 9 pages