

INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT

Serial No.: 80-003072 Instrument Location: GRAF

Reason for Install/Repair: _____ Install After Lab Repair/Inspection
x Other (Specify) New Install

Check When Done:

- x A. Surge Protector Installed/Properly Grounded.
- x B. Telephone Line Connected to Intoxilyzer®.
- x C. Breath Tube Heated.
- x D. Review/Enter Preliminary Data Entry (Level 2, Function E).
- x E. Review/Enter Gas Setup (Level 1, Function S).
- x F. Print Test (Level 1, Function P). (Attach Test Record.)
- x G. ACA Test (Level 1, Function C). (Attach Test Record.)
- x H. RFI Test (Use CMS Mode-First Room Air). (Attach Test Record.)
- I. Repair and/or Maintenance Performed: _____

12/4/09
Date


Field Inspector's Signature

15 Dec 2009
Date

Deb Kashur
Reviewed By

80-003072
P 12/4/09

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003072
Location = GRAF 8164.13.00 06/09
12/04/2009 13:58

***** Printer Test *****

abcdefghijklmnopqrstuvwxy1234567890- = |
ABCDEFGHIJKLMNPNQRSTUUVWXYZ!@#\$%^&*()_+?

abcdefghijklmnopqrstuvwxy1234567890- = |
ABCDEFGHIJKLMNPNQRSTUUVWXYZ!@#\$%^&*()_+?

Current Instrument Setup

Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 659358
Standard Cyl #? 16
Standard Expiration? 09/03/2011
Oper No? 049907

Flow Cal. Date: 06/05/2008
Slope 732
Intercept -720104

IR Calibration Date: 06/05/2008
 3um 9um

0th Coef(*100): -20044 -17692
1st Coef(*100): 264480 139417
2nd Coef(*100): 1329 776
H2O adj(mg/l*10k): 540 318

***** Printer Test End *****



Operator Signature
AARON HUMMEL

Remarks:

Print Test


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003072
Location = GRAF 8164.13.00 06/09
12/04/2009 13:59

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:00
02 Std. Gas	0.082	14:00
03 Room Air	0.000	14:00
04 Std. Gas	0.081	14:01
05 Room Air	0.000	14:01
06 Std. Gas	0.082	14:02
07 Room Air	0.000	14:02

Lot No = 659358
Cyl No = 16
Exp Date = 09/03/2011
County = 50 Oper No. = 049907

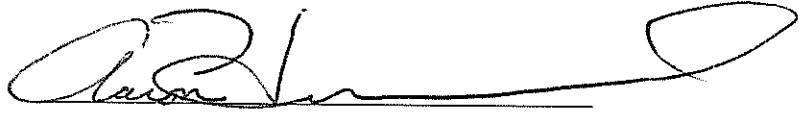

Operator Signature
AARON HUMMEL

Remarks:

Intox Install ACA

Form 106-I8000

*80-003072
p. 3 of 4 rkm*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003072
Location = GRAF 8164.13.00 06/09
12/04/2009 14:18

Test	AC	Time
01 Diagnostic	OK	14:21
02 Room Air	RFI*	14:21
03 Room Air	0.000	14:21

*Invalid Test
Inhibited - RFI

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

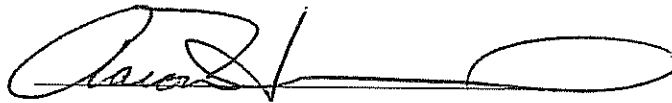

Operator Signature
AARON R. HUMMEL

Remarks:

RFI Test

Form 106-I8000

80-003072
0.4/4 pm