

INTOXILYZER 5000 G and 8000 INSPECTION

Serial No.: 80-004193

Location: TOX LAB

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5190 Lot No. — Rep AC 0.021 AC ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5113 Lot No. — Rep AC 0.285 AC ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.) ✓
Sim SN DR5135 Lot No. — Rep AC 0.10 AC ✓
- J. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 659358 Cyl No. 17 Exp. Date 09/03/2011
Test 1 0.082 Test 4 0.082 Test 7 0.082
Test 2 0.082 Test 5 0.083 Test 8 0.083
Test 3 0.082 Test 6 0.083 Test 9 0.082
Average 0.082

L. Remarks/Maintenance record: Software 864.13 (Annual Inspec)
4 December 2009 Margy Pearson
Date Field Inspector's Signature
08 Dec. 09 Deb Keshu
Date Reviewed by

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 10:52

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-==|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-==|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 659358
Standard Cyl #? 17
Standard Expiration? 09/03/2011
Oper No? 107501

Flow Cal. Date: 06/04/2009
Slope 692
Intercept -611138

IR Calibration Date: 06/03/2009
 3um 9um

0th Coef(*100): -12132 -15777
1st Coef(*100): 261160 129907
2nd Coef(*100): 1958 971
H2O adj(mg/l*10k): 413 335

***** Printer Test End *****

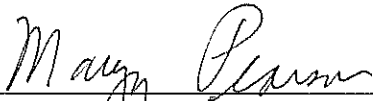

Operator Signature
MARGY PEARSON

Remarks:

Print

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 12:11

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:11
02 Std. Sol.	0.019	12:12
03 Room Air	0.000	12:13
04 Std. Sol.	0.019	12:14
05 Room Air	0.000	12:14
06 Std. Sol.	0.019	12:15
07 Room Air	0.000	12:15


08 Sim Temp = 34.0°C

Simul Ser No = 00000000

Std Sol No = 00000

County = 08

Oper No. = 107501

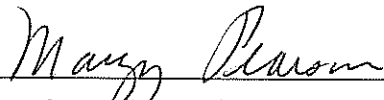

Operator Signature
MARGY PEARSON

Remarks:

Low AC 0.021 AC

Form 106-I8000

80-004193
P. 3 of 9 yzh

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 12:24

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:24
02 Std. Sol.	0.289	12:26
03 Room Air	0.000	12:26
04 Std. Sol.	0.295	12:27
05 Room Air	0.000	12:28
06 Std. Sol.	0.292	12:28
07 Room Air	0.000	12:29

08 Sim Temp = 34.0°C

Simul Ser No = 00000000

Std Sol No = 00000

County = 08

Oper No. = 107501



Operator Signature
MARGY PEARSON

Remarks:

Linearity - 0.285 AC

Form 106-18000

80-004193
P. 4/9/11


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 13:46

Test	AC	Time
01 Room Air	0.000	13:47
02 *Subject Test	INT*	13:47
03 Room Air	0.000	13:48

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = 00000000
Std Sol No = 00000
County = 08 Oper No. = 107501

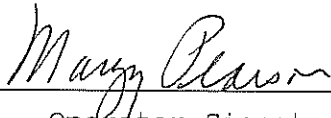

Operator Signature
MARGY PEARSON

Remarks:

Interferent Alc + Acetone

Form 106-I8000

*80-004193
P. 5 of 9, 10/11*

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 13:50

Test	AC	Time
01 Diagnostic	OK	13:51
02 Room Air	RFI*	13:51
03 Room Air	0.000	13:52

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS821456
Simul Ser No = 00000000
Std Sol No = 00000
County = 08 Oper No. = 107501

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

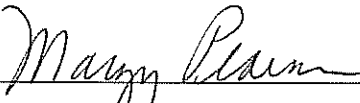

Operator Signature
MARGY PEARSON

Remarks:

RFI Test

Form 106-18000

80-004193
P. 6 of 9 1/24


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 14:06

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:07
02 Std. Gas	0.082	14:07
03 Room Air	0.000	14:08
04 Std. Gas	0.082	14:08
05 Room Air	0.000	14:09
06 Std. Gas	0.082	14:09
07 Room Air	0.000	14:09

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

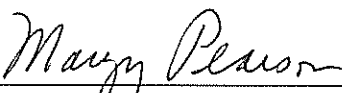

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-18000

*80-004193
p. 2 of 8 ym*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 14:10

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:10
02 Std. Gas	0.082	14:11
03 Room Air	0.000	14:11
04 Std. Gas	0.083	14:11
05 Room Air	0.000	14:12
06 Std. Gas	0.083	14:12
07 Room Air	0.000	14:13

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

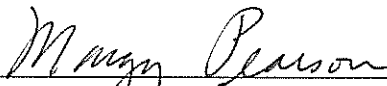

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

*80-004193
P. P. 8 15/11*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004193
Location = TOXL 8164.13.00 06/09
12/04/2009 14:13

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:14
02 Std. Gas	0.082	14:14
03 Room Air	0.000	14:15
04 Std. Gas	0.083	14:15
05 Room Air	0.000	14:16
06 Std. Gas	0.082	14:16
07 Room Air	0.000	14:17

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 107501

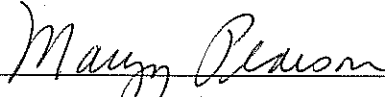

Operator Signature
MARGY PEARSON

Remarks:

ACA

Form 106-I8000

80-004193
P. 9 of 9 JAH