



INTOXILYZER 5000 G and 8000 INSPECTION

Serial No.: 80-004201

Location: TOXL

Check When Done

- A. Is the warm-up time less than 20 minutes? ✓
- B. Is a three-pronged grounded outlet used? ✓
- C. Is the breath tube heated? ✓
- D. Is the diagnostic check complete? ✓
- E. Is the time, date, and year correct (re-set if necessary)? ✓
- F. Print test. (Attach test record.) ✓
- G. Low AC. Use <0.03 AC in ACA mode. (Attach test record.)
Sim SN DR5144 Lot No. — Rep AC 0.019 ✓
- H. Linearity Test. Use ≥ 0.25 AC in ACA mode. (Attach test record.)
Sim SN DR5187 Lot No. — Rep AC 0.285 ✓
- I. Interferent check. Use 0.05% acetone plus 0.10 AC ethanol in ABA mode.
(Attach test record.)
Sim SN DR5114 Lot No. — Rep AC 0.160 ✓
- J. RFI check. Run CMS mode. Key radio on first room air.
(Attach test record.) ✓
- K. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.
(Attach test record.)

Lot No. 659358 Cyl No. 17 Exp. Date 03 Sept 2011

Test 1 0.080 Test 4 0.080 Test 7 0.079

Test 2 0.080 Test 5 0.080 Test 8 0.080

Test 3 0.079 Test 6 0.080 Test 9 0.080

Average 0.080

L. Remarks/Maintenance record: Software upgrade 8/16/13. Annual inspection.

22 Dec 2009
Date

Deb Kashur
Field Inspector's Signature

30 Dec 2009
Date

Maryn Pearson
Reviewed by

80-004201
P. 1 of 9
JAN

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 14:09

***** Printer Test *****

abcdefghijklmnopqrstuvxyz1234567890-|=|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%&^*()_+?

abcdefghijklmnopqrstuvxyz1234567890-|=|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%&^*()_+?

Current Instrument Setup

Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 659358
Standard Cyl #? 17
Standard Expiration? 09/03/2011
Oper No? 777777

Flow Cal. Date: 06/03/2009
Slope 696
Intercept -646527

IR Calibration Date: 12/23/2009

	3um	9um
0th Coef(*100):	-31049	-31538
1st Coef(*100):	280282	134511
2nd Coef(*100):	-770	412
H2O adj(mg/l*10k):	634	579

***** Printer Test End *****

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Print test

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 14:46

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:47
02 Std. Sol.	0.019	14:48
03 Room Air	0.000	14:49
04 Std. Sol.	0.019	14:49
05 Room Air	0.000	14:50
06 Std. Sol.	0.020	14:51
07 Room Air	0.000	14:51

08 Sim Temp = 34.0°C

Simul Ser No = 00000

Std Sol No = 000

County = 08

Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Low AC 0.019 AC

Form 106-I8000

80-004201
P. 3 of 9 Jan

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 14:58

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:59
02 Std. Sol.	0.286	14:59
03 Room Air	0.000	15:00
04 Std. Sol.	0.285	15:01
05 Room Air	0.000	15:01
06 Std. Sol.	0.286	15:02
07 Room Air	0.000	15:03

08 Sim Temp = 34.0°C

Simul Ser No = 00000

Std Sol No = 000

County = 08

Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Linearity test 0.285 AC

Form 106-18000

80-004201
P. 447/42

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 16:01

Test	AC	Time
01 Room Air	0.000	16:02
02 *Subject Test	INT*	16:02
03 Room Air	0.000	16:03

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS921456
Simul Ser No = 00000
Std Sol No = 000
County = 08 Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Interferent check

Form 106-I8000

80-004201
P. Sep 9 2009

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 16:23

Test	AC	Time
01 Diagnostic	OK	16:24
02 Room Air	RFI*	16:24
03 Room Air	0.000	16:24

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = NA
Dr. Lic. = ND/DIS921456
Lot No = 659358
Cyl No = 17
Expiration Date = 09/03/2011
County = 08 Oper No. = 777777

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature
DEB KASHUR

Remarks:

RFI test

Form 106-I8000

*80-004201
P. 6 of 9*

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 16:17

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:17
02 Std. Gas	0.080	16:18
03 Room Air	0.000	16:18
04 Std. Gas	0.080	16:18
05 Room Air	0.000	16:19
06 Std. Gas	0.079	16:19
07 Room Air	0.000	16:20

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 777777

Deb Kashur

Operator Signature
DEB KASHUR

Remarks:

Calibration check

Form 106-I8000

80-004201
P. 7/29/09

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 16:25

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:25
02 Std. Gas	0.080	16:26
03 Room Air	0.000	16:26
04 Std. Gas	0.080	16:26
05 Room Air	0.000	16:27
06 Std. Gas	0.080	16:27
07 Room Air	0.000	16:28

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 777777

Deb Kashur
Operator Signature
DEB KASHUR

Remarks:

Calibration check

Form 106-I8000

80-004201
P. 8 of 9 JTW

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = TOXL 8164.13.00 06/09
12/22/2009 16:28

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:29
02 Std. Gas	0.079	16:29
03 Room Air	0.000	16:30
04 Std. Gas	0.080	16:30
05 Room Air	0.000	16:30
06 Std. Gas	0.080	16:31
07 Room Air	0.000	16:31

Lot No = 659358
Cyl No = 17
Exp Date = 09/03/2011
County = 08 Oper No. = 777777

Deb Kashur
Operator Signature
DEB KASHUR

Remarks:

Form 106-I8000

80-004201
P. 9 of 9 JEM