



ETHANOL BREATH STANDARD CYLINDER REPORT

OFFICE OF ATTORNEY GENERAL

CRIME LABORATORY DIVISION

Toxicology Section/Breath Alcohol Program

SFN 59282 (5-2011)

Chemical Test Operator Name (Print) <i>Chad Harmon</i>		
Location <i>KYPD</i>	Intoxilyzer® Serial Number <i>80-004194</i>	
Gas Lot Number <i>29614080A2</i>	Gas Cylinder Number <i>26</i>	Gas Expiration Date <i>12/05/2016</i>

Check When Done:

1. Scan/Enter Gas Cylinder Information (Level 1, Function S).

2. Perform an ACA Test (Level 1, Function C).

Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).

A. 0. *084* AC

B. 0. *084* AC

C. 0. *083* AC

3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.

4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.

5. Send the Following to the Crime Laboratory:

A. Ethanol Breath Standard Cylinder Report (SFN59282, Form 105-G).

B. ACA Test Record.

Chemical Test Operator Signature <i>[Signature]</i>	Date <i>12-31-15</i> ✓
Reviewed By (Crime Laboratory Use Only) <i>Deb Shanaver</i>	Date <i>19 Feb 16</i>

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004194
Location = FYPD 8164.13.00 06/09
12/31/2015 16:30

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:31
02 Std. Gas	0.084	16:32
03 Room Air	0.000	16:32
04 Std. Gas	0.084	16:33
05 Room Air	0.000	16:33
06 Std. Gas	0.083	16:33
07 Room Air	0.000	16:34

Lot No = 29614080A2
Cyl No = 26
Exp Date = 12/05/2016
County = 61 Oper No. = 010606



Operator Signature
CHAD HARMON

Remarks: