BREATH ALCOHOL EQUIPMENT REPAIR AND MAINTENANCE

Instr	ument: <u>8000 </u>	Serial Number:	30-005938 Loc	cation: ToxL	
LOAT INO	ERATOR FILE), PROBAD DED **. DAT FILE AND TRUMENT WAS PLACE	AMERROR IN SLY DUE TO IN CONLO NOT I D IN CLASSROOM	med: FIELD. DISCOVERED COMPLETE UPLOAD, REPLICATE ERROR ON M FROM OCTOBER TO MA	AT LOCATION: OCT. 12,13,14,2016. ARCH (13 CLASSES).	
BAT	TELY WAS REPLACED	F TRAINES	. CER 3/28/17		
2.	Does the instrument require further testing? If Yes, continue. If No, sign and date below.			Yes or No	
3.	If needed, set the time	If needed, set the time, date, and location.			
4.	Does the flow need to be calibrated? If Yes, attach paperwork.			Yes o No	
5.	Does the optical bench need to be calibrated? If Yes, attach test records. Yes o No				
6.	Does the gas regulator	need to be calib	orated?	Yes or No	
	PSI gauge reading	-750 PSI	lisplay reading <u>77</u> 6	2	
Set in	nstrument to Wet Bath ar	nd run tests:			
7.	Low AC. Use < 0.03 A	C in ACA mode.	Attach test record.		
	Sim SN:	_ Lot #:	AC:		
8.	Linearity Test. Use ≥ 0).25 AC in ACA r	mode. Attach test record		
	Sim SN:	_ Lot #:	AC:		

OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION

Toxicology Section/Breath Alcohol Program

BrW-005

Set in	strument to Gas	and run tests:					
9.	Print test. Attach test record.						
10.	Interferent Check. Use a 0.10 AC ethanol plus 0.05% acetone in ABA mode. Attach test record.						
	Sim SN:	Lot #:	AC:				
11.	RFI Check. Run in CMS mode. Key radio during any room air or subject test. Attach test record.						
12.	Calibration Check. Use a valid 0.080 AC Ethanol Gas Standard. Run three calibration sets in ACA mode. Attach test records.						
Lot No.		Cylinder No	Expiration Date				
	Test 1	Test 1	Test 1				
	Test 2	Test 2					
	Test 3						
Avera	ge						
Field Inspector's Signature 3/28/17 Date							
D	Jeb Shana	ive	28 Mar 17				
Revie	wed by		Date				