



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT

OFFICE OF ATTORNEY GENERAL

CRIME LABORATORY DIVISION

Toxicology Section/Breath Alcohol Program

SFN 59281 (10-2016)

Serial Number 80-005361	Instrument Location Mountrail County Jail
Reason for Install/Repair <input type="checkbox"/> Install After Receiving From Crime Laboratory <input checked="" type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Properly Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. RFI Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if Any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer®8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Gas cyl info is ok DS crime lab

Field Inspector Signature 	Date 4/14/17
------------------------------------------------------------------------------------------------------------------	------------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified by Deb Shanaver	Certified Date 04 May 17
----------------------------------------------	------------------------------------

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005361
Location = MOSO 8164.14.00 09/16
04/14/2017 16:58

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:59
02 Std. Gas	0.080	17:00
03 Room Air	0.000	17:00
04 Std. Gas	0.080	17:00
05 Room Air	0.000	17:01
06 Std. Gas	0.081	17:01
07 Room Air	0.000	17:02

Lot No = 19615080A5
Cyl No = 16
Exp Date = 09/05/2017
County = 31 Oper No. = 120513



Operator Signature
TRAVIS SKAR

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005361
Location = MOSO 8164.14.00 09/16
04/14/2017 17:02

DRY CAL CHECK

Test	AC	Time
01 Room Air	RFI*	17:03
02 Room Air	0.000	17:03

*Invalid Test
Inhibited - RFI

Lot No = 19615080A5
Cyl No = 16
Exp Date = 09/05/2017
County = 31 Oper No. = 120513



Operator Signature
TRAVIS SKAR

Remarks: