

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-003058

Location: TOXL

- |   | Check When Done |
|---|-----------------|
| A. Is the breath tube heated?   | <u>X</u>        |
| B. Is the diagnostic check complete?  | <u>X</u>        |
| C. Is the time, date, and year correct (re-set if necessary)?   | <u>X</u>        |
| D. Print test. (Attach test record.)  | <u>X</u>        |
| E. Low AC. Use $\leq 0.03$ AC in ACA mode. (Attach test record.)<br>Sim SN <u>DR3846</u> Lot No. <u>160200</u> Rep AC <u>0.020</u>  | <u>X</u>        |
| F. High AC. Use $\geq 0.25$ AC in ACA mode. (Attach test record.)<br>Sim SN <u>DR7111</u> Lot No. <u>16060</u> Rep AC <u>0.300</u>  | <u>X</u>        |
| G. Interferent check. Use 0.05% acetone in ethanol solution in ABA mode.<br>(Attach test record.)<br>Sim SN <u>G10329</u> Lot No. <u>1054</u> Solution <u>0.10 AC + 0.05% Acetone</u> | <u>X</u>        |
| H. RFI check. Run CMS mode. Key radio during CMS test.<br>(Attach test record.)   | <u>X</u>        |
| I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.<br>(Attach test record.) <u>0.080</u> AC   |                 |

Lot No. 17316080A2 Cyl No. 27 Exp. Date 5 Aug 18

Test 1 0.080 Test 4 0.080 Test 7 0.080

Test 2 0.080 Test 5 0.080 Test 8 0.080

Test 3 0.081 Test 6 0.080 Test 9 0.080

Average 0.080

J. Tank pressure, Level 3 Func D Sub G

P: 767 psi Reg 750 psi <sup>775 psi</sup>

K. Remarks/Maintenance record: Annual inspection.

18 Apr 17  
Date

Roberta Guzman Nimmo  
Field Inspector's Signature

18 Apr 17  
Date

Db Shanaver  
Reviewed by

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-003058  
Location = TOXL                    8164.14.00 09/16  
04/18/2017                            10:29

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

Current Instrument Setup


Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?        No  
Display Volume?            No  
Disable On Memfull?        Yes  
# of Print Copies?         1  
Select Std (D/W/I)?        Dry  
Standard Value?            0.080  
Standard Lot #?            17316080A2  
Standard Cyl #?            27  
Standard Expiration?       08/05/2018  
Oper No?                    888888

Flow Cal. Date:            04/07/2008  
Slope                      698  
Intercept                  -678970

IR Calibration Date:       04/15/2015  
                                    3um                      9um

-----  
0th Coef(\*100):        -18934                    -10185  
1st Coef(\*100):        272557                    131040  
2nd Coef(\*100):        2331                      1317  
H2O adj(mg/l\*10k):    333                        280

\*\*\*\*\* Printer Test End \*\*\*\*\*

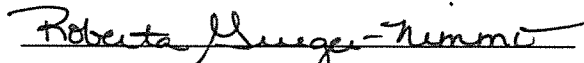

Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks:            *Print Test*

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003058  
Location = TOXL      8164.14.00 09/16  
04/18/2017      10:30

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:31
02 Std. Sol.	0.021	10:31
03 Room Air	0.000	10:32
04 Std. Sol.	0.020	10:33
05 Room Air	0.000	10:33
06 Std. Sol.	0.020	10:34
07 Room Air	0.000	10:34

08 Sim Temp = 34.0°C

Simul Ser No = DR3846

Std Sol No = 16200

County = 08

Oper No. = 888888



Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: Low AC  
0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003058  
Location = TOXL      8164.14.00 09/16  
04/18/2017      10:35

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:36
02 Std. Sol.	0.299	10:37
03 Room Air	0.000	10:37
04 Std. Sol.	0.301	10:38
05 Room Air	0.000	10:38
06 Std. Sol.	0.300	10:39
07 Room Air	0.000	10:40

08 Sim Temp = 34.0°C

Simul Ser No = DR7111  
Std Sol No = 16060  
County = 08      Oper No. = 888888

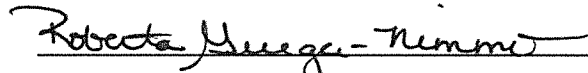  
Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: High AC  
0.300 AC



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer                      Alcohol Analyzer  
North Dakota Model 8000                      SN 80-003058  
Location = TOXL                              8164.14.00 09/16  
04/18/2017    10:45

Test	AC	Time
01 Diagnostic	OK	10:45
02 Room Air	0.000	10:46
03 Subject Test 1	0.000	10:46
04 Room Air	RFI*	10:47
05 Room Air	0.000	10:48

\*Invalid Test  
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT  
Sub DOB = 01/01/1982  
Sub Sex = Female                              Weight = 150  
Test = DUI                                      Cit = RFI CHECK  
Dr. Lic. = ND/DIS821456  
Lot No = 17316080A2  
Cyl No = 27  
Expiration Date = 08/05/2018  
County = 08                                      Oper No. = 888888

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

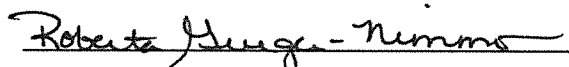

Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: RFI Check


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-003058  
Location = TOXL                    8164.14.00 09/16  
04/18/2017                            10:49

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:49
02 Std. Gas	0.080	10:50
03 Room Air	0.000	10:50
04 Std. Gas	0.080	10:50
05 Room Air	0.000	10:51
06 Std. Gas	0.081	10:51
07 Room Air	0.000	10:52

Lot No = 17316080A2  
Cyl No = 27  
Exp Date = 08/05/2018  
County = 08                            Oper No. = 888888

  
Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check  
                  0.080 AC


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-003058  
Location = TOXL                    8164.14.00 09/16  
04/18/2017                            10:52

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:52
02 Std. Gas	0.080	10:53
03 Room Air	0.000	10:53
04 Std. Gas	0.080	10:54
05 Room Air	0.000	10:54
06 Std. Gas	0.080	10:54
07 Room Air	0.000	10:55

Lot No = 17316080A2  
Cyl No = 27  
Exp Date = 08/05/2018  
County = 08                            Oper No. = 888888

  
Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check  
            0.080 AC




Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-003058  
Location = TOXL      8164.14.00 09/16  
04/18/2017      10:55

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:56
02 Std. Gas	0.080	10:56
03 Room Air	0.000	10:57
04 Std. Gas	0.080	10:57
05 Room Air	0.000	10:57
06 Std. Gas	0.080	10:58
07 Room Air	0.000	10:58

Lot No = 17316080A2  
Cyl No = 27  
Exp Date = 08/05/2018  
County = 08      Oper No. = 888888

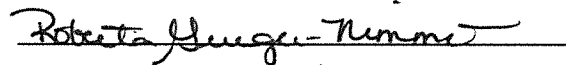  
Operator Signature  
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check  
0.080 AC