

INTOXILYZER® 8000 INSPECTION

Serial No.: 80-005357

Location: UND PD

- Check When Done
- A. Is the breath tube heated? ✓
- B. Is the diagnostic check complete? ✓
- C. Is the time, date, and year correct (re-set if necessary)? ✓
- D. Print test. (Attach test record.) ✓
- E. Low AC. Use  $\leq 0.03$  AC in ACA mode. (Attach test record.)  
Sim SN DR5113 Lot No. 201507 F Rep AC 0.020 ✓
- F. High AC. Use  $\geq 0.25$  AC in ACA mode. (Attach test record.)  
Sim SN DR3378 Lot No. 16060 Rep AC 0.300 ✓
- G. Interferent check. Use 0.05% acetone in ethanol solution in ABA mode.  
(Attach test record.)  
Sim SN DR3803 Lot No. 1C55 Solution 0.10 AC + 0.05% acetone ✓
- H. RFI check. Run CMS mode. Key radio during CMS test.  
(Attach test record.) ✓
- I. Calibration Check. Use Ethanol Breath Standard cylinder. Do three sets in ACA mode.  
(Attach test record.) 0.080 AC

Lot No. 19615080AS Cyl No. 21 Exp. Date 09/05/2017

Test 1 0.080 Test 4 0.081 Test 7 0.080

Test 2 0.081 Test 5 0.080 Test 8 0.081

Test 3 0.081 Test 6 0.080 Test 9 0.081

Average 0.081

J. Tank pressure, Level 3 Func D Sub G P: 504 psi Reg 500 psi

K. Remarks/Maintenance record: annual inspection

25 May 17  
Date

30 MAY 2017  
Date

Deb Shanaver  
Field Inspector's Signature

Charles E. E...  
Reviewed by

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-005357  
Location = UNDP                8164.14.00 09/16  
05/25/2017                      08:29

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

Current Instrument Setup

Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?        No  
Display Volume?            No  
Disable On Memfull?        Yes  
# of Print Copies?         1  
Select Std (D/W/I)?        Dry  
Standard Value?            0.080  
Standard Lot #?            19615080A5  
Standard Cyl #?            21  
Standard Expiration?       09/05/2017  
Oper No?                    777777

Flow Cal. Date:            08/17/2011  
Slope                      663  
Intercept                  -589414

IR Calibration Date:       04/07/2015

	3um	9um
0th Coef(*100):	-16389	-26777
1st Coef(*100):	269467	141875
2nd Coef(*100):	1088	917
H2O adj(mg/l*10k):	159	330

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
DEB SHANAVER

Remarks:

*Print test*

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005357  
Location = UNDP      8164.14.00 09/16  
05/25/2017      08:48

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:49
02 Std. Sol.	0.018	08:50
03 Room Air	0.000	08:51
04 Std. Sol.	0.018	08:51
05 Room Air	0.000	08:52
06 Std. Sol.	0.017	08:52
07 Room Air	0.000	08:53

08 Sim Temp = 34.0°C

Simul Ser No = DR5113  
Std Sol No = 201507F  
County = 18      Oper No. = 777777

*Deb Shanaver*

Operator Signature  
DEB SHANAVER

Remarks:

*Low AC      0.020AC*

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-005357  
Location = UNDP                8164.14.00 09/16  
05/25/2017                        08:55

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:55
02 Std. Sol.	0.305	08:56
03 Room Air	0.000	08:57
04 Std. Sol.	0.301	08:57
05 Room Air	0.000	08:58
06 Std. Sol.	0.300	08:59
07 Room Air	0.000	08:59

08 Sim Temp = 34.0°C

Simul Ser No = DR3378  
Std Sol No = 16060  
County = 18                        Oper No. = 777777



Operator Signature  
DEB SHANAVER

Remarks:

*High AC 0.300AC*

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005357  
Location = UNDP      8164.14.00 09/16  
05/25/2017      09:00

Test	AC	Time
01 Room Air	0.000	09:02
02 *Subject Test	INT*	09:02
03 Room Air	0.000	09:03

\*Invalid Test  
Interferent Detected

Sub Name = ~~RPI, TEST NA~~ <sup>DS</sup> Interferent  
Sub DOB = 09/23/1980  
Sub Sex = Unknown      Weight = NA  
Test = DUI      Cit = NA  
Dr. Lic. = ND/NA  
Simul Ser No = DR3378  
Std Sol No = 16060  
County = 18      Oper No. = 777777

Deb Shanaver

Operator Signature  
DEB SHANAVER

Remarks:

Interferent detected check

Form 106-I8000

0.10AC + 0.05% acetone

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-005357  
Location = UNDP                8164.14.00 09/16  
05/25/2017                        09:05

Test	AC	Time
01 Diagnostic	OK	09:06
02 Room Air	0.000	09:06
03 *Subject Test	RFI*	09:07
04 Room Air	0.000	09:07

\*Invalid Test  
  Inhibited - RFI

Sub Name = RFI, TEST NA  
Sub DOB = 03/27/1998  
Sub Sex = Unknown                Weight = NA  
Test = DUI                         Cit = NA  
Dr. Lic. = ND/NA  
Simul Ser No = DR3378  
Std Sol No = 16060  
County = 18                        Oper No. = 777777

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

Deb Shanaver

Operator Signature  
DEB SHANAVER

Remarks:

RFI test

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005357  
Location = UNDP      8164.14.00 09/16  
05/25/2017      08:32

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:33
02 Std. Gas	0.080	08:33
03 Room Air	0.000	08:34
04 Std. Gas	0.081	08:34
05 Room Air	0.000	08:35
06 Std. Gas	0.081	08:35
07 Room Air	0.000	08:36

Lot No = 19615080A5  
Cyl No = 21  
Exp Date = 09/05/2017  
County = 18      Oper No. = 777777



Operator Signature  
DEB SHANAVER

Remarks:

Calibration check 0.080AC

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005357  
Location = UNDP      8164.14.00 09/16  
05/25/2017      08:37

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:37
02 Std. Gas	0.081	08:38
03 Room Air	0.000	08:38
04 Std. Gas	0.080	08:39
05 Room Air	0.000	08:39
06 Std. Gas	0.080	08:40
07 Room Air	0.000	08:40

Lot No = 19615080A5  
Cyl No = 21  
Exp Date = 09/05/2017  
County = 18      Oper No. = 777777

Deb Shanaver  
Operator Signature  
DEB SHANAVER

Remarks:

Calibration check 0.080AC



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005357  
Location = UNDP      8164.14.00 09/16  
05/25/2017      08:40

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:41
02 Std. Gas	0.080	08:41
03 Room Air	0.000	08:42
04 Std. Gas	0.081	08:42
05 Room Air	0.000	08:43
06 Std. Gas	0.081	08:43
07 Room Air	0.000	08:44

Lot No = 19615080A5  
Cyl No = 21  
Exp Date = 09/05/2017  
County = 18      Oper No. = 777777



Operator Signature  
DEB SHANAVER

Remarks:

*Calibration check 0.080AC*

Form 106-I8000