

INTOXILYZER® 8000 INSPECTION

Intoxilyzer® 8000 Serial Number: 80-004184

Location: TOXL

A. General Setup and Checks:

1. Diagnostic Tests Pass and Instrument in Ready Mode
2. Breath tube heated
3. Date, time and location code (Level 2,E). Re-set if necessary.
4. Print test (Level 1,P). Sign and attach test record.
5. Tank monitor (Level 3,D,G). Display and Regulator within 50 psi.
 - a. Display: 402 psi Regulator: 400 psi

B. Tests (Sign and attach test records):

Configure simulator for the following tests (Level 1,S).

1. Wet Calibration Check

- a. Low AC (Level 1,C): Known Value ≤ 0.03 AC: 0.030 AC
Sim. Ser #: MP3060 Lot #: 16200 Exp. Date: 17 Aug 18
- b. High AC (Level 1,C): Known Value ≥ 0.25 AC: 0.300 AC
Sim. Ser #: MP3062 Lot #: 17350 Exp. Date: 11 Oct 19

Configure dry gas standard for the following tests (Level 1,S).

2. Interferent Check (Level 1,B): Known Value: 0.10 AC + 0.05% Acetone
Sim. Ser #: DR5144 Lot #: 1CS5 Exp. Date: 15 May 19
 Display reads "Interferent Detect"
3. RFI Check (CMS Mode) Display reads "RFI Detect"
4. Dry Calibration Check (Level 1,C): Known Value 0.080 AC
Gas Cyl Lot #: 20016080A Cyl #: 48 Exp. Date: 5 Sep 18

Test 1 <u>0.081</u>	Test 4 <u>0.080</u>	Test 7 <u>0.081</u>
Test 2 <u>0.081</u>	Test 5 <u>0.080</u>	Test 8 <u>0.081</u>
Test 3 <u>0.080</u>	Test 6 <u>0.081</u>	Test 9 <u>0.081</u>
Average <u>0.081</u>		

C. Remarks/Maintenance: Annual inspection. Battery was replaced and trained. Calibration ~~test~~ completed take of tank sensor. Replaced O-rings on simulator return, breath hose

Instrument is acceptable to be used in the field.

Roberto Inequi-Tummo
Breath Analyst Signature

9 Apr 18
Date

Charles E. Edr
Reviewed by

09 APR 2018
Date

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:18

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 20016080A2
Standard Cyl #? 48
Standard Expiration? 09/05/2018
Oper No? 888888

Flow Cal. Date: 07/30/2010
Slope 804
Intercept -925008

IR Calibration Date: 12/07/2015
3um 9um

	3um	9um
0th Coef(*100):	-15679	-18680
1st Coef(*100):	263558	136948
2nd Coef(*100):	2594	1146
H2O adj(mg/l*10k):	405	327

***** Printer Test End *****

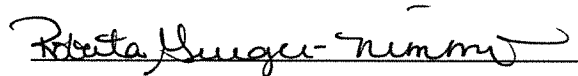

Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Print Test

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:50

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:50
02 Std. Sol.	0.021	10:51
03 Room Air	0.000	10:52
04 Std. Sol.	0.021	10:52
05 Room Air	0.000	10:53
06 Std. Sol.	0.021	10:54
07 Room Air	0.000	10:54

08 Sim Temp = 34.0°C

Simul Ser No = MP3060
Std Sol No = 16200
County = 08 Oper No. = 888888

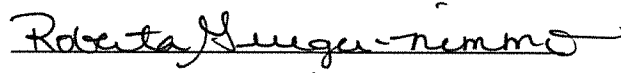
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: low AC
 0.020 AC

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:56

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:57
02 Std. Sol.	0.297	10:58
03 Room Air	0.000	10:58
04 Std. Sol.	0.297	10:59
05 Room Air	0.000	11:00
06 Std. Sol.	0.299	11:01
07 Room Air	0.000	11:01

08 Sim Temp = 34.0°C

Simul Ser No = MP3062
Std Sol No = 17350
County = 08 Oper No. = 888888

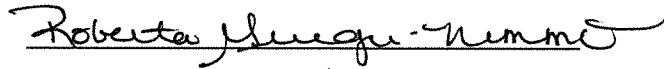
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: High AC
 0.0 Res 0.300 AC


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 11:08

Test	AC	Time
01 Room Air	0.000	11:09
02 *Subject Test	INT*	11:09
03 Room Air	0.000	11:10

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS821456
Lot No = 20016080A2
Cyl No = 48
Expiration Date = 09/05/2018
County = 08 Oper No. = 888888

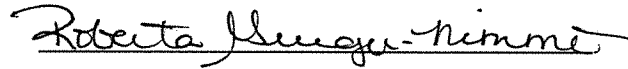

Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Interferent Check
0.080 AC RW
0.10 AC + 0.05% Acetone
Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 11:11

Test	AC	Time
01 Diagnostic	OK	11:11
02 Room Air	RFI*	11:12
03 Room Air	0.000	11:12

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = RFI CHECK
Dr. Lic. = ND/DIS821456
Lot No = 20016080A2
Cyl No = 48
Expiration Date = 09/05/2018
County = 08 Oper No. = 888888

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

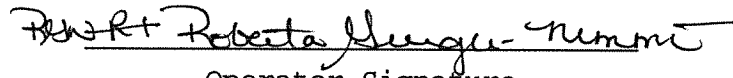

Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: RFI Check


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:20

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:21
02 Std. Gas	0.081	10:21
03 Room Air	0.000	10:22
04 Std. Gas	0.081	10:22
05 Room Air	0.000	10:23
06 Std. Gas	0.080	10:23
07 Room Air	0.000	10:24

Lot No = 20016080A2
Cyl No = 48
Exp Date = 09/05/2018
County = 08 Oper No. = 888888

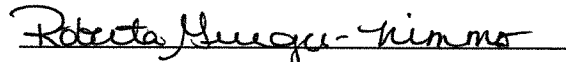
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Calibration Check
0.080 AC


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:26

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:27
02 Std. Gas	0.080	10:27
03 Room Air	0.000	10:28
04 Std. Gas	0.080	10:28
05 Room Air	0.000	10:29
06 Std. Gas	0.081	10:29
07 Room Air	0.000	10:30

Lot No = 20016080A2
Cyl No = 48
Exp Date = 09/05/2018
County = 08 Oper No. = 888888

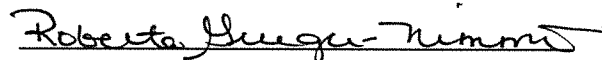
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: *Calibration Check*
0.080 AC


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004184
Location = TOXL 8164.14.00 09/16
04/09/2018 10:30

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:31
02 Std. Gas	0.081	10:31
03 Room Air	0.000	10:32
04 Std. Gas	0.081	10:32
05 Room Air	0.000	10:33
06 Std. Gas	0.081	10:33
07 Room Air	0.000	10:34

Lot No = 20016080A2
Cyl No = 48
Exp Date = 09/05/2018
County = 08 Oper No. = 888888

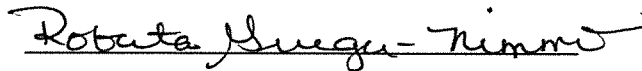

Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: *Calibration Check*
0.080 AC

Form 106-I8000