

INTOXILYZER® 8000 CALIBRATION ADJUSTMENT

Intoxilyzer® 8000 Serial Number: 80-002666 Location: TOXL

- A. Flow Sensor Calibration and Verification Check (Level 3,M,C,F)
1. Replaced o-rings if damaged
 2. Flow Meter Serial Number: 40655
 3. Air Supplied to Intoxilyzer® 8000 at:
 - a. 5 L/min 15 L/min 30 L/min
 4. Flow Rate Calibration Printout Attached
 - a. Correlation ≥ 0.99000
 5. Flow Sensor Calibration Verification (Level 3,D,F)
 - a. 10 L/min: 0. 16.0 L/S X 60 Sec/min = 9.60 L/min
 - b. 20 L/min: 0. 33.2 L/S X 60 Sec/min = 19.92 L/min
 - c. Flow Rates within ± 1 L/min of Expected Value
- B. Gas Tank Sensor Check (Level 3,D,G)
1. Display: 1027 psi Regulator: 1000 psi
 2. Display and Regulator within 50 psi
 3. Completed tare of tank sensor if needed (Level 3,M,C,G)
- C. Optical Bench Calibration and Verification Check (Level 3,M,C,O)
1. Autocalibration Printout Attached
 - a. Max Power Res Value ≥ 10
 - b. Auto Range Res Value ≥ 4
 2. Simulator Solutions for Optical Bench Calibration Adjustment
 - a. Set # Solutions to Run at 5

Soln.	g/210 L	Lot No.	Exp. Date	Simulator SN
1	0.000	NA – MilliQ H ₂ O	NA – MilliQ H ₂ O	MP3002
2	0.040	201706B	6 Jun 19	MP3057
3	0.080	201707E	25 Jul 19	MP3058
4	0.150	201705D	24 May 19	MP3059
5	0.300	201803H	22 Mar 20	MP3061

3. 0.100 AC Calibration Gas for H₂O Adjustment
 - a. Lot No. 13518100A3 Cyl No. 6 Exp. Date: 5 Aug 20
4. Atmospheric Pressure
 - a. 962 mbar Displayed by Intoxilyzer® 8000
 - b. 962 mbar Adjusted to using barometer
 - c. 961 mbar on Auto Calibration Report printout
5. Screen displayed "Calibration Success"

6. Calibration Adjustment Printout Attached
- Solution 1 Avg % Abs ≤ 0.2500
 - Solution 2-5 REL STD DEV ≤ 3.000
 - Residual (g/210 L) Values for Solutions 1-5 ≤ 0.0020 for 3 μm and 9 μm channels
 - Dry Gas H₂O Adjustment Sum for 3 μm and 9 μm channels within ± 10

	Average ^{3μm}	H ₂ O Adjust	
3 μm	<u>4276.6665</u> + <u>961485</u>	= <u>5237.6665</u>	²⁵⁰ 4761.6665
9 μm	<u>4507.0000</u> + <u>254</u>	= <u>4761</u>	

7. Optical Bench Calibration Verification (Level 1, S and C)
- Wet Calibration Check
 - Low AC Known Value ≤ 0.03 AC: 0.030 AC
 Sim. SN: DR5188 Lot No.: 18020 Exp. Date: 9 Jan 20
 - High AC Known Value ≥ 0.25 AC: 0.250 AC
 Sim. SN: DR3374 Lot No.: 2018036 Exp. Date: 22 Mar 20
 - Dry Calibration Check: Known Value 0.08 AC
 Lot No. 34A17080A3 Cyl No. 28 Exp. Date: 5 Feb 20
 Test 1 0.079 AC Test 4 0.079 AC Test 7 0.079 AC
^{Res} AC
 Test 2 0.078 AC Test 5 0.078 AC Test 8 0.079 AC
^{Res} AC
 Test 3 0.079 AC Test 6 0.079 AC Test 9 0.079 AC
 Average 0.079 AC
 - Wet Calibration Check and Dry Calibration Check AC results are within ± 0.005 or $\pm 5\%$ (whichever is greater) of stated value.

D. Remarks/Maintenance: Changed o-ring on breath hose, and simulator return. Completed calibration adjustment to check steps for calibration adjustment SOP. Display intensity is decreasing and barcode scanner has failed. Due to extent of repairs, instrument will be disposed of.

Instrument is acceptable to be used in the field.

Roberta Muegge-Nimmo
 Breath Analyst Signature
Charles E. Ehr
 Reviewed by

5 Feb 19
 Date
2/5/2019
 Date

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 10:28

Flow Rate Calibration*****

- 1: Rate (Liters/min) = 5
 SQRT(Diff)) = 6.082
- 2: Rate (Liters/min) = 15
 SQRT(Diff)) = 11.223
- 3: Rate (Liters/min) = 30
 SQRT(Diff)) = 22.133

Dependent Data Scale Factor = 100000 L/min

Independent Data Scale Factor = 256

Rounded Slope = 597

Rounded Intercept = -343366

Correlation = 0.99598

TOXL
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-002666
02/05/2019 12:07:27

Auto Calibration
Max Power Res Value = 45
Auto Range Res Value = 29

TOXL
 Intoxilyzer - Alcohol Analyzer
 Model 8000 SN 80-002666
 02/05/2019 12:07:27

Auto Calibration

pg 1 of 2

<<<<< 3um >>>>>			<<<<< 9um >>>>>		

Solution = 0.000 g/210L or 0.0000 mg/l, Samples = 4, Discarded = 1					
Sample	% Abs	(% Abs Ref)	% Abs	(% Abs Ref)	
Sample #1	0.1420	(-0.0170)	0.1240	(-0.0070)	
Sample #2	0.0970	(0.0380)	0.1170	(0.0020)	
Sample #3	0.0790	(0.0850)	0.1130	(0.0180)	
Sample #4	0.1060	(0.0860)	0.1040	(0.0200)	
Avg % Abs	0.0940	(0.0697)	0.1113	(0.0133)	
STD DEV	0.0137	(0.0274)	0.0067	(0.0099)	
REL STD DEV	14.625	(39.371)	5.981	(73.993)	

Solution = 0.040 g/210L or 0.1905 mg/l, Samples = 4, Discarded = 1					
Sample	% Abs	(% Abs Ref)	% Abs	(% Abs Ref)	
Sample #1	0.7620	(0.0060)	1.4670	(0.0100)	
Sample #2	0.8020	(0.0060)	1.4920	(0.0140)	
Sample #3	0.7910	(0.0260)	1.4980	(0.0160)	
Sample #4	0.7980	(0.0250)	1.5090	(0.0140)	
Avg % Abs	0.7970	(0.0190)	1.4997	(0.0147)	
STD DEV	0.0056	(0.0113)	0.0086	(0.0012)	
REL STD DEV	0.699	(59.313)	0.575	(7.873)	

Solution = 0.080 g/210L or 0.3810 mg/l, Samples = 4, Discarded = 1					
Sample	% Abs	(% Abs Ref)	% Abs	(% Abs Ref)	
Sample #1	1.4570	(0.0120)	2.8510	(0.0060)	
Sample #2	1.5070	(0.0090)	2.8840	(0.0000)	
Sample #3	1.4640	(0.0340)	2.8830	(0.0060)	
Sample #4	1.5100	(0.0220)	2.8900	(0.0060)	
Avg % Abs	1.4937	(0.0217)	2.8857	(0.0040)	
STD DEV	0.0257	(0.0125)	0.0038	(0.0035)	
REL STD DEV	1.723	(57.708)	0.131	(86.603)	

Solution = 0.150 g/210L or 0.7143 mg/l, Samples = 4, Discarded = 1					
Sample	% Abs	(% Abs Ref)	% Abs	(% Abs Ref)	
Sample #1	2.6040	(0.0020)	5.0660	(-0.0010)	
Sample #2	2.6460	(0.0070)	5.1240	(0.0050)	
Sample #3	2.6810	(-0.0040)	5.1490	(-0.0010)	
Sample #4	2.6660	(0.0080)	5.1590	(-0.0010)	
Avg % Abs	2.6643	(0.0037)	5.1440	(0.0010)	
STD DEV	0.0176	(0.0067)	0.0180	(0.0035)	
REL STD DEV	0.659	(181.591)	0.350	(346.410)	

Solution = 0.300 g/210L or 1.4286 mg/l, Samples = 4, Discarded = 1					
Sample	% Abs	(% Abs Ref)	% Abs	(% Abs Ref)	
Sample #1	5.0460	(-0.0140)	9.6820	(-0.0090)	
Sample #2	5.0190	(0.0210)	9.6960	(0.0140)	
Sample #3	5.0750	(0.0040)	9.7240	(0.0150)	
Sample #4	5.0850	(0.0040)	9.7490	(0.0140)	
Avg % Abs	5.0597	(0.0097)	9.7230	(0.0143)	
STD DEV	0.0356	(0.0098)	0.0265	(0.0006)	
REL STD DEV	0.703	(101.534)	0.273	(4.028)	

TOXL
 Intoxilyzer - Alcohol Analyzer
 Model 8000 SN 80-002666
 02/05/2019 12:07:27

Auto Calibration

pg 2 of 2

<<<<< 3um >>>>>			<<<<< 9um >>>>>		
-----			-----		
Zero Order Coef	-249.62		Zero Order Coef	-147.94	
First Order Coef	2661.62		First Order Coef	1336.08	
Second Order Coef	41.75		Second Order Coef	15.28	
-----			-----		
Act	Fit	Residual	Act	Fit	Residual
(g/210L)	(g/210L)	(g/210L)	(g/210L)	(g/210L)	(g/210L)
0.000	0.000	-0.0000	0.000	0.000	-0.0000
0.040	0.040	0.0001	0.040	0.040	0.0003
0.080	0.080	-0.0002	0.080	0.081	-0.0005
0.150	0.150	0.0001	0.150	0.150	0.0003
0.300	0.300	-0.0000	0.300	0.300	-0.0000
-----			-----		

<<<<< 3um >>>>>		<<<<< 9um >>>>>	
-----		-----	
Solution = 0.100 g/210L or 0.4762 mg/l, Samples = 4, Discarded = 1			
Sample			
Sample #1	4295.00	4475.00	
Sample #2	4206.00	4475.00	
Sample #3	4353.00	4516.00	
Sample #4	4271.00	4530.00	
Avg	4276.6665	4507.0000	
STD DEV	73.6637	28.5832	
REL STD DEV	1.722	0.634	
H2O adjust (mg/l*10k)	485	254	

Atmospheric Pressure = 961

*****CALIBRATION SUCCESSFUL*****

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 12:52

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:53
02 Std. Sol.	0.021	12:54
03 Room Air	0.000	12:54
04 Std. Sol.	0.021	12:55
05 Room Air	0.000	12:56
06 Std. Sol.	0.021	12:56
07 Room Air	0.000	12:57

08 Sim Temp = 34.0°C

Simul Ser No = DR5188

Std Sol No = 18020

County = 08

Oper No. = 888888

Roberta Grieger-Nimmo

Operator Signature

ROBERTA GRIEGER-NIMMO

Wet calibration Check

Remarks: Low AC

0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 12:58

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:59
02 Std. Sol.	0.250	13:00
03 Room Air	0.000	13:01
04 Std. Sol.	0.250	13:01
05 Room Air	0.000	13:02
06 Std. Sol.	0.251	13:03
07 Room Air	0.000	13:03


08 Sim Temp = 34.0°C

Simul Ser No = DR3374

Std Sol No = 201803G

County = 08

Oper No. = 888888

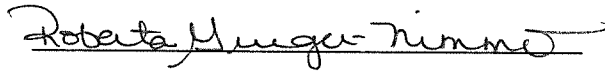

Operator Signature

ROBERTA GRIEGER-NIMMO

Wet Calibration Check

Remarks: High AC

0.250 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 13:05

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:06
02 Std. Gas	0.079	13:06
03 Room Air	0.000	13:06
04 Std. Gas	0.078	13:07
05 Room Air	0.000	13:07
06 Std. Gas	0.079	13:08
07 Room Air	0.000	13:08

Lot No = 34917080A3
Cyl No = 028
Exp Date = 02/05/2020
County = 08 Oper No. = 888888

Roberta Grieger-Nimmo
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: *Dry calibration check*
0.080AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 13:08

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:09
02 Std. Gas	0.079	13:10
03 Room Air	0.000	13:10
04 Std. Gas	0.078	13:10
05 Room Air	0.000	13:11
06 Std. Gas	0.079	13:11
07 Room Air	0.000	13:12

Lot No = 34917080A3
Cyl No = 028
Exp Date = 02/05/2020
County = 08 Oper No. = 888888

Roberta Grieger-Nimmo
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: *Dry Calibration Check*
0.080 AC

Form 106-I8000


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-002666
Location = TOXL 8164.14.00 09/16
02/05/2019 13:12

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:13
02 Std. Gas	0.079	13:13
03 Room Air	0.000	13:14
04 Std. Gas	0.079	13:14
05 Room Air	0.000	13:15
06 Std. Gas	0.079	13:15
07 Room Air	0.000	13:16

Lot No = 34917080A3
Cyl No = 028
Exp Date = 02/05/2020
County = 08 Oper No. = 888888

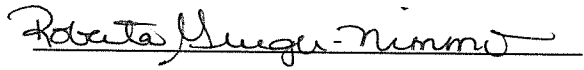
Operator Signature
ROBERTA GRIEGER-NIMMO

Remarks: Dry Calibration Check
0.080AC

Form 106-I8000