

ETHANOL GAS STANDARD CYLINDER REPORT NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM SFN 59282 (08/2018)

| Chemical Test Operator Name (Print) GALLAGHER, JONATHAN | | | | | |
|--|---------------------|----------------------------|---------------------|--|--|
| Location | | Intoxilyzer® Serial Number | | | |
| HARV | | 80-006500 | | | |
| Gas Lot Number | Gas Cylinder Number | | Gas Expiration Date | | |
| 34418080A2 | 061 | | 02/05/2021 | | |

Check When Done:

- X 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- X 2. Perform an ACA Test (Level 1, Function C).

Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).

- A. 0,083 AC
- B. 0.082 AC
- C. 0.082 AC

3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.

- X 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- S. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

| Chemical Test Operator Signature | Date |
|---|------------------|
| the defen | 26 FEB19 |
| Reviewed By (Crime Laboratory Use Only) | Date 27 FFB 2019 |
| | |

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

| CMI, Inc. Intoxilyzer | Alcohol Analyzer |
|-------------------------|------------------|
| North Dakota Model 8000 | SN 80-006500 |
| Location = HARU | 8164.13.00 06/09 |
| 02/26/2019 | 20:03 |

| | DRY CAL CHECK |] |
|---|---|--|
| Test | AC | Time |
| 01 Room Air 02 Std. Gas 03 Room Air 04 Std. Gas 05 Room Air 06 Std. Gas 07 Room Air | 0.000 0.083 0.000 0.082 0.000 0.082 0.000 | 20:03 20:04 20:04 20:04 20:05 20:05 20:05 20:06 |

Lot No = 34418080A2 Cyl No = 61Exp Date = 02/05/2021 County = 52Oper No. = 131749 Operator Signature JONATHAN GALLAGHER

Remarks: