

INTOXILYZER® 8000 INSPECTION

Intoxilyzer® 8000 Serial Number: 80-005951

Location: HEART OF AMERICA CORRECTIONS (RUGB)

A. General Setup and Checks:

1. Diagnostic Tests Pass and Instrument in Ready Mode
2. Breath tube heated
3. Date, time and location code (Level 2,E). Re-set if necessary.
4. Print test (Level 1,P). Sign and attach test record.
5. Tank monitor (Level 3,D,G). Display and Regulator within 50 psi.
 - a. Display: 941 psi Regulator: 950 psi

B. Tests (Sign and attach test records):

Configure simulator for the following tests (Level 1,S).

1. Wet Calibration Check
 - a. Low AC (Level 1,C): Known Value ≤ 0.03 AC: 0.020 AC
Sim. Ser #: DR3375 Lot #: 17160 Exp. Date: 6.8.19
 - b. High AC (Level 1,C): Known Value ≥ 0.25 AC: 0.300 AC
Sim. Ser #: PR7111 Lot #: 17350 Exp. Date: 10.11.19

Configure dry gas standard for the following tests (Level 1,S).

2. Interferent Check (Level 1,B): Known Value: 0.10 AC + 0.05% Acetone
Sim. Ser #: DR7009 Lot #: ICS 5 Exp. Date: NA
 Display reads "Interferent Detect"
3. RFI Check (CMS Mode) Display reads "RFI Detect"
4. Dry Calibration Check (Level 1,C): Known Value 0.080 AC
Gas Cyl Lot #: 13518080A6 Cyl #: 023 Exp. Date: 8.5.20

Test 1 <u>0.082</u>	Test 4 <u>0.081</u>	Test 7 <u>0.081</u>
Test 2 <u>0.082</u>	Test 5 <u>0.081</u>	Test 8 <u>0.082</u>
Test 3 <u>0.081</u>	Test 6 <u>0.081</u>	Test 9 <u>0.081</u>
Average <u>0.081</u>		

C. Remarks/Maintenance: REPLACED BREATH HOSE O-RING
ANNUAL INSPECTION

Instrument is acceptable to be used in the field.

Breath Analyst Signature

NA

Date

5.6.19
NA

Reviewed by

Date

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 10:57

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^*()_+?

Current Instrument Setup

Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 13518080A6
Standard Cyl #? 023
Standard Expiration? 08/05/2020
Oper No? 666666

Flow Cal. Date: 06/21/2013
Slope 693
Intercept -721277

IR Calibration Date: 05/08/2017
 3um 9um

0th Coef(*100): -17201 -17496
1st Coef(*100): 267144 137813
2nd Coef(*100): 2948 1389
H2O adj(mg/l*10k): 438 428

***** Printer Test End *****



Operator Signature
CHARLES EDER

Remarks: PRINT TEST

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:12

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:13
02 Std. Sol.	0.020	11:13
03 Room Air	0.000	11:14
04 Std. Sol.	0.020	11:15
05 Room Air	0.000	11:15
06 Std. Sol.	0.020	11:16
07 Room Air	0.000	11:16


08 Sim Temp = 34.0°C

Simul Ser No = DR3375

Std Sol No = 17160

County = 35

Oper No. = 666666

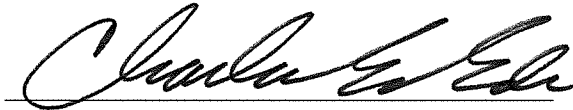

Operator Signature
CHARLES EDER

Remarks:

Low AC
0.020 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:23

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:24
02 Std. Sol.	0.300	11:24
03 Room Air	0.000	11:25
04 Std. Sol.	0.302	11:26
05 Room Air	0.000	11:26
06 Std. Sol.	0.303	11:27
07 Room Air	0.000	11:27

08 Sim Temp = 34.0°C

Simul Ser No = DR7111
Std Sol No = 17350
County = 35 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: *HIGH AC*
0.300 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:08

Test	AC	Time
01 Room Air	0.000	11:09
02 *Subject Test	INT*	11:09
03 Room Air	0.000	11:10

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS821456
Lot No = 13518080A6
Cyl No = 023
Expiration Date = 08/05/2020
County = 35 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: *Interferent Check*
0.10d AC + 0.05% ACETONE
CEE
Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:02

Test	AC	Time
01 Diagnostic	OK	11:03
02 Room Air	0.000	11:03
03 *Subject Test	RFI*	11:03
04 Room Air	0.000	11:04

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 01/01/1982
Sub Sex = Female Weight = 150
Test = DUI Cit = RFI CHECK
Dr. Lic. = ND/DIS821456
Lot No = 13518080A6
Cyl No = 023
Expiration Date = 08/05/2020
County = 35 Oper No. = 666666

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.

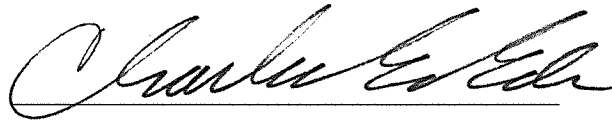

Operator Signature
CHARLES EDER

Remarks:

RFI CHECK

Form 106-I8000


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 10:58

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:59
02 Std. Gas	0.082	10:59
03 Room Air	0.000	11:00
04 Std. Gas	0.082	11:00
05 Room Air	0.000	11:00
06 Std. Gas	0.081	11:01
07 Room Air	0.000	11:01

Lot No = 13518080A6
Cyl No = 023
Exp Date = 08/05/2020
County = 35 Oper No. = 666666

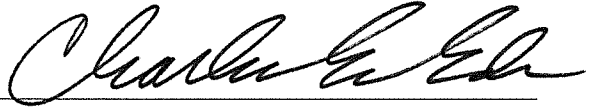

Operator Signature
CHARLES EDER

Remarks: **CALIBRATION CHECK**
0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:04

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:04
02 Std. Gas	0.081	11:05
03 Room Air	0.000	11:05
04 Std. Gas	0.081	11:06
05 Room Air	0.000	11:06
06 Std. Gas	0.081	11:07
07 Room Air	0.000	11:07

Lot No = 13518080A6
Cyl No = 023
Exp Date = 08/05/2020
County = 35 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005951
Location = RUGB 8164.14.00 09/16
05/06/2019 11:29

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:30
02 Std. Gas	0.081	11:30
03 Room Air	0.000	11:31
04 Std. Gas	0.082	11:31
05 Room Air	0.000	11:32
06 Std. Gas	0.081	11:32
07 Room Air	0.000	11:33

Lot No = 13518080A6
Cyl No = 023
Exp Date = 08/05/2020
County = 35

Oper No. = 666666

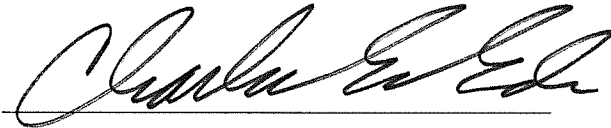

Operator Signature
CHARLES EDER

Remarks: CALIBRATION CHECK
0.080 AC

Form 106-I8000