

INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT

NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM SFN 59281 (06/2018)

Serial Number 86 - 005940	Instrument Location				
Reason for Install/Repair					
☐ Install After Receiving From Crime Laboratory ☐ Install After Location Change					
Other (Specify) Muchine uplate					
Check When Done:					
1. Surge Protector Installed/Property Grounded.					
2. Telephone Line Connected to Intoxilyzer® 8000.					
☑ 3. Breath Tube Heated.					
4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).					
5. Scan/Enter Gas Cylinder Information (Level 1, Function S).					
A. Print Test (Level1, Function P).					
C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).					
7. Repair and/or Maintenance Performed (if any):	/A				
8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.					
9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.					
10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.					
Field Inspector Signature		Date			
Jac Gan Store	Z	2-1-21			
Crime Laboratory Use Only					
This installation has been reviewed and the instrument is approved to be the date the Field Inspector performed the installation. This record on for County of Burleigh, North Dakota, is certified to be a true and correct county.	ile at the Office of Attorney General, Crin	mine alcohol concentration from ne Laboratory Division, in the			

FORM 104-G

Reviewed/Certified By

Certified Date 2 · 3 · 2 /

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005940
Location = NEWT 8164.14.00 09/16
02/01/2021 19:50

******** Printer Test ********

abcdefghijklmnopqrstuvwxyz1234567890-=|
ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=|
ABCDEFGHIJKLMNOPQRSTUVWXYZ!@#\$%^&*()_+?

Current Instrument Setup Enabled Data Entry Mode: DABACABA Start Test Sequence: Display Prelim Rslt? Yes Display Third Digit? Yes Inhib Printer(Y/N)? No Display Volume? No Disable On Memfull? Yes # of Print Copies? 1 Select Std (D/W/I)? Dry Standard Value? 0.080 Standard Lot #? 24119080A1 Standard Cyl #? 26 Standard Expiration? 11/05/2021 132358 Oper No?

Flow Cal. Date: 06/21/2013 Slope 688 Intercept -659053

IR Calibration Date: 04/02/2015

3um 9um

Oth Coef(*100): -18917 -25088

1st Coef(*100): 267616 136158

2nd Coef(*100): 3374 1493

H2O adj(mq/l*10k): 427 417

****** Printer Test End *******

Operator Signature DEMETRIUS JAMERSON

Remarks:

Post test - Install

Form 106-I8000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer
North Dakota Model 8000
Location = NEWT
02/01/2021

Alcohol Analyzer SN 80-005940 8164.14.00 09/16 19:52

DRY CAL CHECK

Τe	est		AC	Time
01	Room	Air	0.000	19:53
02	Std.	Gas	0.080	19:53
03	Room	Air	0.000	19:54
04	Std.	Gas	0.080	19:54
05	Room	Air	0.000	19:55
06	Std.	Gas	0.080	19:55
07	Room	Air	0.000	19:55

Lot No = 24119080A1

Cyl No = 26

Exp Date = 11/05/2021

County = 60

Oper No. = 132358

DEMETRIUS

Operator Signature DEMETRIUS JAMERSON

Remarks:

ACA TEST - Instill

Form 106-I8000

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer
North Dakota Model 8000
Location = NEWT
02/01/2021

Alcohol Analyzer SN 80-005940 8164.14.00 09/16 20:00

Test	AC	Time
01 Room Air	RFI*	20:01
02 Room Air	0.000	20:02

*Invalid Test Inhibited - RFI

Sub Name = JAMERSON, DEMETRIUS JAMAUL

Sub DOB = 03/19/1994

Sub Sex = Male

Weight = 230

Test = DUI

Cit = 00000

Dr. Lic. = ND/JAM944020

Lot No = 24119080A1

Cyl No = 26

Expiration Date = 11/05/2021

County = 60

Oper No. = 132358

Operator Signature DEMETRIUS JAMERSON

Remarks:

RFI test - Install

Form 106-I8000