

INTOXILYZER® 8000 INSPECTION

Intoxilyzer® 8000 Serial Number: 80-005939 Location: TOXL

A. General Setup and Checks:

1. Diagnostic Tests Pass and Instrument in Ready Mode
2. Breath tube heated
3. Date, time and location code (Level 2,E). Re-set if necessary.
4. Print test (Level 1,P). Sign and attach test record.
5. Tank monitor (Level 3,D,G). Display and Regulator within 50 psi.
 - a. Display: 724 psi Regulator: 725 psi

B. Tests (Sign and attach test records):

Configure simulator for the following tests (Level 1,S).

1. Wet Calibration Check
 - a. Low AC (Level 1,C): Known Value ≤ 0.03 AC: 0.030 AC
Sim. Ser #: MP5289 Lot #: 201911E Exp. Date: 11.19.21
 - b. High AC (Level 1,C): Known Value ≥ 0.25 AC: 0.300 AC
Sim. Ser #: MP3069 Lot #: 20030 Exp. Date: 1.21.22

Configure dry gas standard for the following tests (Level 1,S).

2. Interferent Check (Level 1,B): Known Value: 0.10AC + 0.05% Acetone
Sim. Ser #: DR3847 Lot #: ICS6 Exp. Date: NA
 Display reads "Interferent Detect"
3. RFI Check (CMS Mode) Display reads "RFI Detect"
4. Dry Calibration Check (Level 1,C): Known Value 0.080 AC
Gas Cyl Lot #: 05620080A1 Cyl #: 6 Exp. Date: 4.5.22

Test 1	<u>0.080</u>	Test 4	<u>0.079</u>	Test 7	<u>0.080</u>
Test 2	<u>0.079</u>	Test 5	<u>0.080</u>	Test 8	<u>0.079</u>
Test 3	<u>0.079</u>	Test 6	<u>0.079</u>	Test 9	<u>0.079</u>
Average	<u>0.079</u>				

C. Remarks/Maintenance: INSPECTION, AFTER REPAIR, TO BE PLACED ON SUPPLEMENTAL DEVICE LIST.

ANNUAL INSPECTION

Instrument is acceptable to be used in the field.

Charles E. Ehr
Breath Analyst Signature

8.20.2021
Date

NA

NA

Reviewed by

Date

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 15:41

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

Current Instrument Setup

Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 05620080A1
Standard Cyl #? 6
Standard Expiration? 04/05/2022
Oper No? 666666

Flow Cal. Date: 06/18/2020
Slope 611
Intercept -592495

IR Calibration Date: 06/18/2020
 3um 9um

0th Coef(*100): -15726 -19510
1st Coef(*100): 296627 143165
2nd Coef(*100): -369 494
H2O adj(mg/l*10k): 220 240

***** Printer Test End *****



Operator Signature
CHARLES EDER

Remarks:

PRINT TEST

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 15:49

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	15:50
02 Std. Sol.	0.030	15:51
03 Room Air	0.000	15:51
04 Std. Sol.	0.030	15:52
05 Room Air	0.000	15:52
06 Std. Sol.	0.029	15:53
07 Room Air	0.000	15:54

08 Sim Temp = 34.0°C

Simul Ser No = MP5289

Std Sol No = 201911E

County = 08

Oper No. = 666666



Operator Signature

CHARLES EDER

Remarks:

LOW AC

*~~0.30~~ 0.030 AC
CEE*

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:02

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:03
02 Std. Sol.	0.295	16:04
03 Room Air	0.000	16:04
04 Std. Sol.	0.296	16:05
05 Room Air	0.000	16:05
06 Std. Sol.	0.296	16:06
07 Room Air	0.000	16:07


08 Sim Temp = 34.0°C

Simul Ser No = MP3069

Std Sol No = 20030

County = 08

Oper No. = 666666

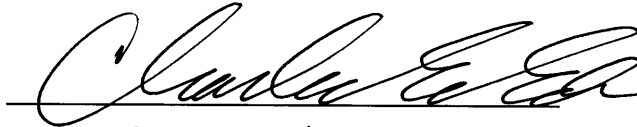

Operator Signature

CHARLES EDER

Remarks:

HIGH AK
0.300 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:09

Test	AC	Time
01 Room Air	0.000	16:10
02 *Subject Test	INT*	16:10
03 Room Air	0.000	16:11

*Invalid Test
Interferent Detected

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = INTERFERENT
Dr. Lic. = ND/DIS921456
Lot No = 05620080A1
Cyl No = 6
Expiration Date = 04/05/2022
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

Interferent Check
0.10AC + 0.05% ACETONE

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:11

Test	AC	Time
01 Diagnostic	OK	16:12
02 Room Air	0.000	16:13
03 *Subject Test	RFI*	16:13
04 Room Air	0.000	16:13

*Invalid Test
Inhibited - RFI

Sub Name = DISCOVER, THE SPIRIT
Sub DOB = 02/01/1992
Sub Sex = Female Weight = 150
Test = DUI Cit = RFI CHECK
Dr. Lic. = ND/DIS921456
Lot No = 05620080A1
Cyl No = 6
Expiration Date = 04/05/2022
County = 08 Oper No. = 666666

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature
CHARLES EDER

Remarks: RFI CHECK

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:18

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:19
02 Std. Gas	0.080	16:19
03 Room Air	0.000	16:19
04 Std. Gas	0.079	16:20
05 Room Air	0.000	16:20
06 Std. Gas	0.079	16:21
07 Room Air	0.000	16:21

Lot No = 05620080A1
Cyl No = 6
Exp Date = 04/05/2022
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

CALIBRATION CHECK
0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:22

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:22
02 Std. Gas	0.079	16:22
03 Room Air	0.000	16:23
04 Std. Gas	0.080	16:23
05 Room Air	0.000	16:24
06 Std. Gas	0.079	16:24
07 Room Air	0.000	16:25

Lot No = 05620080A1
Cyl No = 6
Exp Date = 04/05/2022
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

CALIBRATION CHECK
0.080 AC

Form 106-I8000

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005939
Location = TOXL 8164.14.00 09/16
08/20/2021 16:25

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:26
02 Std. Gas	0.080	16:26
03 Room Air	0.000	16:27
04 Std. Gas	0.079	16:27
05 Room Air	0.000	16:27
06 Std. Gas	0.079	16:28
07 Room Air	0.000	16:28

Lot No = 05620080A1
Cyl No = 6
Exp Date = 04/05/2022
County = 08 Oper No. = 666666



Operator Signature
CHARLES EDER

Remarks:

CALIBRATION CHECK
0.080 AC

Form 106-I8000