



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59281 (06/2018)

Serial Number 80-004201	Instrument Location CARS
Reason for Install/Repair	
<input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level 1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): Exchange
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature 	Date 11-22-2022
-------------------------------	---------------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By 	Certified Date 13 Dec 2022
---------------------------	--------------------------------------

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501


CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = CARS 8164.14.00 09/16
11/22/2022 14:28

DRY CAL CHECK

Test	AC	Time
01 Room Air	RFI*	14:29
02 Room Air	0.000	14:29

*Invalid Test
Inhibited - RFI

Lot No = 26021080A1
Cyl No = 6
Exp Date = 10/05/2023
County = 19 Oper No. = 132318

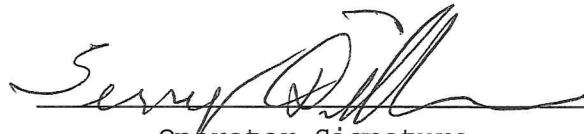

Operator Signature
TERRY DILLON

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = CARS 8164.14.00 09/16
11/22/2022 14:23

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:23
02 Std. Gas	0.079	14:24
03 Room Air	0.000	14:24
04 Std. Gas	0.080	14:25
05 Room Air	0.000	14:25
06 Std. Gas	0.080	14:26
07 Room Air	0.000	14:26

Lot No = 26021080A1
Cyl No = 6
Exp Date = 10/05/2023
County = 19 Oper No. = 132318


Operator Signature
TERRY DILLON

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004201
Location = CARS 8164.14.00 09/16
11/22/2022 14:10

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=_|
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 26021080A1
Standard Cyl #? 6
Standard Expiration? 10/05/2023
Oper No? 132318

Flow Cal. Date: 04/25/2018
Slope 580
Intercept -164747

IR Calibration Date: 04/26/2018

	3um	9um
0th Coef(*100):	-26833	-28328
1st Coef(*100):	272526	132392
2nd Coef(*100):	3966	1397
H2O adj (mg/l*10k):	561	514

***** Printer Test End *****

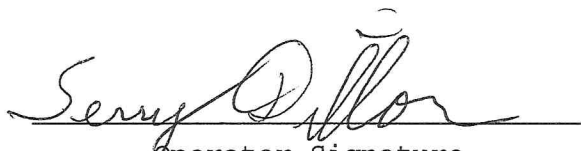
Operator Signature
TERRY DILLON

Remarks:



INTOXILYZER® RECORD
ND OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION
SFN 50496 (05/2018)

Intoxilyzer® Serial Number <i>80-004201</i>		Location <i>CARS</i>
Gas Lot Number <i>26021080 A1</i>	Gas Cylinder Number <i>6</i>	Gas Expiration Date <i>10/05/23</i>

Test Date	Chemical Test Operator Number	Operator's Name (PRINT Last Name, First, MI)
<i>11-22-22</i>	<i>132318</i>	<i>Dillon, Terry Print Test</i>
<i>11-22-22</i>	<i>132318</i>	<i>Dillon, Terry Calibration</i>
<i>11-22-22</i>	<i>132318</i>	<i>Dillon, Terry RFI</i>