



ETHANOL GAS STANDARD CYLINDER REPORT
 OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION
 Toxicology Section/Breath Alcohol Program
 SFN 59282 (10-2013)

Chemical Test Operator Name (Print) <i>Melby, Robert M.</i>		
Location <i>Cross Divide Co. S.O.</i>	Intoxilyzer® Serial Number <i>80-005957</i>	
Gas Lot Number <i>19615080A5</i>	Gas Cylinder Number <i>23</i>	Gas Expiration Date <i>9/05/2017</i>

Check When Done:

- 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 2. Perform an ACA Test (Level 1, Function C).
 Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 A. *0.078* AC
 B. *0.078* AC
 C. *0.079* AC
- 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

Chemical Test Operator Signature <i>[Signature]</i> #4242	Date <i>06-20-2016</i> ✓
Reviewed By (Crime Laboratory Use Only) <i>[Signature]</i>	Date <i>23 June 16</i>


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005957
Location = CROS 8164.13.00 06/09
06/20/2016 16:40

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:41
02 Std. Gas	0.078	16:41
03 Room Air	0.000	16:42
04 Std. Gas	0.078	16:42
05 Room Air	0.000	16:43
06 Std. Gas	0.079	16:43
07 Room Air	0.000	16:43

Lot No = 19615080A5
Cyl No = 23
Exp Date = 09/05/2017
County = 12 Oper No. = 100011

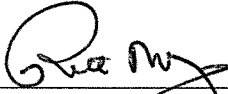 #4242


Operator Signature
ROBERT MELBY

Remarks:

Test OK

Form 106-I8000

I, undersigned, having custody of the original record, certifies that the contents of this true and correct copy of document maintained as part of the files of this office.

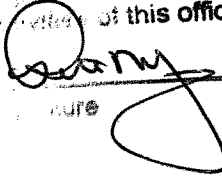

Signature

6-20-16
Date