



NORTH DAKOTA OFFICE OF ATTORNEY GENERAL  
CRIME LABORATORY DIVISION

INTOXILYZER® 8000 ANNUAL INSPECTION

Intoxilyzer® 8000 Serial Number: 80-00 499 Inspection Location: TOXL

A. Pre-Inspection

1. Items with Instrument:
  - Gas Cylinder Yes or No (If Yes, Lot # \_\_\_\_\_ Cyl. # \_\_\_\_\_)
  - Keys Yes or No
  - Power Cord Yes or No
2.  Download Data
3.  Upload Operator File
4.  Current Location Code: FAPD
5.  Battery Check
  - Was the external battery pack replaced? Yes or No
6.  O-Rings
  - Replaced Simulator O-Ring Yes or No
  - Replaced Breath Tube O-Ring Yes or No

B. General Setup and Checks:

1.  Diagnostics passed and instrument in "Ready" mode
2.  Breath tube heated
3.  Date, time and location code (Level 2,E). Re-set if necessary.
  - Time Zone: CST or MDT (Time on test records will be in time zone circled).
4.  Print test (Level 1,P). Sign and attach test record.
5.  Tank monitor (Level 3,D,G).
  - Display: 902 psi Regulator: 900 psi
  - Display and Regulator ± 50 psi of each other Yes or No
  - Gas tank tare necessary? Yes or No
  - If Yes, display readings after tare (Level 3,M,C,G):
  - Display: \_\_\_\_\_ psi Regulator: \_\_\_\_\_ psi

C. Tests (Sign and attach test records):

1.  Configure simulator for the following test (Level 1,S).
  - Wet Calibration Check - Low AC (Level 1,C)
  - Known Value ≤ 0.03 AC: 0.020 AC
  - Sim. Ser #: 20 MP5318
  - Lot #: 202204
  - Exp. Date: 1Jan24
  - Results ± 0.005 of known AC

Toxicology Unit - Breath Alcohol Section	Qualtrax ID: 11698 Revision: 1
Intoxilyzer 8000 Annual Inspection	Status: Published Published: 03/24/2023
UNCONTROLLED WHEN PRINTED	Page 1 of 2



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer           Alcohol Analyzer  
North Dakota Model 8000       SN 80-004199  
Location = TOXL                8164.14.00 09/16  
05/18/2023                      11:38

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxy1234567890-=\_|  
ABCDEFGHIJKLMN0PQRSTUVWXYZ!@#%&^\*()\_+?

abcdefghijklmnopqrstuvwxy1234567890-=\_|  
ABCDEFGHIJKLMN0PQRSTUVWXYZ!@#%&^\*()\_+?

Current Instrument Setup

Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?         No  
Display Volume?             No  
Disable On Memfull?         Yes  
# of Print Copies?          1  
Select Std (D/W/I)?         Dry  
Standard Value?             0.080  
Standard Lot #?             02621080A1  
Standard Cyl #?             13  
Standard Expiration?        03/05/2023  
Oper No?                    133237

Flow Cal. Date:             06/17/2019  
Slope                        593  
Intercept                    -481610

IR Calibration Date:        06/17/2019  
                              3um                    9um

-----  
0th Coef(\*100):            -17512            -19237  
1st Coef(\*100):            277207            136841  
2nd Coef(\*100):            1620               1027  
H2O adj(mg/l\*10k):        211                 375

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
ANNA NAREHOOD

Remarks: Print Test

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004199  
Location = TOXL      8164.14.00 09/16  
05/18/2023      11:55

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:56
02 Std. Sol.	0.022	11:57
03 Room Air	0.000	11:57
04 Std. Sol.	0.020	11:58
05 Room Air	0.000	11:59
06 Std. Sol.	0.022	11:59
07 Room Air	0.000	12:00


08 Sim Temp = 34.0°C

Simul Ser No = MP5318

Std Sol No = 202201A

County = 08

Oper No. = 133237

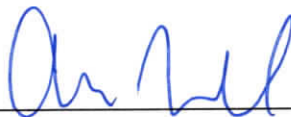

Operator Signature  
ANNA NAREHOOD

Remarks: LOW AC Check-0.020AC

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004199  
Location = TOXL      8164.14.00 09/16  
05/18/2023      12:01

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:01
02 Std. Sol.	0.300	12:02
03 Room Air	0.000	12:03
04 Std. Sol.	0.301	12:03
05 Room Air	0.000	12:04
06 Std. Sol.	0.302	12:05
07 Room Air	0.000	12:05

08 Sim Temp = 34.0°C

Simul Ser No = MP3067

Std Sol No = 202201F

County = 08

Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks:

High Ac Check -0.300Ac

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-004199  
Location = TOXL                      8164.14.00 09/16  
05/18/2023                              12:06

Test	AC	Time
01 Room Air	0.000	12:07
02 *Subject Test	INT*	12:07
03 Room Air	0.000	12:08

\*Invalid Test  
Interferent Detected

Sub Name = TEST, DONOR2 NONE  
Sub DOB = 07/25/1998  
Sub Sex = Male                              Weight = NA  
Test = OTH                              Cit = INTERFERENT CK  
Dr. Lic. = ND/TES989643  
Lot No = 26021080A1  
Cyl No = 3  
Expiration Date = 10/05/2023  
County = 08                              Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks: Interferent Check

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer                      Alcohol Analyzer  
North Dakota Model 8000                      SN 80-004199  
Location = TOXL                              8164.14.00 09/16  
05/18/2023    12:08

Test	AC	Time
01 Diagnostic	OK	12:09
02 Room Air	0.000	12:10
03 *Subject Test	RFI*	12:10
04 Room Air	0.000	12:10

\*Invalid Test  
Inhibited - RFI

Sub Name = TEST, DONOR2 NONE  
Sub DOB = 07/25/1998  
Sub Sex = Male                                      Weight = NA  
Test = OTH    Cit = RFI CHECK  
Dr. Lic. = ND/TES989643  
Lot No = 26021080A1  
Cyl No = 3  
Expiration Date = 10/05/2023  
County = 08    Oper No. = 133237

I followed the Approved Method and the instructions displayed by the Intoxilyzer in conducting this test.



Operator Signature  
ANNA NAREHOOD

Remarks: RFI check

Form 106-I8000


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004199  
Location = TOXL      8164.14.00 09/16  
05/18/2023      12:11

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:12
02 Std. Gas	0.081	12:12
03 Room Air	0.000	12:13
04 Std. Gas	0.081	12:13
05 Room Air	0.000	12:14
06 Std. Gas	0.081	12:14
07 Room Air	0.000	12:15

Lot No = 26021080A1  
Cyl No = 3  
Exp Date = 10/05/2023  
County = 08      Oper No. = 133237

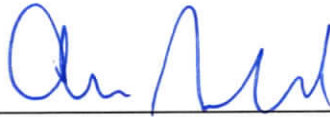

Operator Signature  
ANNA NAREHOOD

Remarks: Dry Cal check - 0. #1-3

Form 106-I8000



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004199  
Location = TOXL      8164.14.00 09/16  
05/18/2023      12:15

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:16
02 Std. Gas	0.082	12:16
03 Room Air	0.000	12:16
04 Std. Gas	0.082	12:17
05 Room Air	0.000	12:17
06 Std. Gas	0.082	12:18
07 Room Air	0.000	12:18

Lot No = 26021080A1  
Cyl No = 3  
Exp Date = 10/05/2023  
County = 08      Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks: Dry Cal Check #4-6

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004199  
Location = TOXL      8164.14.00 09/16  
05/18/2023      12:18

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:19
02 Std. Gas	0.081	12:19
03 Room Air	0.000	12:20
04 Std. Gas	0.081	12:20
05 Room Air	0.000	12:21
06 Std. Gas	0.081	12:21
07 Room Air	0.000	12:22

Lot No = 26021080A1  
Cyl No = 3  
Exp Date = 10/05/2023  
County = 08      Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks: Dry Cal check #7-9

Form 106-I8000