



ETHANOL GAS STANDARD CYLINDER REPORT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59282 (08/2018)

Chemical Test Operator Name (Print) <i>Allen Kathryn</i>		
Location <i>Cass County Jail</i>	Intoxilyzer® Serial Number <i>80-005944</i>	
Gas Lot Number <i>34921080A1</i>	Gas Cylinder Number <i>50</i>	Gas Expiration Date <i>2/5/24</i>

Check When Done:

- 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 2. Perform an ACA Test (Level 1, Function C).
 Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 A. 0. *082* AC
 B. 0. *081* AC
 C. 0. *082* AC
- 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G)
 - B. ACA Test Record.

Chemical Test Operator Signature <i>Kathryn Allen 367H</i>	Date <i>5-26-23</i>
Reviewed By (Crime Laboratory Use Only) <i>Aime Nahl</i>	Date <i>26May2023</i>

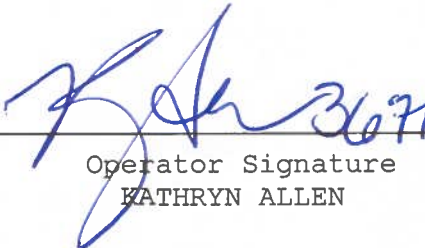
Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005944
Location = CASS 8164.14.00 09/16
05/26/2023 08:22

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	08:22
02 Std. Gas	0.082	08:23
03 Room Air	0.000	08:23
04 Std. Gas	0.081	08:24
05 Room Air	0.000	08:24
06 Std. Gas	0.082	08:25
07 Room Air	0.000	08:25

Lot No = 34921080A1
Cyl No = 50
Exp Date = 02/05/2024
County = 09 Oper No. = 090801


Operator Signature
KATHRYN ALLEN

Remarks:

Acceptable Range!
Form 106-I8000