



NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION

INTOXILYZER® 8000 REPAIR AND MAINTENANCE

Intoxilyzer® 8000 Serial Number: 80-00 4939 Maintenance Location: TOXL

1. Reported Issue and Issue Found:

High dry gas calibration values.

2. Repair/Maintenance Performed:

Calibration adjustment.

3. Does the instrument flow sensor need to be adjusted? Yes or No

- If Yes, complete Flow Sensor Adjustment and Calibration (Qualtrax ID: 11872), then go to step 4.
- If No, go to step 4.

4. Does the instrument calibration need to be adjusted? Yes or No

- If Yes, complete Intoxilyzer® 8000 Calibration Adjustment (Qualtrax ID: 11859), then go to step 5.
- If No, go to step 5.

5. Was the instrument cover removed for the repair? Yes or No

- If Yes, complete Intoxilyzer® 8000 Initial Inspection (Qualtrax ID: 11784), steps 6-7 are not needed, and sign below.
- If No, go to step 6.

6. Is an Annual Inspection due on the Intoxilyzer® 8000? Yes or No

- If Yes, complete Intoxilyzer® 8000 Annual Inspection (Qualtrax ID: 11698), step 7 is not needed, and sign below.
- If No, go to step 7.

Intoxilyzer 8000 Repair and Maintenance

Laboratory Unit: Toxicology Unit - Breath Alcohol Section

Approved By: Laboratory Director

UNCONTROLLED WHEN PRINTED

Qualtrax ID: 11860 Revision: 2

Status: Published

Date Approved: 06/21/2023

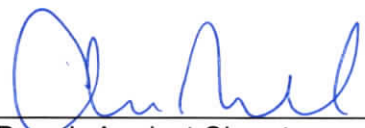
Page 1 of 2

7. Complete Intoxilyzer® 8000 Calibration (Qualtrax ID: 11871) and sign below.

Remarks/Notes: NIA

Instrument is acceptable to be used in the field. Yes or No

If No, state reason(s) why:


Breath Analyst Signature

22 June 2023
Date


Reviewer

22 June 2023
Date