




**INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT**  
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL  
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM  
 SFN 59281 (06/2018)

Serial Number <b>80-004944</b>	Instrument Location <b>Will</b>
Reason for Install/Repair	
<input type="checkbox"/> Install After Receiving From Crime Laboratory <input checked="" type="checkbox"/> Install After Location Change <input checked="" type="checkbox"/> Other (Specify) <b>Moved it</b>	

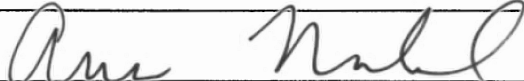
Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
  - A. Print Test (Level1, Function P).
  - B. ACA Test (Level 1, Function C).
  - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): **N/A**
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature 	Date <b>8-3-23</b>
--	-----------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By 	Certified Date <b>04Aug2023</b>
--	------------------------------------

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000            SN 80-004944  
Location = WILL                    8164.14.00 09/16  
08/03/2023                            16:28

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-=\_|  
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#\$\$%^&\*()\_+?

Current Instrument Setup

Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?        No  
Display Volume?            No  
Disable On Memfull?        Yes  
# of Print Copies?          1  
Select Std (D/W/I)?        Dry  
Standard Value?            0.080  
Standard Lot #?            26021080A1  
Standard Cyl #?            13  
Standard Expiration?       10/05/2023  
Oper No?                    130739

Flow Cal. Date:            07/01/2011  
Slope                        685  
Intercept                   -634740

IR Calibration Date:       07/01/2011  
                                  3um                    9um

-----  
0th Coef(\*100):            -22591                -17234  
1st Coef(\*100):            265915                132691  
2nd Coef(\*100):            2841                    1562  
H2O adj(mg/l\*10k):        559                    425

\*\*\*\*\* Printer Test End \*\*\*\*\*

  
Operator Signature  
MICHELLE ROMANS

Remarks:

Print test

Form 106-I8000

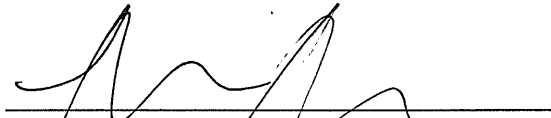
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-004944  
Location = WILL      8164.14.00 09/16  
08/03/2023      16:29

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	16:29
02 Std. Gas	0.080	16:30
03 Room Air	0.000	16:30
04 Std. Gas	0.080	16:31
05 Room Air	0.000	16:31
06 Std. Gas	0.079	16:31
07 Room Air	0.000	16:32

Lot No = 26021080A1  
Cyl No = 13  
Exp Date = 10/05/2023  
County = 53      Oper No. = 130739

  
Operator Signature  
MICHELLE ROMANS

Remarks:

ACA Test

Form 106-I8000


Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer            Alcohol Analyzer  
North Dakota Model 8000        SN 80-004944  
Location = WILL                 8164.14.00 09/16  
08/03/2023                        16:33

Test	AC	Time
01 Room Air	RFI*	16:34
02 Room Air	0.000	16:34

\*Invalid Test  
Inhibited - RFI

Sub Name = TEST, TEST TEST  
Sub DOB = 01/01/1999  
Sub Sex = Male                    Weight = 200  
Test = DUI                         Cit = 99999  
Dr. Lic. = ND/999999  
Lot No = 26021080A1  
Cyl No = 13  
Expiration Date = 10/05/2023  
County = 53                        Oper No. = 130739



Operator Signature  
MICHELLE ROMANS

Remarks:

RFI TEST

Form 106-I8000