



**INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT**  
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL  
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM  
 SFN 59281 (06/2018)

Serial Number <b>80-006491</b>	Instrument Location <b>W.CSO.</b>
--------------------------------	-----------------------------------

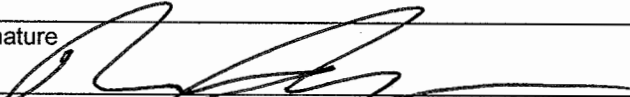
Reason for Install/Repair

Install After Receiving From Crime Laboratory     Install After Location Change

Other (Specify) Annual inspection at crime Laboratory

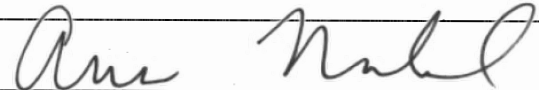
Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
  - A. Print Test (Level1, Function P).
  - B. ACA Test (Level 1, Function C).
  - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): N/A
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature 	Date <b>05-Dec-23</b>
---	-----------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By 	Certified Date <b>06Dec2023</b>
---	---------------------------------

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer           Alcohol Analyzer  
North Dakota Model 8000        SN 80-006491  
Location = WESO                 8164.14.00 09/16  
12/05/2023                       12:35

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxyz1234567890-|=|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

abcdefghijklmnopqrstuvwxyz1234567890-|=|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#\$\$%^&\*()\_+?

Current Instrument Setup

Data Entry Mode:                Enabled  
Start Test Sequence:            DABACABA  
Display Prelim Rslt?            Yes  
Display Third Digit?            Yes  
Inhib Printer(Y/N)?            No  
Display Volume?                 No  
Disable On Memfull?            Yes  
# of Print Copies?              1  
Select Std (D/W/I)?            Dry  
Standard Value?                 0.080  
Standard Lot #?                 14323080A1  
Standard Cyl #?                 19  
Standard Expiration?            06/05/2025  
Oper No?                         131495

Flow Cal. Date:                 05/22/2018  
Slope                            639  
Intercept                       -618311

IR Calibration Date:            05/23/2018

	3um	9um
0th Coef(*100):	-23641	-17066
1st Coef(*100):	264855	137703
2nd Coef(*100):	2663	1471
H2O adj(mg/l*10k):	556	451

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
MICHAEL BEAVER

Remarks:

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-006491  
Location = WESO      8164.14.00 09/16  
12/05/2023      14:07

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	14:07
02 Std. Gas	0.082	14:07
03 Room Air	0.000	14:08
04 Std. Gas	0.082	14:08
05 Room Air	0.000	14:09
06 Std. Gas	0.082	14:09
07 Room Air	0.000	14:10

Lot No = 14323080A1  
Cyl No = 19  
Exp Date = 06/05/2025  
County = 52      Oper No. = 131495



Operator Signature  
MICHAEL BEAVER

Remarks:

Form 106-I8000

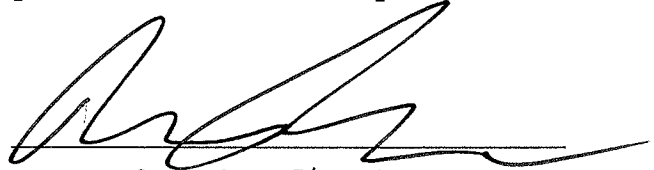
Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-006491  
Location = WESO      8164.14.00 09/16  
12/05/2023      14:10

Test	AC	Time
01 Room Air	RFI*	14:11
02 Room Air	0.000	14:12

\*Invalid Test  
Inhibited - RFI

Sub Name = TEST, TEST TEST  
Sub DOB = 01/01/1999  
Sub Sex = Male      Weight = 120  
Test = DUI      Cit = 123  
Dr. Lic. = ND/123  
Lot No = 14323080A1  
Cyl No = 19  
Expiration Date = 06/05/2025  
County = 52      Oper No. = 131495



Operator Signature  
MICHAEL BEAVER

Remarks:

Form 106-I8000