



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59281 (06/2018)

Serial Number <i>80-003067</i>	Instrument Location <i>Grafton PD</i>
Reason for Install/Repair	
<input type="checkbox"/> Install After Receiving From Crime Laboratory <input checked="" type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>David Nishro</i>	Date <i>1/26/2024</i>
--	--------------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By <i>Aime Nahl</i>	Certified Date <i>29Jan2024</i>
---	------------------------------------

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003067
Location = GFTN 8164.14.00 09/16
01/26/2024 12:04

***** Printer Test *****

abcdefghijklmnopqrstuvwxy1234567890-|=|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^*()_+?

abcdefghijklmnopqrstuvwxy1234567890-|=|
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 01923080A3
Standard Cyl #? 16
Standard Expiration? 02/05/2025
Oper No? 132597

Flow Cal. Date: 04/17/2020
Slope 609
Intercept -236750

IR Calibration Date: 04/17/2020
 3um 9um

0th Coef(*100): -19307 -18350
1st Coef(*100): 282277 142540
2nd Coef(*100): 1257 744
H2O adj(mg/l*10k): 302 214

***** Printer Test End *****

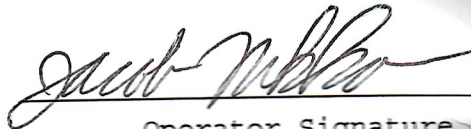

Operator Signature
JACOB NIKKO

Remarks:

Form 106-I8000

The undersigned, having custody of the original public record, certifies that the information hereon is a true and correct copy of the original document maintained as a part of the activities of this agency.

Signature/Title *Jacob Nikko / Police Officer*
Agency/Date *Crafton PD / 1/26/2024*


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003067
Location = GFTN 8164.14.00 09/16
01/26/2024 12:05

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	12:05
02 Std. Gas	0.080	12:06
03 Room Air	0.000	12:06
04 Std. Gas	0.080	12:07
05 Room Air	0.000	12:07
06 Std. Gas	0.080	12:08
07 Room Air	0.000	12:08

Lot No = 01923080A3
Cyl No = 16
Exp Date = 02/05/2025
County = 50 Oper No. = 132597



Operator Signature
JACOB NIKKO

Remarks:

Form 106-I8000

The undersigned, having custody of the original public record, certifies that the information hereon is a true and correct copy of the original document maintained as a part of the activities of this agency.

Signature/Title Jacob Nikko / Police Officer
Agency/Date Grafton PD / 1/25/24


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-003067
Location = GFTN 8164.14.00 09/16
01/26/2024 12:10

Test	AC	Time
01 Room Air	RFI*	12:11
02 Room Air	0.000	12:12

*Invalid Test
Inhibited - RFI

Sub Name = TEST, RFI NA
Sub DOB = 11/11/2000
Sub Sex = Male Weight = 180
Test = OTH Cit = TEST
Dr. Lic. = ND/TEST
Lot No = 01923080A3
Cyl No = 16
Expiration Date = 02/05/2025
County = 50 Oper No. = 132597

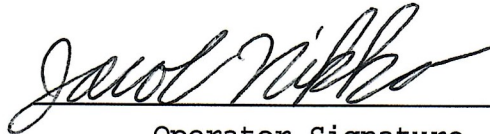

Operator Signature
JACOB NIKKO

Remarks:

Form 106-I8000

The undersigned, having custody of the original public record, certifies that the information hereon is a true and correct copy of the original document maintained as a part of the activities of this agency.

Signature/Title Jacob Nikko / Police Officer

Agency/Date Grafton PD / 1/26/24