



NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
CRIME LABORATORY DIVISION

INTOXILYZER® 8000 REPAIR AND MAINTENANCE

Intoxilyzer® 8000 Serial Number: 80-00 5937 Maintenance Location: TOXL

1. Reported Issue and Issue Found:

Modem port was pushed into the Intoxilyzer.

2. Repair/Maintenance Performed:

Opened Intoxilyzer to reposition modem port.
Replaced the external snap-on part to the modem.
Started the initial inspection but the pump wasn't working
correctly. Replaced the pump + performed a new initial inspection.

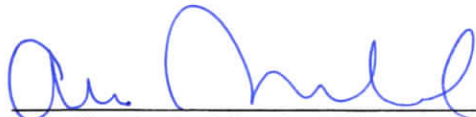
3. Does the instrument flow sensor need to be adjusted? Yes or No
- If Yes, complete Flow Sensor Adjustment and Calibration (Document ID: 11872), then go to step 4.
 - If No, go to step 4.
4. Does the instrument calibration need to be adjusted? Yes or No
- If Yes, complete Intoxilyzer® 8000 Calibration Adjustment (Document ID: 11859), then go to step 5.
 - If No, go to step 5.
5. Was the instrument cover removed for the repair? Yes or No
- If Yes, complete Intoxilyzer® 8000 Initial Inspection (Document ID: 11784), then go to step 6.
 - If No, go to step 6.
6. Is an Annual Inspection due on the Intoxilyzer® 8000? Yes or No
- If Yes, complete Intoxilyzer® 8000 Annual Inspection (Document ID: 11698), step 7 is not needed, and sign below.
 - If No, go to step 7.

7. Was an Intoxilyzer® 8000 Initial Inspection (Document ID: 11784) completed in step 5? Yes or No
- If Yes, no further action is needed and sign below.
 - If No, Complete Intoxilyzer® 8000 Calibration (Document ID: 11871) and sign below.

Remarks/Notes: NIA

Instrument is acceptable to be used in the field. Yes or No

If No, state reason(s) why:


Breath Alcohol Analyst Signature

15 May 2024
Date


Reviewer Signature

16 May 2024
Date