



NORTH DAKOTA OFFICE OF ATTORNEY GENERAL  
CRIME LABORATORY DIVISION

INTOXILYZER® 8000 ANNUAL INSPECTION

Intoxilyzer® 8000 Serial Number: 80-00 5958 Inspection Location: TOXL

A. Pre-Inspection

1. Items with Instrument:
  - Gas Cylinder Yes or No (If Yes, Lot # \_\_\_\_\_ Cyl. # \_\_\_\_\_)
  - Keys Yes or No
  - Power Cord Yes or No
2.  Download Data
3.  Upload Operator File
4.  Current Location Code: 011V
5.  Battery Check
  - Was the external battery pack replaced? Yes or No
6.  O-Rings
  - Replaced Simulator O-Ring Yes or No
  - Replaced Breath Tube O-Ring Yes or No

B. General Setup and Checks:

1.  Diagnostics passed and instrument in "Ready" mode
2.  Breath tube heated
3.  Date, time and location code (Level 2,E). Re-set if necessary.  
Time Zone: CST or MDT (Time on test records will be in time zone circled).
4.  Print test (Level 1,P). Sign and attach test record.
5.  Tank monitor (Level 3,D,G).  
Display: 974 psi Regulator: 1000 psi  
Display and Regulator ± 50 psi of each other Yes or No  
Gas tank tare necessary? Yes or No  
If Yes, display readings after tare (Level 3,M,C,G):  
Display: \_\_\_\_\_ psi Regulator: \_\_\_\_\_ psi

C. Tests (Sign and attach test records):

1.  Configure simulator for the following test (Level 1,S).  
Wet Calibration Check - Low AC (Level 1,C)  
Known Value ≤ 0.03 AC: 0.020 AC  
Sim. Ser #: MP3002  
Lot #: 202301E  
Exp. Date: 17Jan25  
 Results ± 0.005 of known AC



Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer           Alcohol Analyzer  
North Dakota Model 8000       SN 80-005958  
Location = TOXL                8164.14.00 09/16  
06/25/2024                      09:42

\*\*\*\*\* Printer Test \*\*\*\*\*

abcdefghijklmnopqrstuvwxy1234567890-|=|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

abcdefghijklmnopqrstuvwxy1234567890-|=|  
ABCDEFGHIJKLMNopqrstuvwxyz!@#%&^\*()\_+?

Current Instrument Setup

Data Entry Mode:            Enabled  
Start Test Sequence:        DABACABA  
Display Prelim Rslt?        Yes  
Display Third Digit?        Yes  
Inhib Printer(Y/N)?         No  
Display Volume?             No  
Disable On Memfull?         Yes  
# of Print Copies?           1  
Select Std (D/W/I)?         Dry  
Standard Value?             0.080  
Standard Lot #?             28423080A3  
Standard Cyl #?             41  
Standard Expiration?        11/05/2025  
Oper No?                    133237

Flow Cal. Date:             06/21/2013  
Slope                        716  
Intercept                   -760436

IR Calibration Date:        04/07/2015  
                                  3um                    9um  
-----  
0th Coef(\*100):            -17179            -17694  
1st Coef(\*100):            259328            130890  
2nd Coef(\*100):            3473                1654  
H2O adj(mg/l\*10k):        425                 422

\*\*\*\*\* Printer Test End \*\*\*\*\*



Operator Signature  
ANNA NAREHOOD

Remarks: Print Test

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005958  
Location = TOXL      8164.14.00 09/16  
06/25/2024      09:48

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	09:48
02 Std. Sol.	0.019	09:51
03 Room Air	0.000	09:51
04 Std. Sol.	0.020	09:52
05 Room Air	0.000	09:52
06 Std. Sol.	0.020	09:53
07 Room Air	0.000	09:54

08 Sim Temp = 34.0°C

Simul Ser No = MP3002

Std Sol No = 202301E

County = 08

Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks: LOW AC Check 0.020AE

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005958  
Location = TOXL      8164.14.00 09/16  
06/25/2024      10:45

WET CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:46
02 Std. Sol.	0.393	10:47
03 Room Air	0.000	10:48
04 Std. Sol.	0.394	10:48
05 Room Air	0.000	10:49
06 Std. Sol.	0.394	10:50
07 Room Air	0.000	10:50


08 Sim Temp = 34.0°C

Simul Ser No = MP6036

Std Sol No = 202212C

County = 08

Oper No. = 133237

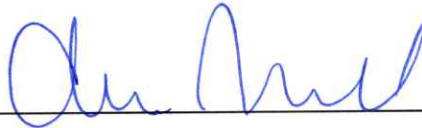

Operator Signature  
ANNA NAREHOOD

Remarks: High AC Check -0.400 AC

Form 106-I8000






Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005958  
Location = TOXL      8164.14.00 09/16  
06/25/2024      10:28

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:29
02 Std. Gas	0.079	10:29
03 Room Air	0.000	10:30
04 Std. Gas	0.079	10:30
05 Room Air	0.000	10:30
06 Std. Gas	0.079	10:31
07 Room Air	0.000	10:31

Lot No = 01923080A3  
Cyl No = 18  
Exp Date = 02/05/2025  
County = 08      Oper No. = 133237

  
\_\_\_\_\_  
Operator Signature  
ANNA NAREHOOD

Remarks: Dry cal check #1-3

Form 106-I8000




Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005958  
Location = TOXL      8164.14.00 09/16  
06/25/2024      10:32

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:33
02 Std. Gas	0.080	10:33
03 Room Air	0.000	10:34
04 Std. Gas	0.079	10:34
05 Room Air	0.000	10:34
06 Std. Gas	0.079	10:35
07 Room Air	0.000	10:35

Lot No = 01923080A3  
Cyl No = 18  
Exp Date = 02/05/2025  
County = 08      Oper No. = 133237

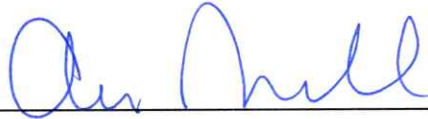

Operator Signature  
ANNA NAREHOOD

Remarks: Dry cal check #4-6

Form 106-I8000

Intoxilyzer Test Record and Checklist  
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer      Alcohol Analyzer  
North Dakota Model 8000      SN 80-005958  
Location = TOXL      8164.14.00 09/16  
06/25/2024      10:36

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:36
02 Std. Gas	0.080	10:37
03 Room Air	0.000	10:37
04 Std. Gas	0.080	10:37
05 Room Air	0.000	10:38
06 Std. Gas	0.079	10:38
07 Room Air	0.000	10:39

Lot No = 01923080A3  
Cyl No = 18  
Exp Date = 02/05/2025  
County = 08      Oper No. = 133237



Operator Signature  
ANNA NAREHOOD

Remarks: Dry Cal check #7-9

Form 106-I8000