

ETHANOL GAS STANDARD CYLINDER REPORT NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM SFN 59282 (08/2018)

Chemical Test Operator Name (Print) Jacob Nikko 132547	1		
Location GRAF	ļ.	ntoxilyzer® Serial Nur	nber 80-00 5357
Gas Lot Number 01529080A4	Gas Cylinder Number 3		Gas Expiration Date 03/05/2026

Check When Done:

- ✓ 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 2. Perform an ACA Test (Level 1, Function C).

Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).

- A. 0.<u>0 \$2</u> AC
- B. 0.<u>∅ ⊉ 2</u> AC
- C. 0.0 \$ 2 AC

3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.

- ☑ 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

Chemical Test Operator Signature	~	2 . 2	Date 1/17/2025
Reviewed By (Crime Laboratory Use Only) Anna Narehood	An	Mall	^{Date} 17Jan2025

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer	Alcohol Analyzer
North Dakota Model 8000	SN 80-005357
Location = GRAF	8164.14.00 09/16
01/17/2025	05:08

				_
			DRY CAL CHECK	
q	lest		۲	Dime
T			AC	Time
01	Room	Air	0.000	05:09
02	Std.	Gas	0.082	05:09
03	Room	Air	0.000	05 : 10
04	Std.	Gas	0.082	05:10
05	Room	Air	Ó.000	05:11
06	Std.	Gas	0.082	05:11
07	Room	Air	0.000	05:12

Lot No = 01524080A4 Cyl No = 30 Exp Date = 03/05/2026 County = 50 Oper M

Oper No. = 132597

Operator Signature JACOB NIKKO

Remarks:

Form 106-18000

The Undersigned, having custody of the original public record, certifies that the information hereon is a true and correct copy of the original document maintained as a part of the activities of this agency.

Signature/Title Agency/Date Valsh Lo So