

NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION

INTOXILYZER® 8000 REPAIR AND MAINTENANCE

Intoxilyzer® 8000 Serial Number: 80-00 7086 Maintenance Location: 10x

- 1. Reported Issue and Issue Found: High dy calibration Values during Annual Inspection.
- 2. Repair/Maintenance Performed: Calibration Adjustment.
- 3. Does a Calibration Adjustment need to be performed? Yes or No
 - If Yes, perform a calibration check.

Configure dry gas standard (Level 1,S). Known Value : 0.080 AC Gas Cylinder Lot #: 14323080A4Cylinder #: 42Exp. Date: 405125Dry Calibration Check (Level 1,C) Test 1 0.083 Test 2 0.083 Test 3 0.083

- If No, go to step 4.
- 4. Does the instrument flow sensor need to be adjusted? Yes or (No)
 - If Yes, complete Flow Sensor Adjustment and Calibration (Document ID: 11872), then go to step 5.
 - If No, go to step 5.

Intoxilyzer 8000 Repair and Maintenance Laboratory Unit: Toxicology Unit - Breath Alcohol Section Approved By: Laboratory Director UNCONTROLLED WHEN PRINTED

- 5. Does the instrument calibration need to be adjusted? Yes or No
 - If Yes, complete Intoxilyzer® 8000 Calibration Adjustment (Document ID: 11859), then go to step 6.
 - If No, go to step 6.
- 6. Was the instrument cover removed for the repair? Yes or No
 - If Yes, complete Intoxilyzer® 8000 Initial Inspection (Document ID: 11784), then go to step 7.
 - If No, go to step 7.
- 7. Is an Annual Inspection due on the Intoxilyzer® 8000? Yes or No
 - If Yes, complete Intoxilyzer® 8000 Annual Inspection (Document ID: 11698), step 8 is not needed, and sign below.
 - If No, go to step 8.
- 8. Was an Intoxilyzer® 8000 Initial Inspection (Document ID: 11784) completed in step 6? Yes or No
 - If Yes, no further action is needed and sign below.
 - If No, Complete Intoxilyzer® 8000 Calibration (Document ID: 11871) and sign below.

Remarks/Notes: Found during Annual Inspection

Instrument is acceptable to be used in the field.

es or

No

Date

If No, state reason(s) why:

Breath Alcohol Analyst Signature

eviewer Signature

Intoxilyzer 8000 Repair and Maintenance Laboratory Unit: Toxicology Unit - Breath Alcohol Section Approved By: Laboratory Director UNCONTROLLED WHEN PRINTED

Document ID: 11860 Revision: 5 Status: Published Date Approved: 03/20/2025 Page 2 of 2

1

04June 2025

05 June 2025

Uploaded 10June2025

AEN

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. IntoxilyzerAlcohol AnalyzerNorth Dakota Model 8000SN 80-007086Location = TOXL8164.16.00 09/1804/29/202510:49

	DRY CAL CHECK	
Test	AC	Time
01 Room Air	0.000	10:49
02 Std. Gas	0.083	10:50
03 Room Air	0.000	10:50
04 Std. Gas	0.083	10:51
05 Room Air	0.000	10:51
06 Std. Gas	0.083	10:52
07 Room Air	0.000	10:52

Lot No = 14323080A4 Cyl No = 42 Exp Date = 06/05/2025 County = 08 Oper No. = 133237

Operator Signature ANNA NAREHOOD

Dry Cal check - cal adj. needed. Remarks:

Form 106-18000