

NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION

INTOXILYZER® 8000 REPAIR AND MAINTENANCE

Intoxilyzer® 8000 Serial Number: 80-00 5362 Maintenance Location: ToxL			
1.	Reported Issue and Issue Found: During annual inspection, high dry calibration		
	values were produced.		
2.	Repair/Maintenance Performed: (alibration adjustment		
 3. Does a Calibration Adjustment need to be performed? Yes or No If Yes, perform a calibration check. Configure dry gas standard (Level 1,S). 			
	Known Value : 0.000 AC Gas Cylinder Lot #: 28433080A3 Cylinder #: 4		
	Exp. Date: 11 5125		
	ADry Calibration Check (Level 1,C) Test 1 <u>0.083</u> Test 2 <u>0.082</u> Test 3 <u>0.083</u>		
	If No, go to step 4.		
4.	 Does the instrument flow sensor need to be adjusted? Yes or No If Yes, complete Flow Sensor Adjustment and Calibration (Document ID: 11872), then go to step 5. If No, go to step 5. 		

Intoxilyzer 8000 Repair and Maintenance

Laboratory Unit: Toxicology Unit - Breath Alcohol Section

Approved By: Laboratory Director

UNCONTROLLED WHEN PRINTED

Document ID: 11860 Revision: 5

Status: Published

Date Approved: 03/20/2025

Page 1 of 2

 5. Does the instrument calibration need to be adjusted? If Yes, complete Intoxilyzer® 8000 Calibration A 11859), then go to step 6. If No, go to step 6. 				
 Was the instrument cover removed for the repair? Yes or No If Yes, complete Intoxilyzer® 8000 Initial Inspection (Document ID: 11784), then go to step 7. If No, go to step 7. 				
 7. Is an Annual Inspection due on the Intoxilyzer® 8000? Yes or No If Yes, complete Intoxilyzer® 8000 Annual Inspection (Document ID: 11698), step 8 is not needed, and sign below. If No, go to step 8. 				
 8. Was an Intoxilyzer® 8000 Initial Inspection (Document ID: 11784) completed in step 6? Yes or No If Yes, no further action is needed and sign below. If No, Complete Intoxilyzer® 8000 Calibration (Document ID: 11871) and sign below. 				
Remarks/Notes: NIA				
Instrument is acceptable to be used in the field. Ves or No If No, state reason(s) why:				
Breath Alcohol Analyst Signature Janelle Futschmur Reviewer Signature	05 June 2025 Date 65 June 2025 Date			

Laboratory Unit: Toxicology Unit - Breath Alcohol Section

Approved By: Laboratory Director UNCONTROLLED WHEN PRINTED

Document ID: 11860 Revision: 5

Status: Published

Date Approved: 03/20/2025

Page 2 of 2

Intoxilyzer Test Record and Checklist NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer North Dakota Model 8000 Location = TOXL 8164.14.00 09/16 04/04/2025

SN 80-005362 13:45

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	13:45
02 Std. Gas	0.083	13:46
03 Room Air	0.000	13:46
04 Std. Gas	0.082	13:47
05 Room Air	0.000	13:47
06 Std. Gas	0.083	13:47
07 Room Air	0.000	13:48

Lot No = 28423080A3

Cyl No = 41

Exp Date = 11/05/2025

County = 08

Oper No. = 133237

Operator Signature ANNA NAREHOOD

calibration adjustment nuded.

Form 106-I8000