

## NORTH DAKOTA OFFICE OF ATTORNEY GENERAL CRIME LABORATORY DIVISION

## INTOXILYZER® 8000 REPAIR AND MAINTENANCE

Int	ntoxilyzer® 8000 Serial Number: 80-00 2667 Maintenance Lo	ocation: TOXL
	Reported Issue and Issue Found:  upon Start-up, the instrument Nad no date	
2.	Repair/Maintenance Performed:  Real-Time Clock Wattery was replaced.	
3.	Does a Calibration Adjustment need to be performed? Yes or  • If Yes, perform a calibration check.  □ Configure dry gas standard (Level 1,S).	No
	Known Value :AC Gas Cylinder Lot #: Cylinder #: Exp. Date: Dry Calibration Check (Level 1,C) Test 1 Test 2 Test 3	
	If No, go to step 4.	
4.	<ul> <li>Does the instrument flow sensor need to be adjusted? Yes or</li> <li>If Yes, complete Flow Sensor Adjustment and Calibration 11872), then go to step 5.</li> <li>If No, go to step 5.</li> </ul>	No (Document ID:
into	oxilyzer 8000 Repair and Maintenance	
	boratory Unit: Toxicology Unit - Breath Alcohol Section	cument ID: 11860 Revision: 5 Status: Published

Page 1 of 2

Date Approved: 03/20/2025

Page 1 of 2

**AEN** 

Approved By: Laboratory Director

Uploaded 09June2025

UNCONTROLLED WHEN PRINTED

<ul> <li>Does the instrument calibration need to be adjusted? Yes or No</li> <li>If Yes, complete Intoxilyzer® 8000 Calibration Adjustment (Document ID: 11859), then go to step 6.</li> <li>If No, go to step 6.</li> </ul>		
<ul> <li>Was the instrument cover removed for the repair? Yes or No</li> <li>If Yes, complete Intoxilyzer® 8000 Initial Inspection (Document ID: 11784), then go to step 7.</li> <li>If No, go to step 7.</li> </ul>		
<ul> <li>7. Is an Annual Inspection due on the Intoxilyzer® 8000? Yes or No</li> <li>If Yes, complete Intoxilyzer® 8000 Annual Inspection (Document ID: 11698), step 8 is not needed, and sign below.</li> <li>If No, go to step 8.</li> </ul>		
<ul> <li>8. Was an Intoxilyzer® 8000 Initial Inspection (Document ID: 11784) completed in step 6? Yes or No</li> <li>If Yes, no further action is needed and sign below.</li> <li>If No, Complete Intoxilyzer® 8000 Calibration (Document ID: 11871) and sign below.</li> </ul>		
Remarks/Notes: NA		
Instrument is acceptable to be used in the field. Yes or No		
If No, state reason(s) why:  From previous annual inspections, this instrument is classroom.  Only because the fan motor doesn't turn on.		
Breath Alcohol Analyst Signature  Objune 2025  Date		
Reviewer Signature  07 Sine 2025  Date		

Laboratory Unit: Toxicology Unit - Breath Alcohol Section

Approved By: Laboratory Director

**UNCONTROLLED WHEN PRINTED** 

Document ID: 11860 Revision: 5

Status: Published

Date Approved: 03/20/2025

Page 2 of 2