



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59281 (06/2018)

Serial Number <u>80-005952</u>	Instrument Location <u>Wisher PD</u>
Reason for Install/Repair	
<input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): N/A AEI 10Feb2026
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature	Date <u>2-2-2026</u>
---------------------------	----------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By <u>Anna Ingemansen</u>	Certified Date <u>10Feb2026</u>
--	---------------------------------

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005952
Location = WISH 8164.14.00 09/16
02/02/2026 10:50

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890- = |
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%&^* () _ + ?

abcdefghijklmnopqrstuvwxyz1234567890- = |
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%&^* () _ + ?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 19825080A2
Standard Cyl #? 25
Standard Expiration? 09/05/2027
Oper No? 010912

Flow Cal. Date: 04/16/2021
Slope 657
Intercept -744615

IR Calibration Date: 04/16/2021

	3um	9um
0th Coef(*100):	-24450	-14639
1st Coef(*100):	266154	132500
2nd Coef(*100):	2284	1091
H2O adj(mg/l*10k):	516	323

***** Printer Test End *****

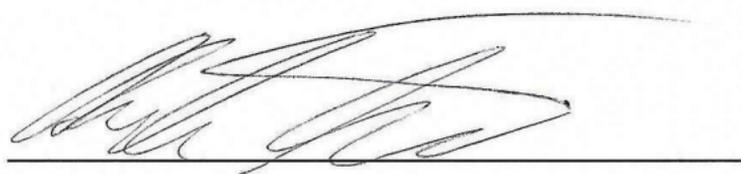

Operator Signature
ANDREW TUNHEIM

Remarks:


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005952
Location = WISH 8164.14.00 09/16
02/02/2026 10:50

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	10:51
02 Std. Gas	0.080	10:51
03 Room Air	0.000	10:52
04 Std. Gas	0.081	10:52
05 Room Air	0.000	10:53
06 Std. Gas	0.080	10:53
07 Room Air	0.000	10:54

Lot No = 19825080A2
Cyl No = 25
Exp Date = 09/05/2027
County = 26 Oper No. = 010912

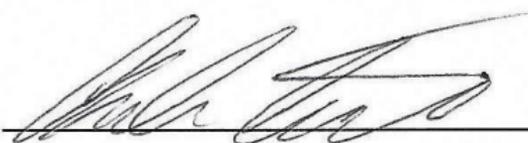

Operator Signature
ANDREW TUNHEIM

Remarks:

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005952
Location = WISH 8164.14.00 09/16
02/02/2026 10:57

DRY CAL CHECK

Test	AC	Time
01 Room Air	RFI*	10:57
02 Room Air	0.000	10:57

*Invalid Test
Inhibited - RFI

Lot No = 19825080A2
Cyl No = 25
Exp Date = 09/05/2027
County = 26 Oper No. = 010912



Operator Signature
ANDREW TUNHEIM

Remarks: